A1A
人体解剖模型 M-100型
Human Anatomy Torso (dual sex)

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
English
Instruction
Manual

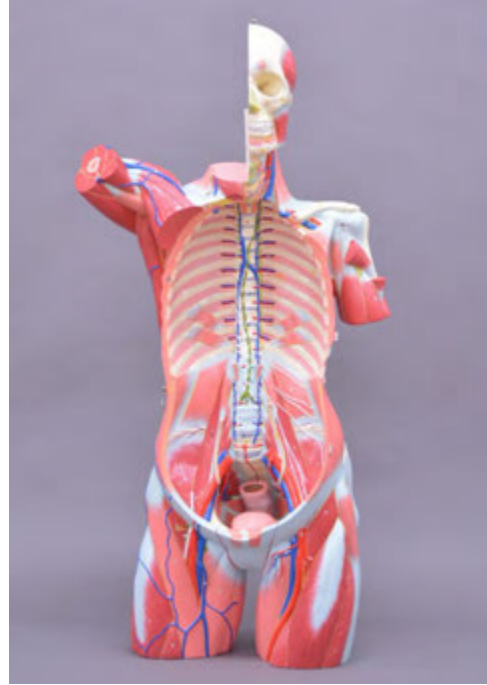
1. Start


2. Lower and fit the reproductive organs. For the female model, please fit the ovaries towards the back. (See point 1)

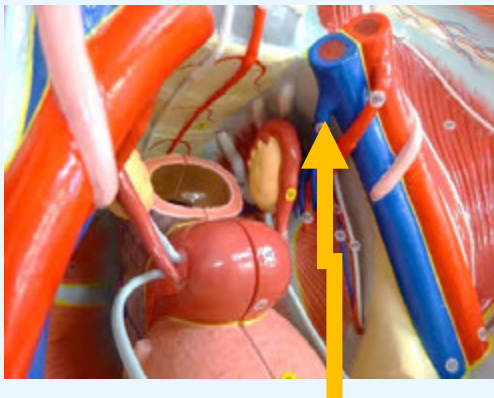Point 1 : For the female model, confirm that the ovaries are attached towards the back before fitting the large intestines.

Behind this blood vessel

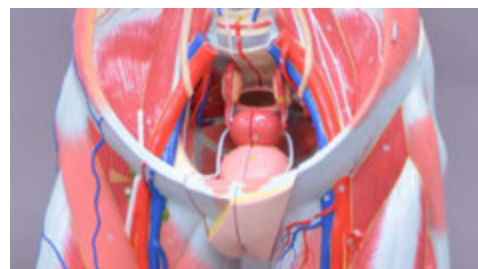▲ Female


▲ Male

3. Attach the lower back of the diaphragm, then the kidneys using the pinholes. (See point 2)

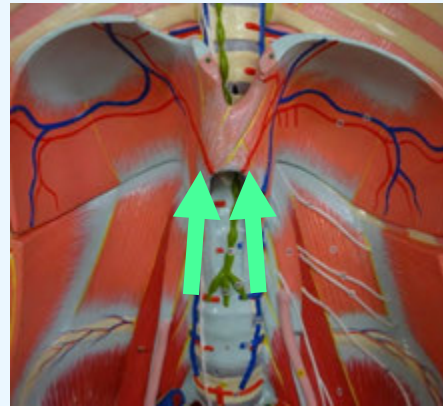
Point 2: Position of the pin on the back of the diaphragm


4. Hang the large intestine on the metal hooks.

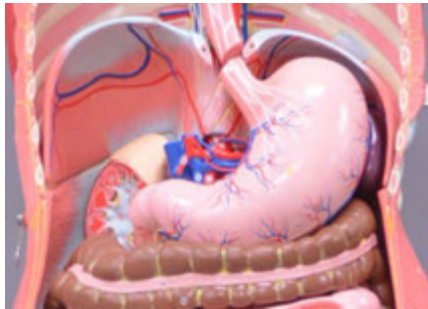5. Attach the small intestine, then esophagus. Place the stomach above the large intestine, matching the with esophagus.

▲ close

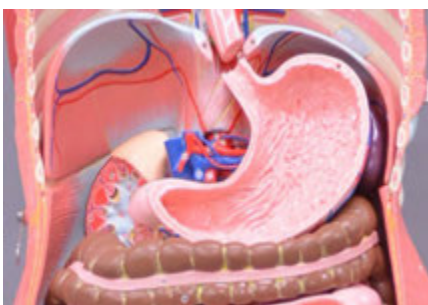▲ open

6. Place the liver on the stomach and the large intestine.

7. Attach the lower front of the diaphragm using the pin.


8. Momentarily remove the esophagus and insert the back of the lungs. Reposition the esophagus after.


9. Attach the front of the lungs to the back of lungs using the pinholes.

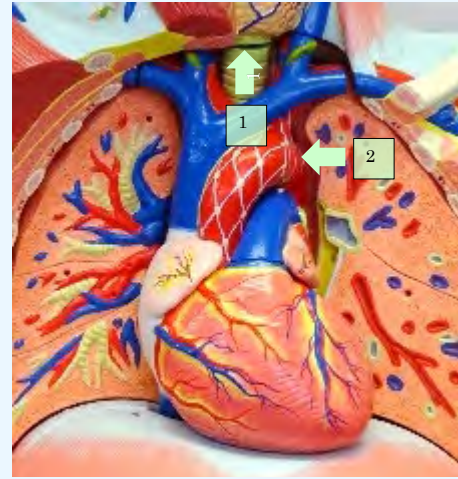
Point 3: For the heart, insert the upper blood vessels towards the throat first, then the aorta via the pinholes. (Fits like a groove)


10. Attach the front of the lungs to the back of lungs using the pinholes.

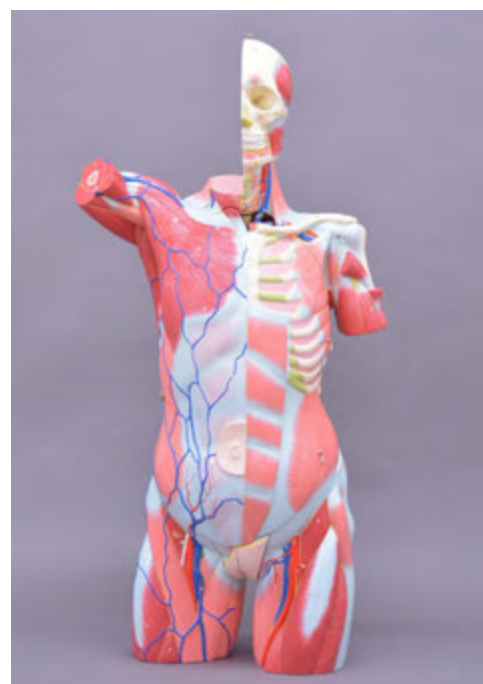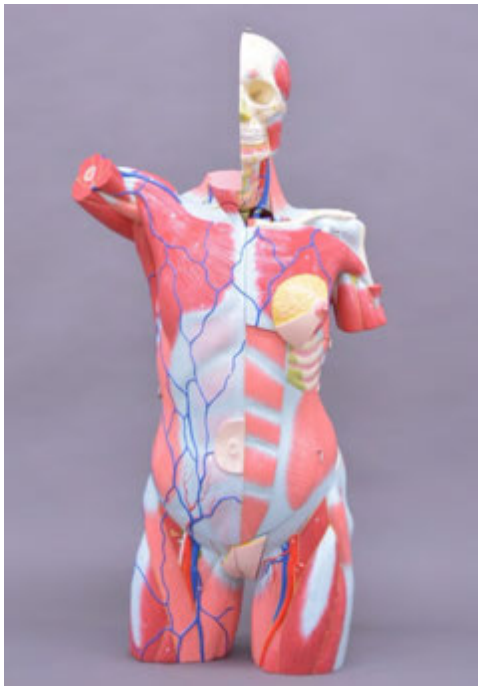
11. Place the abdominal cover over.


12. Attach the muscle piece and breast.

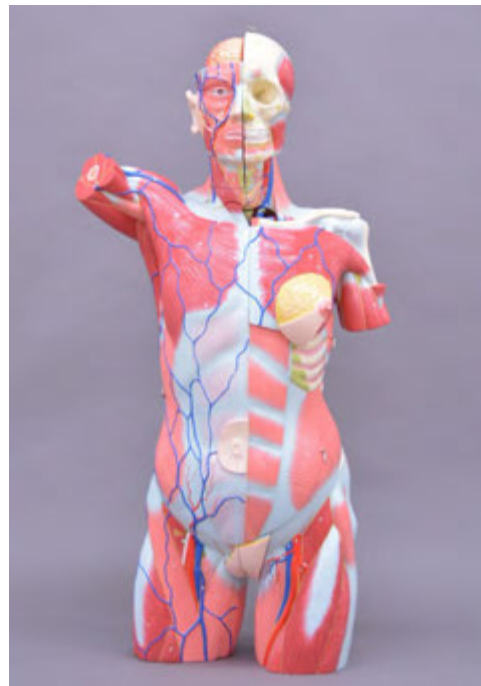13. Attach the right lower head using the pinholes, set the right brain to the head.

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
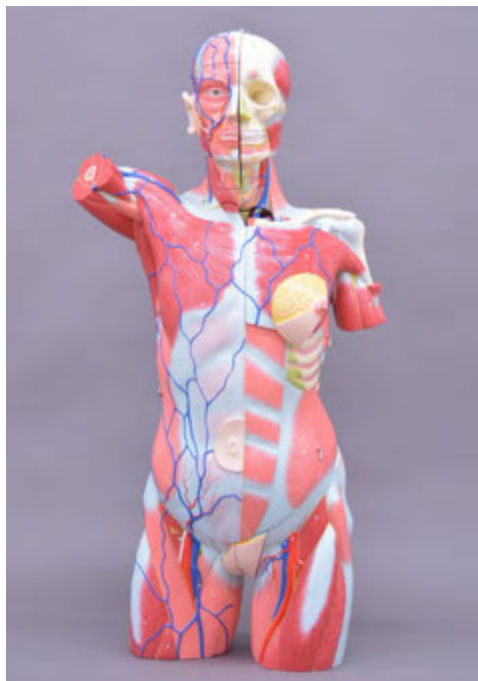
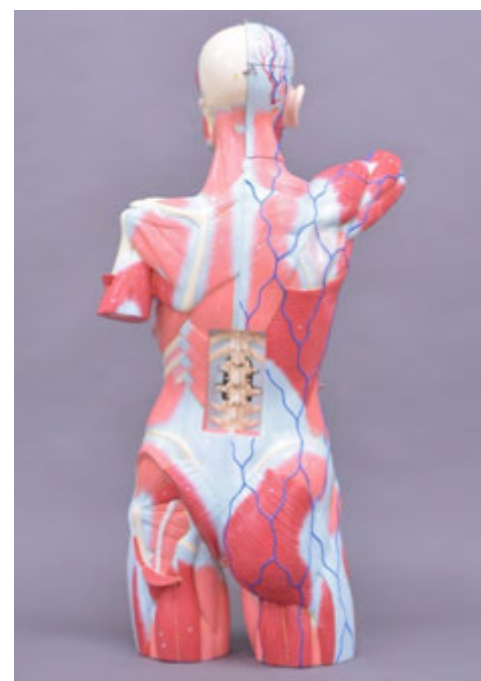
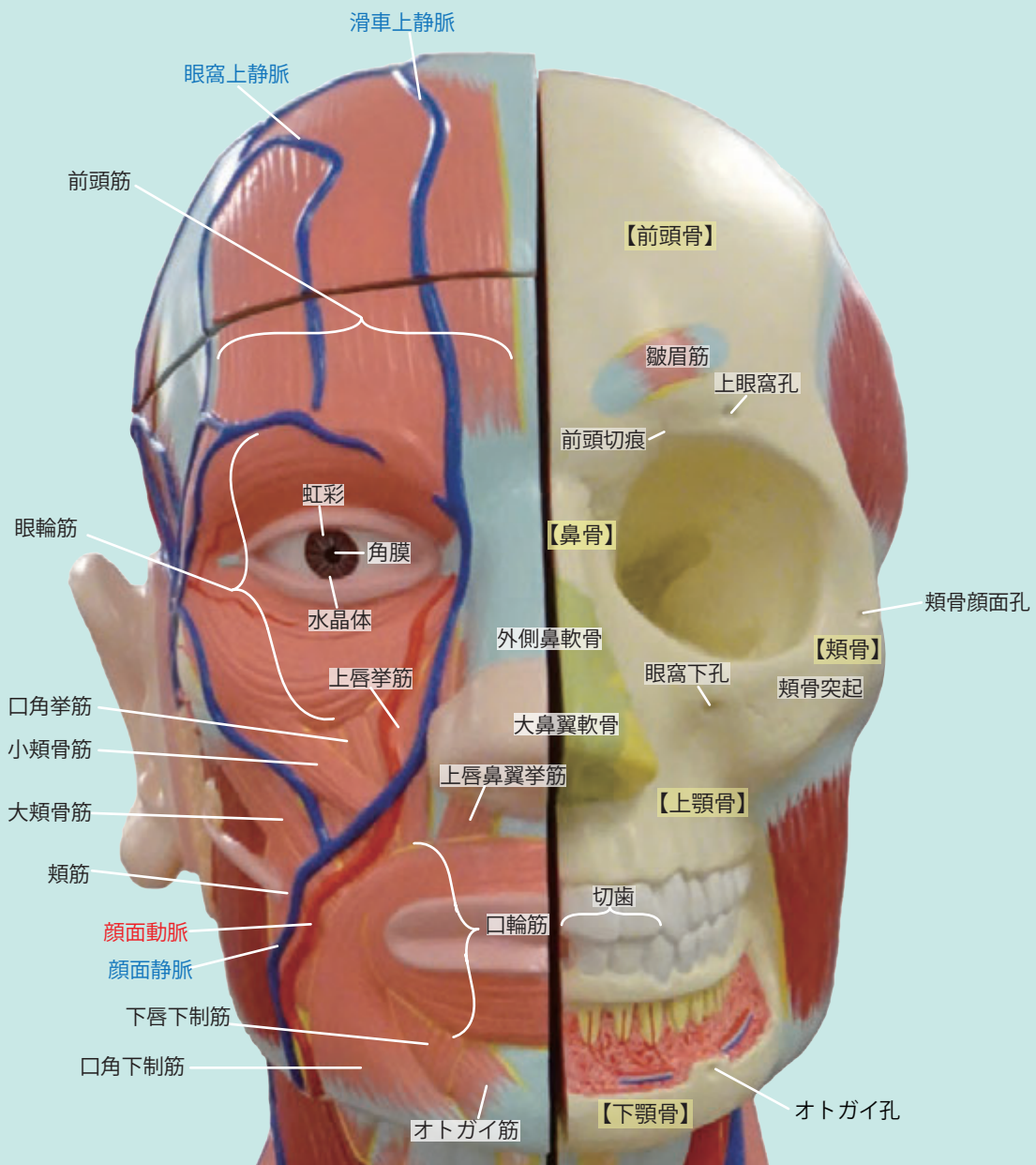
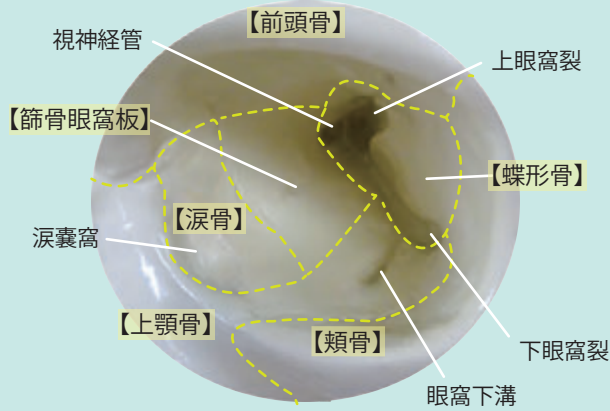
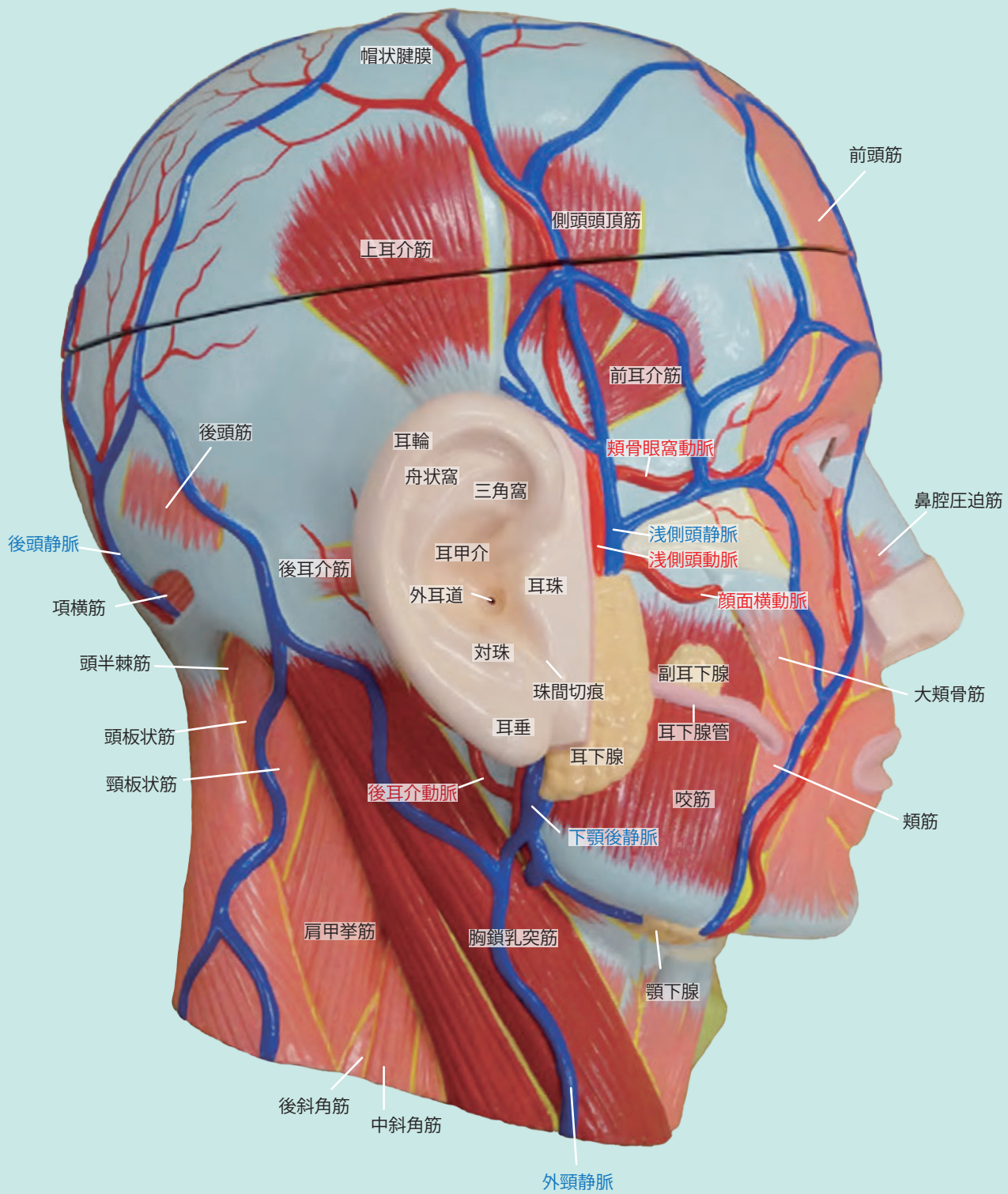
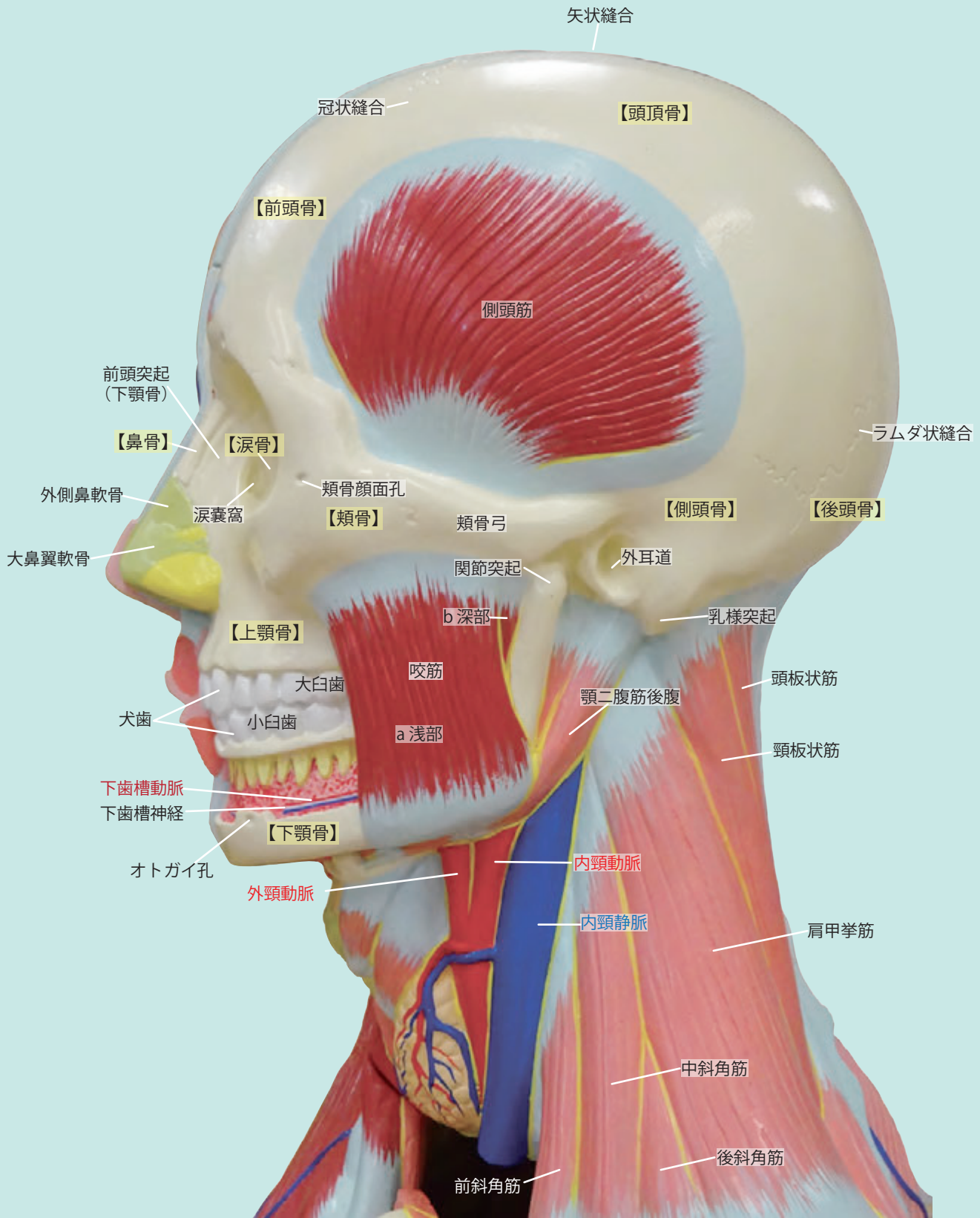
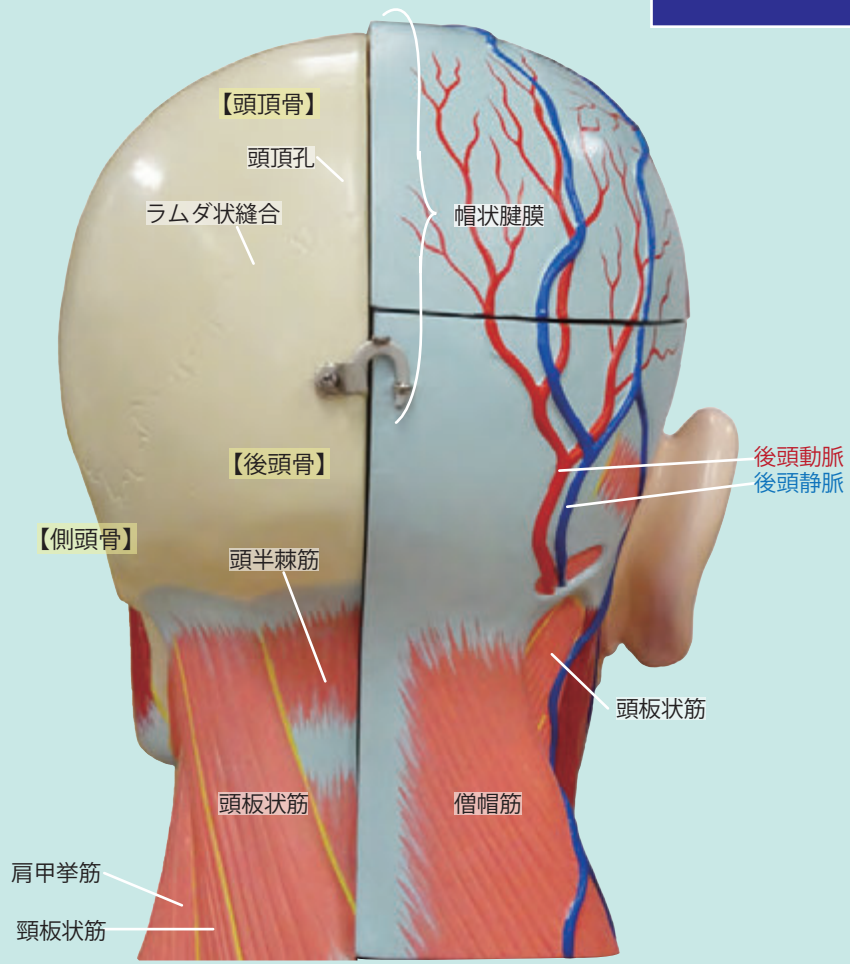
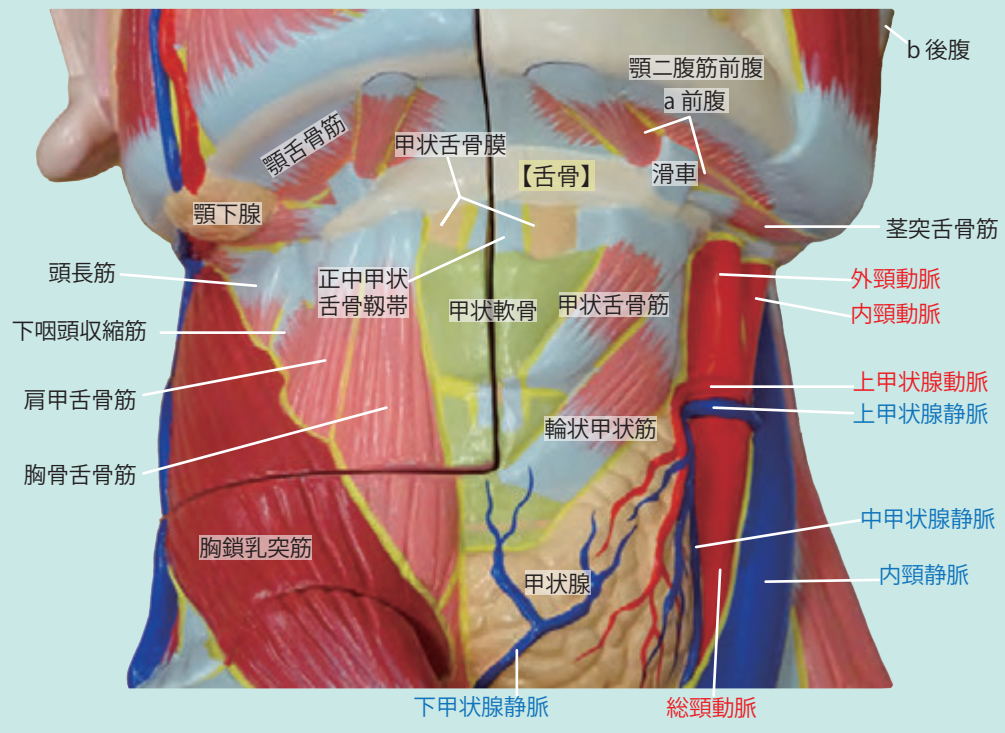
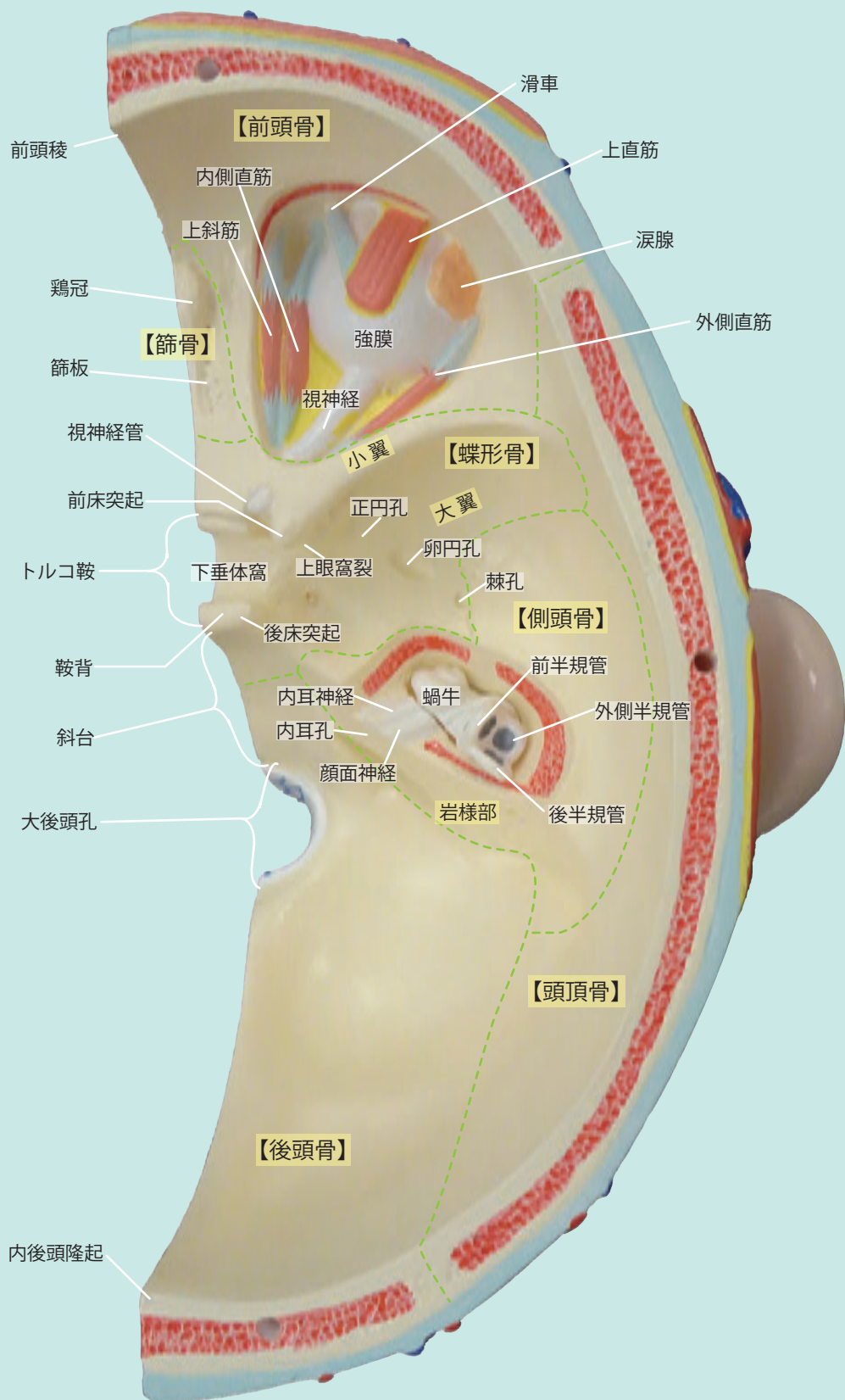
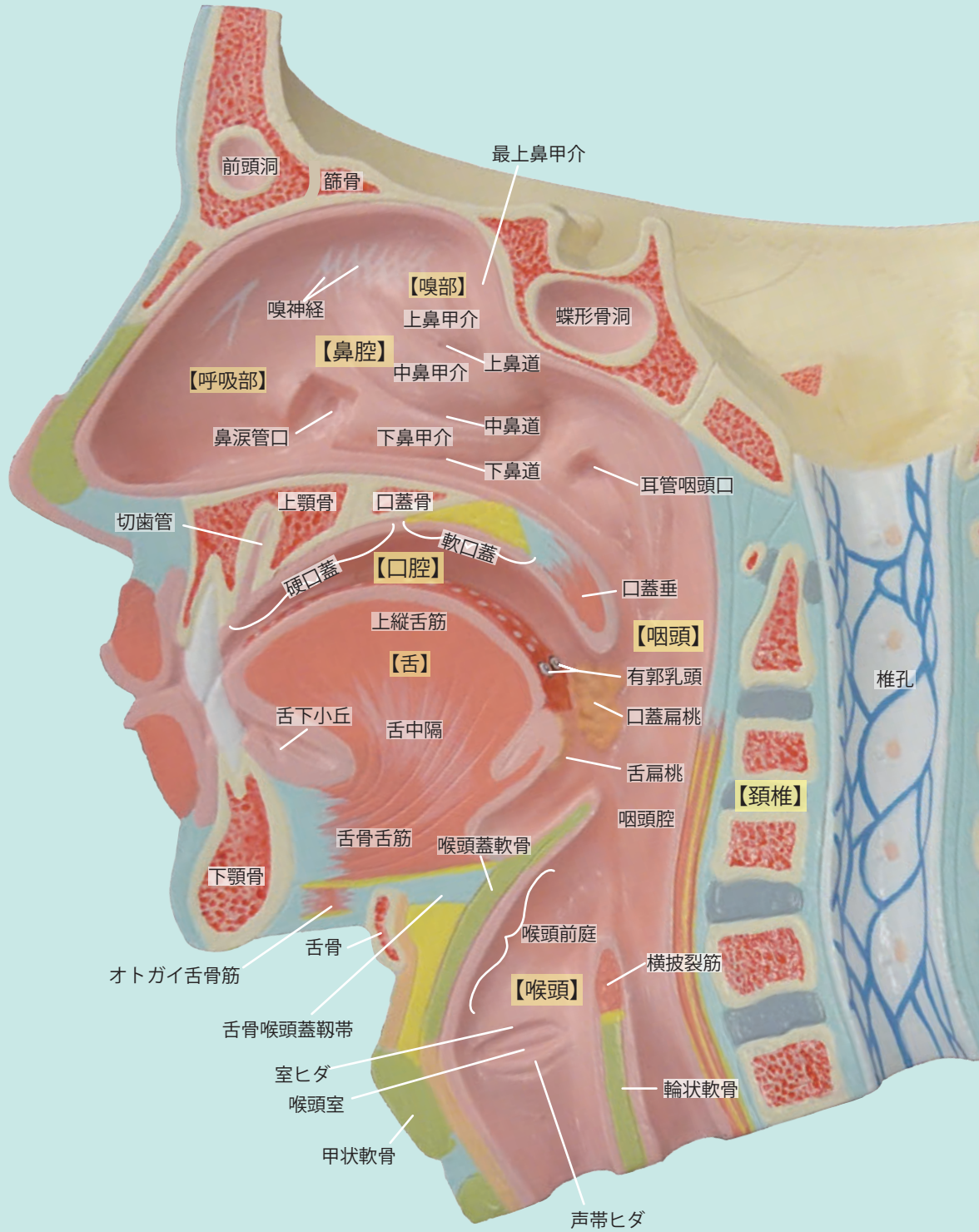
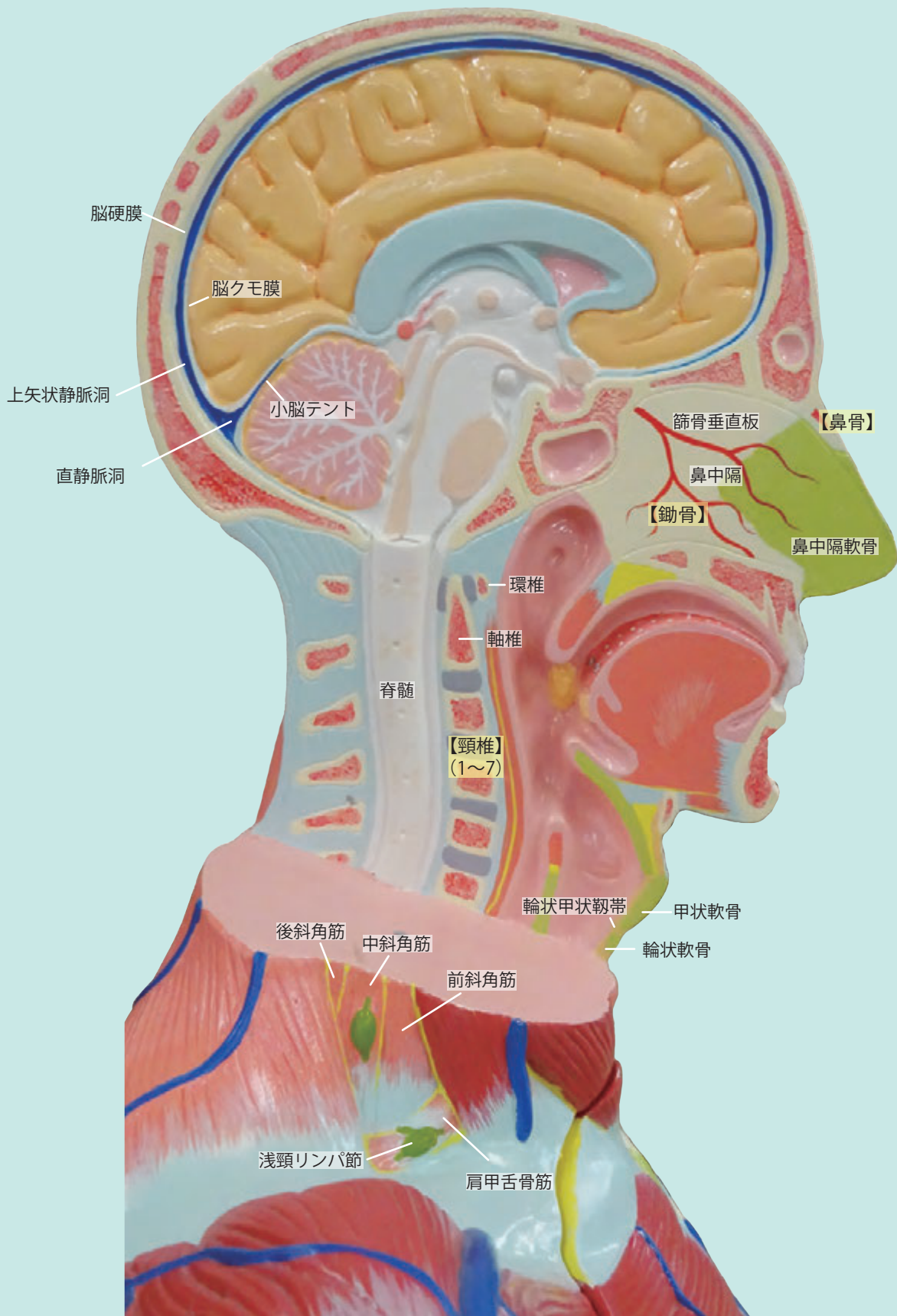
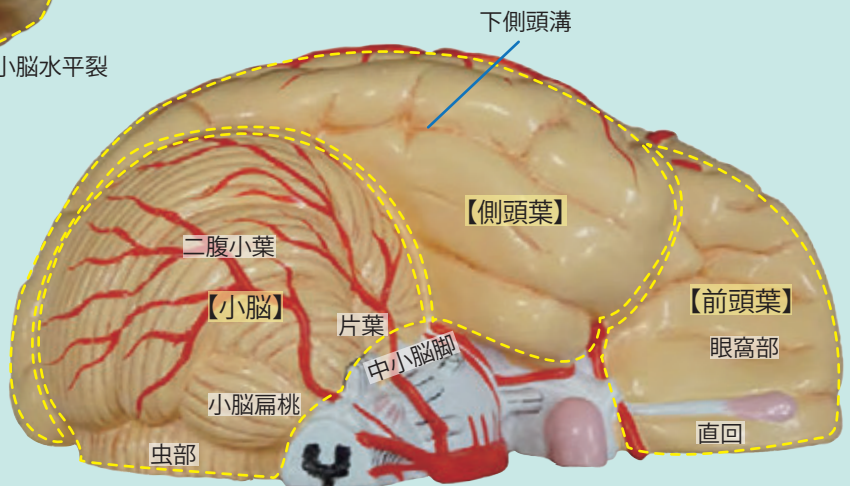
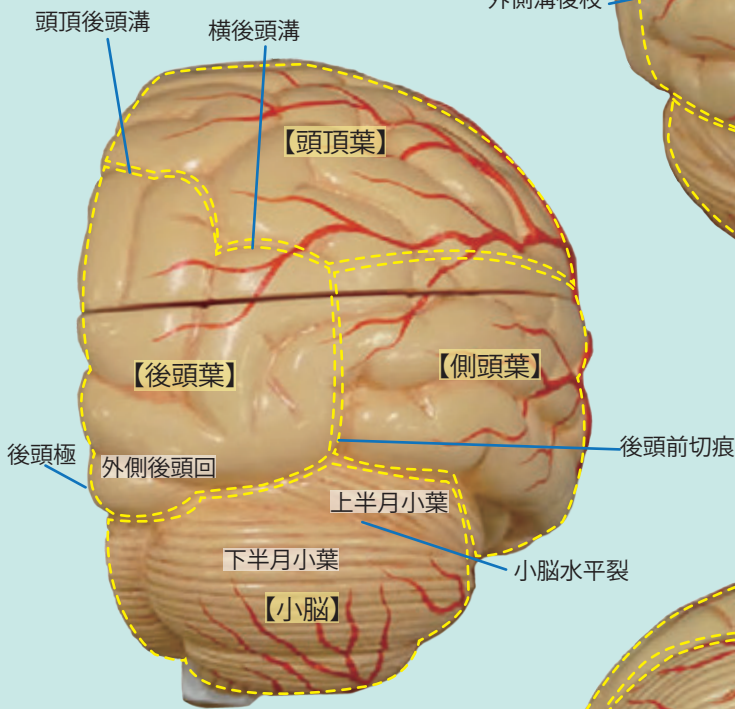
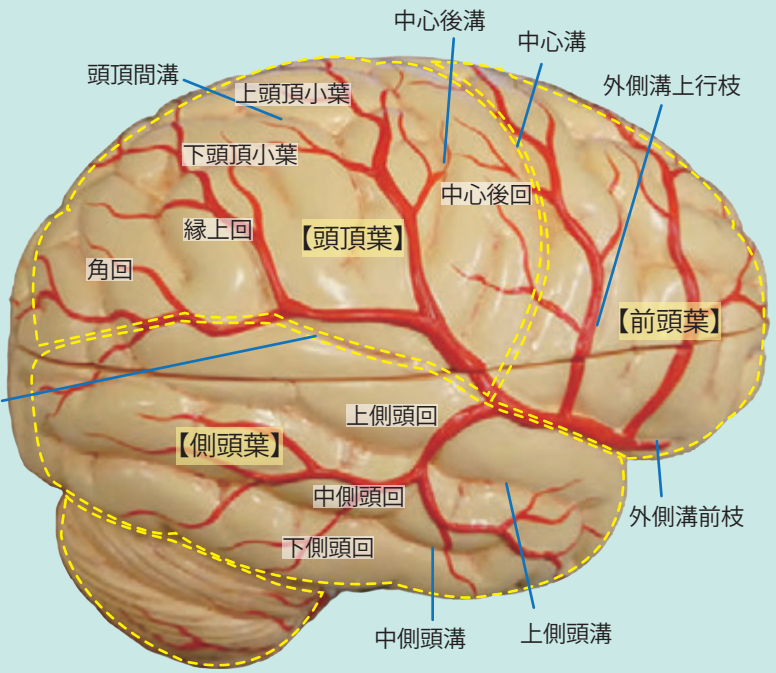
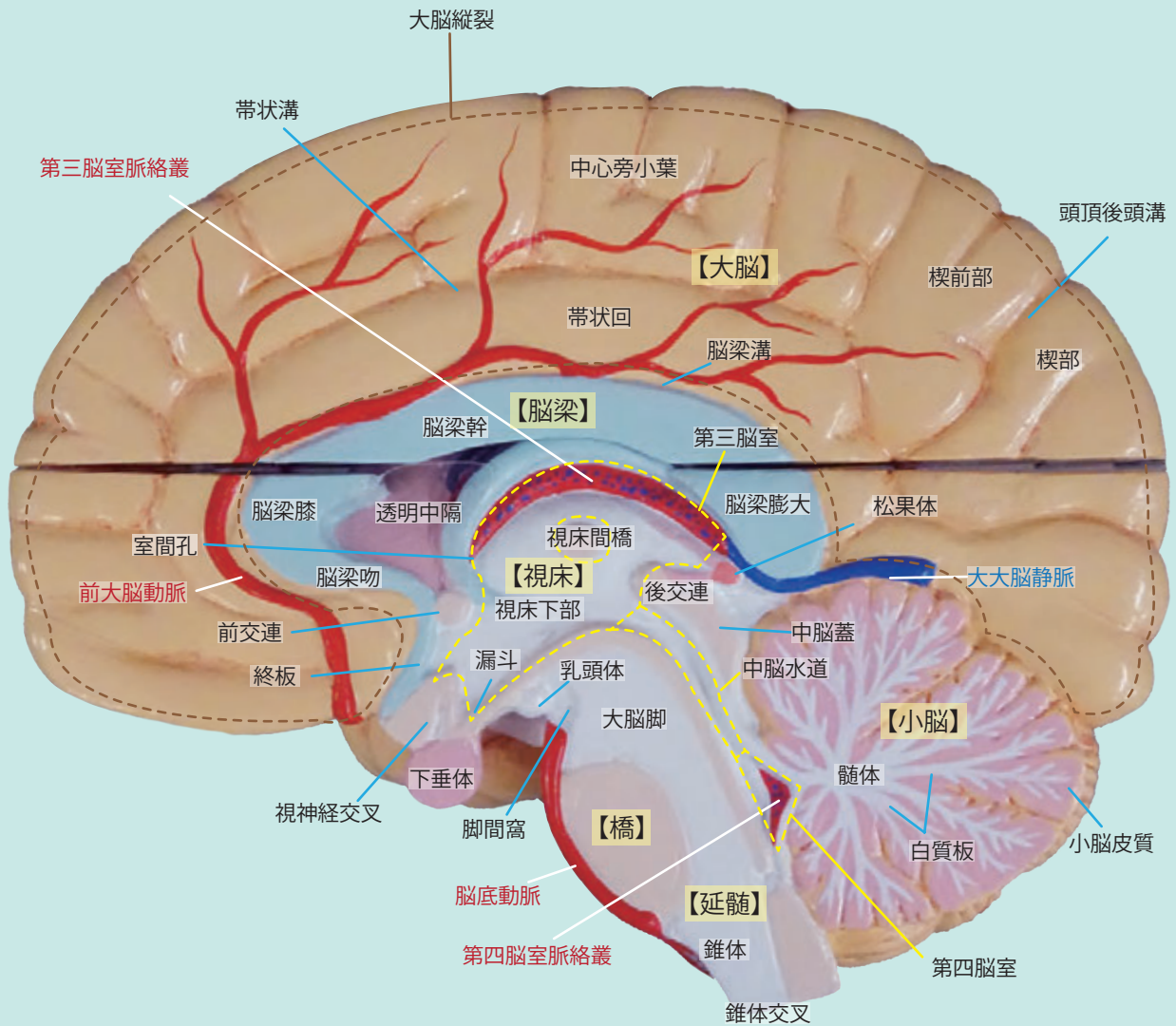
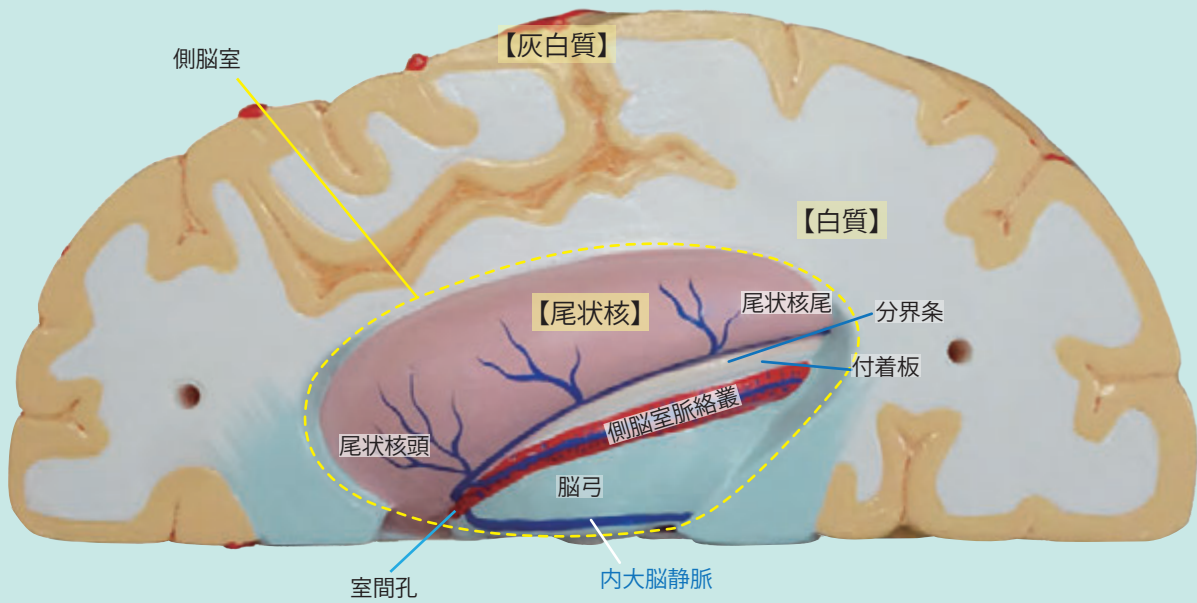
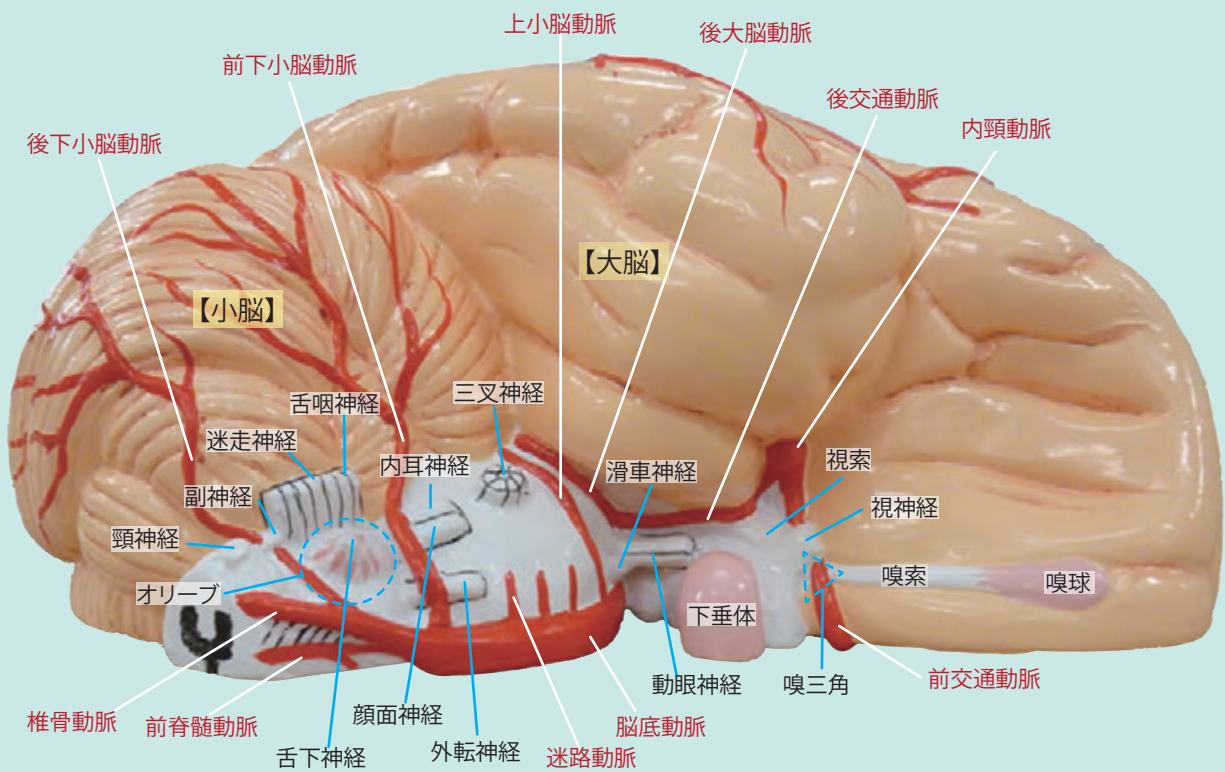
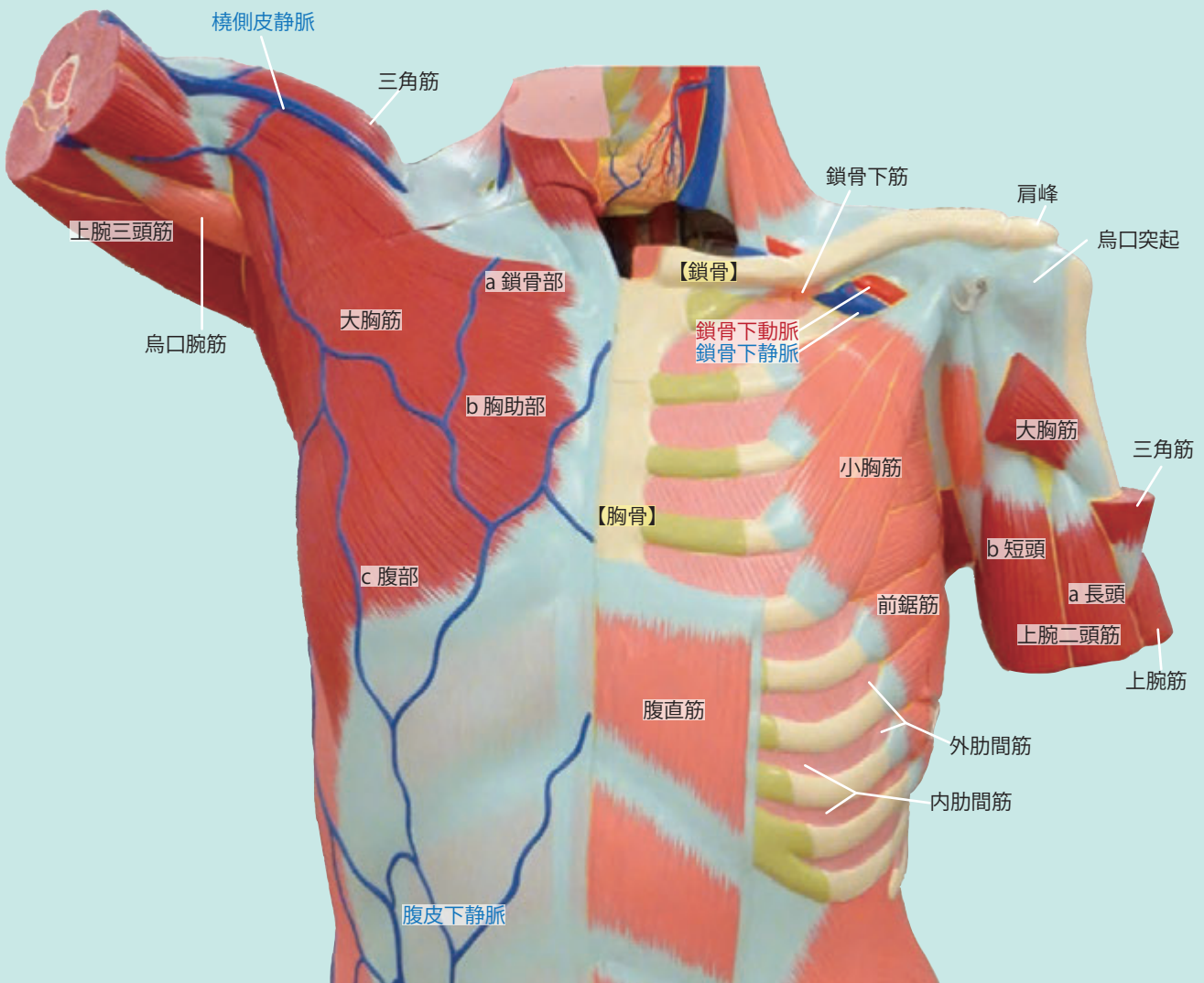
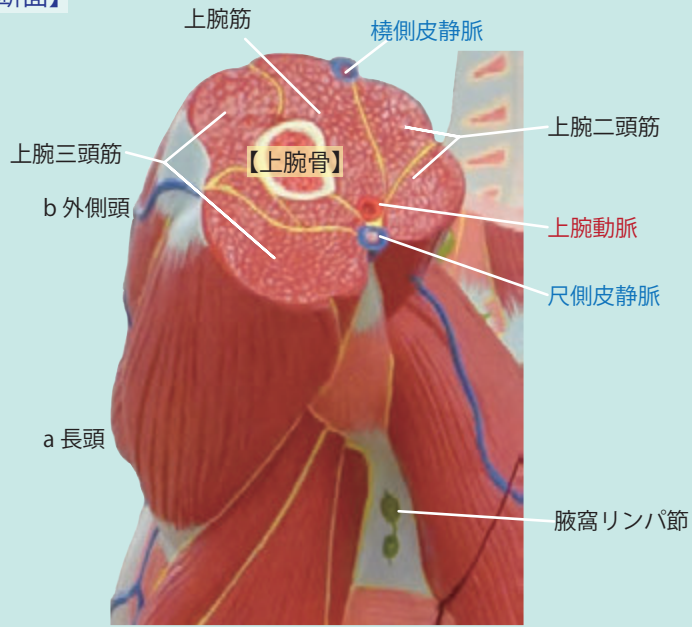
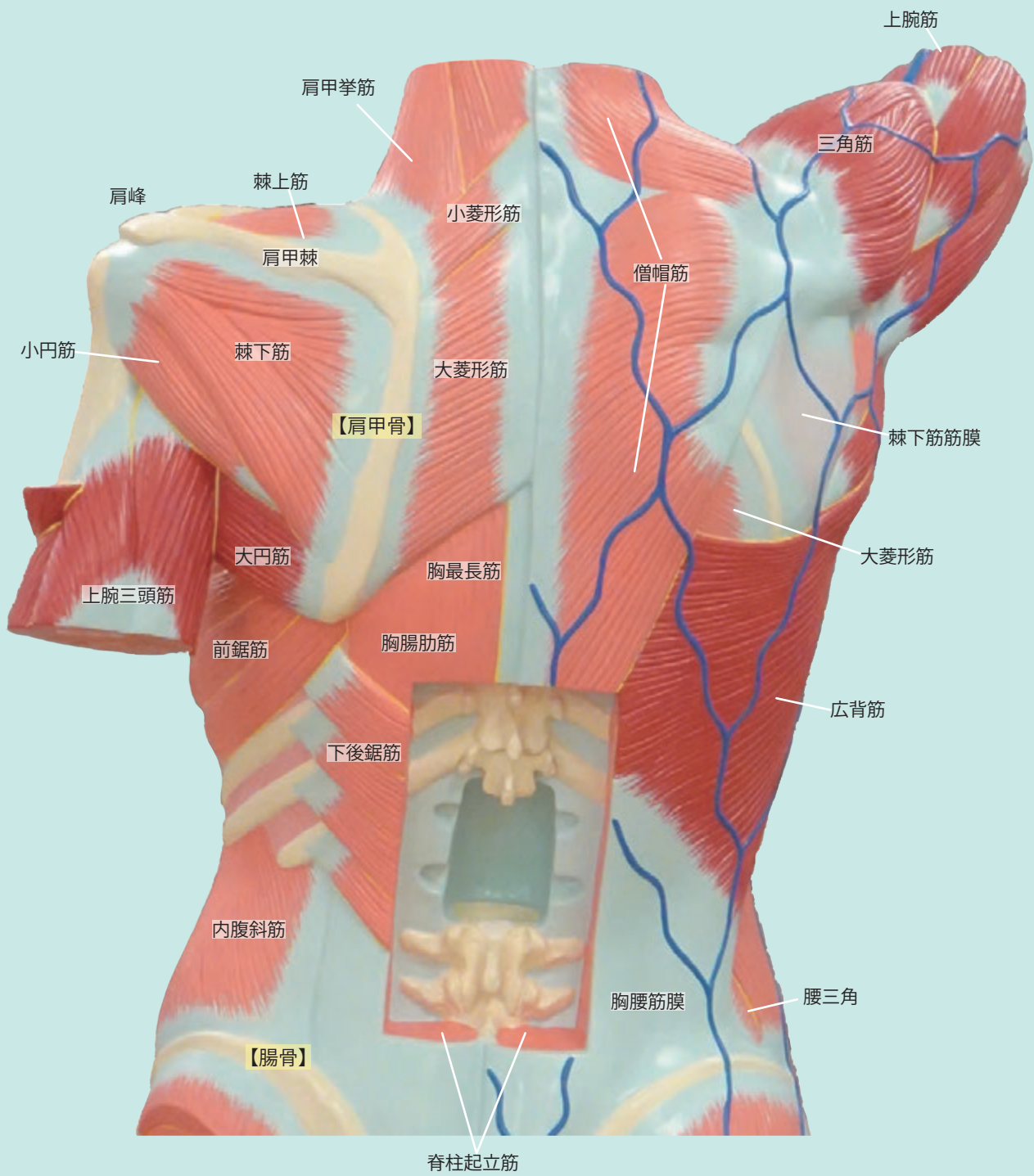
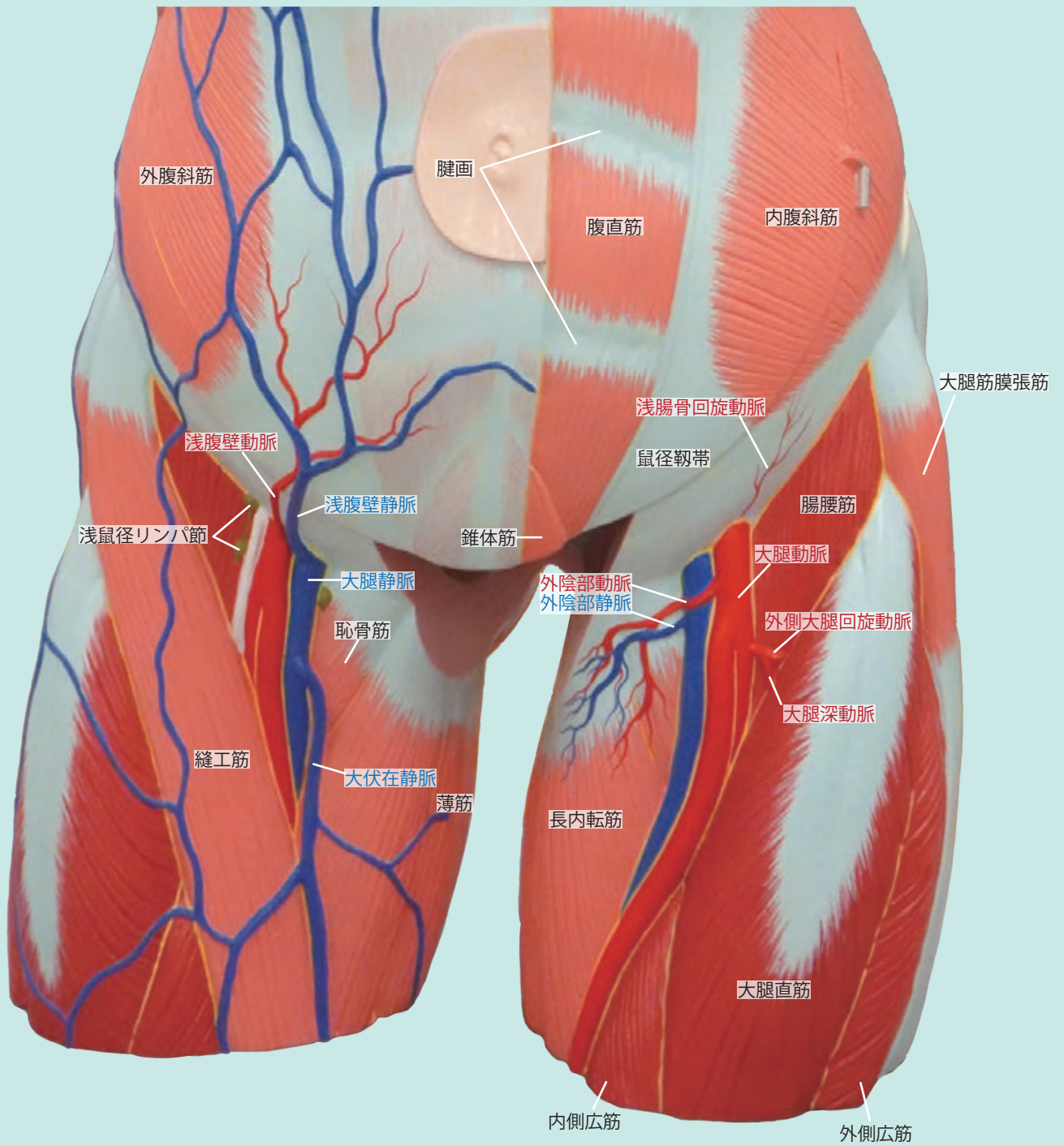
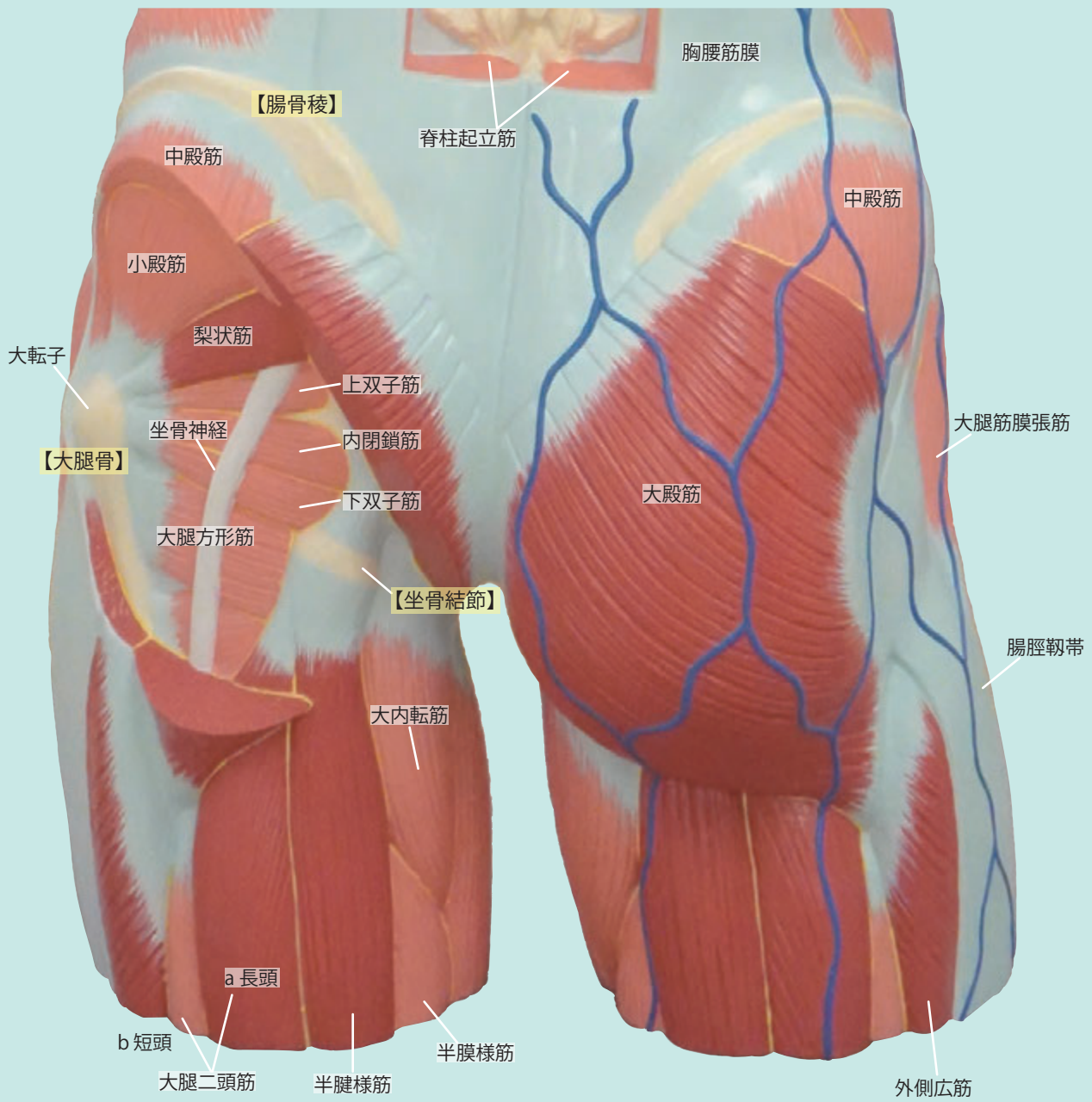
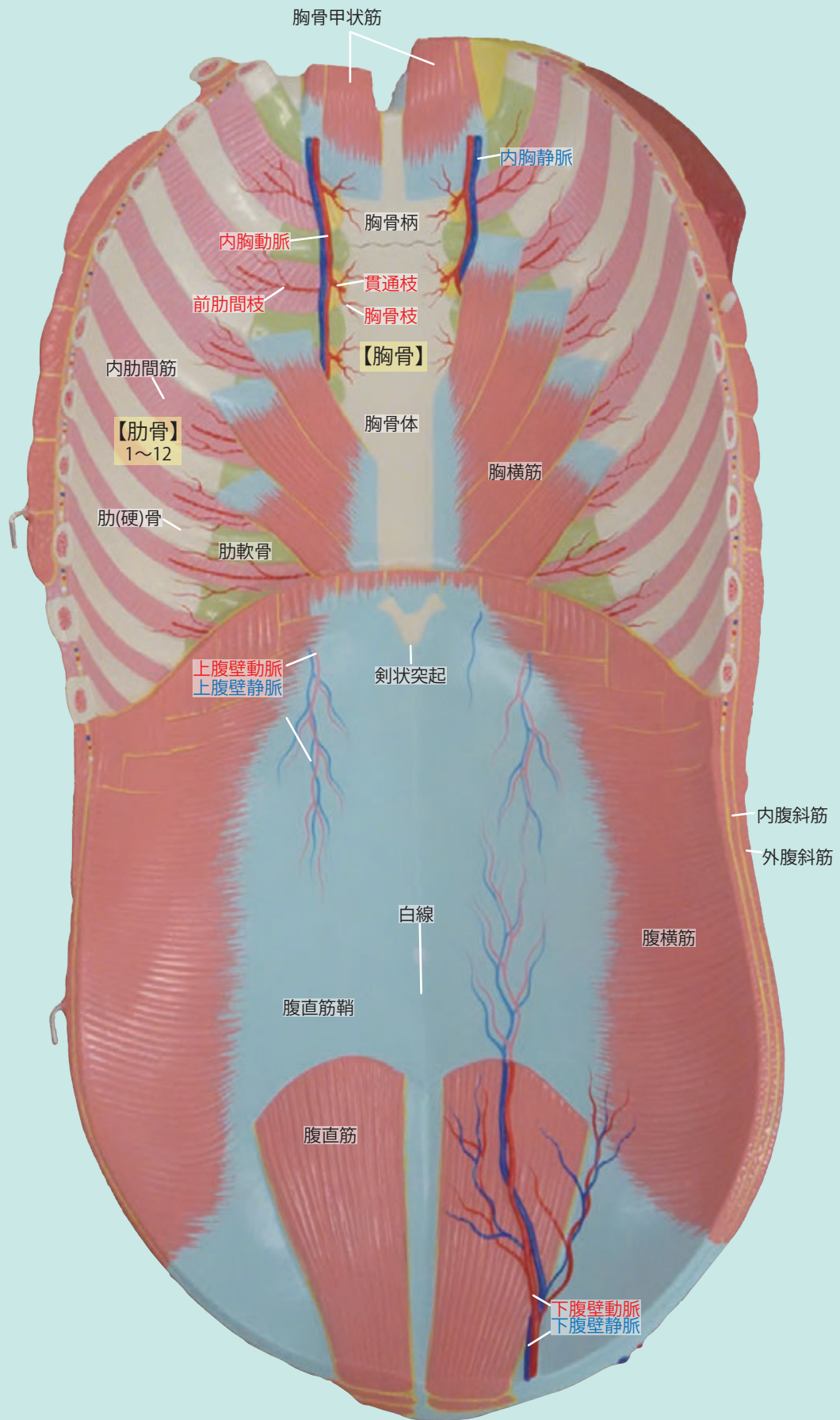
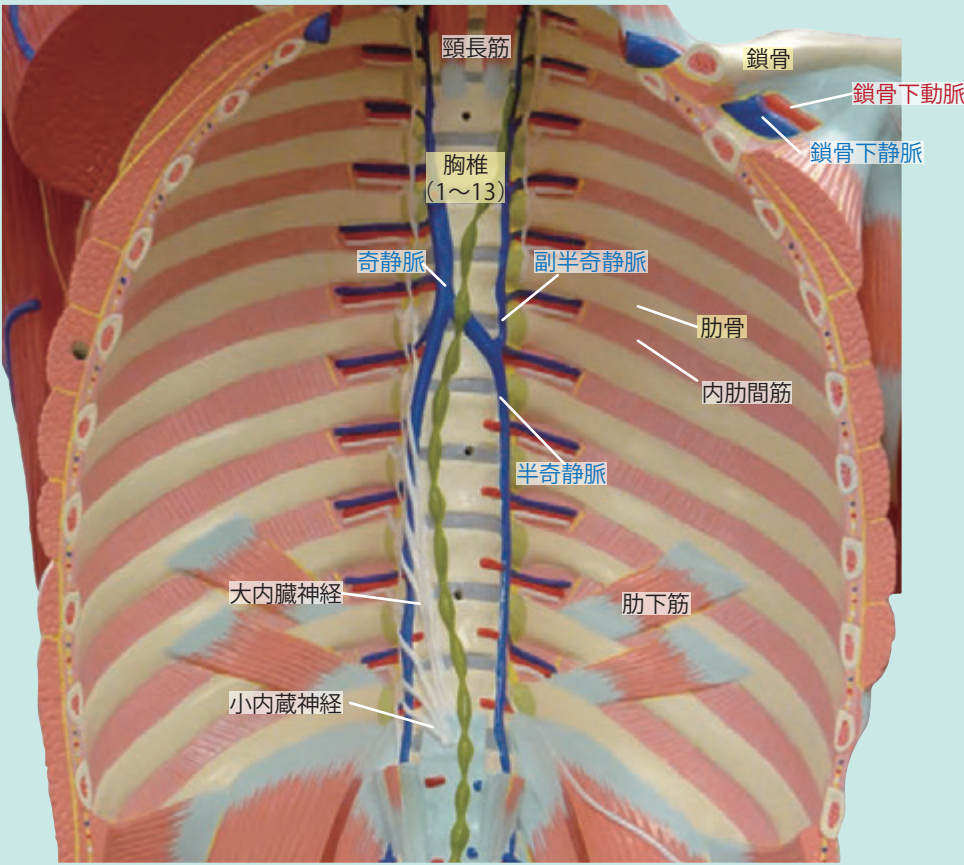
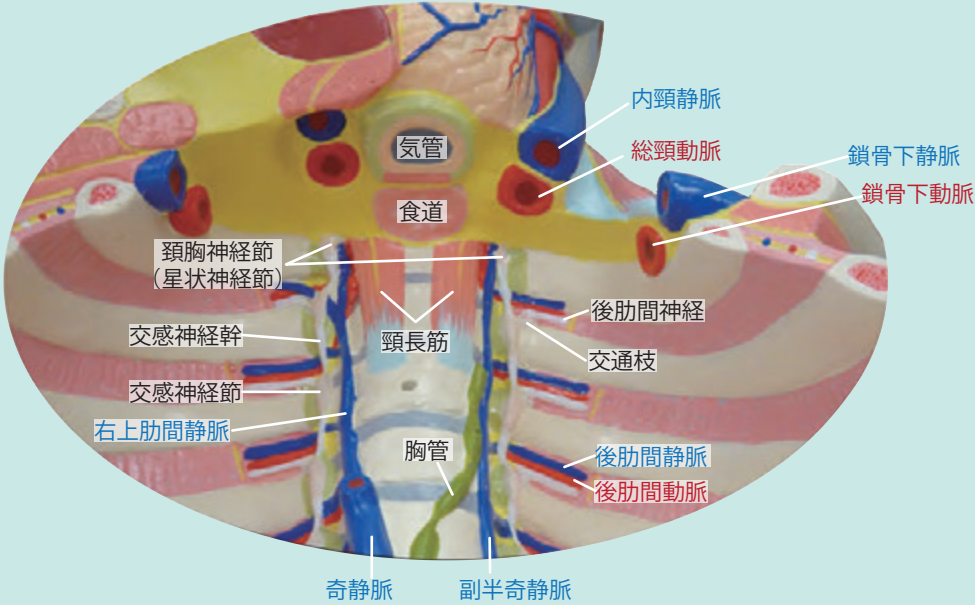
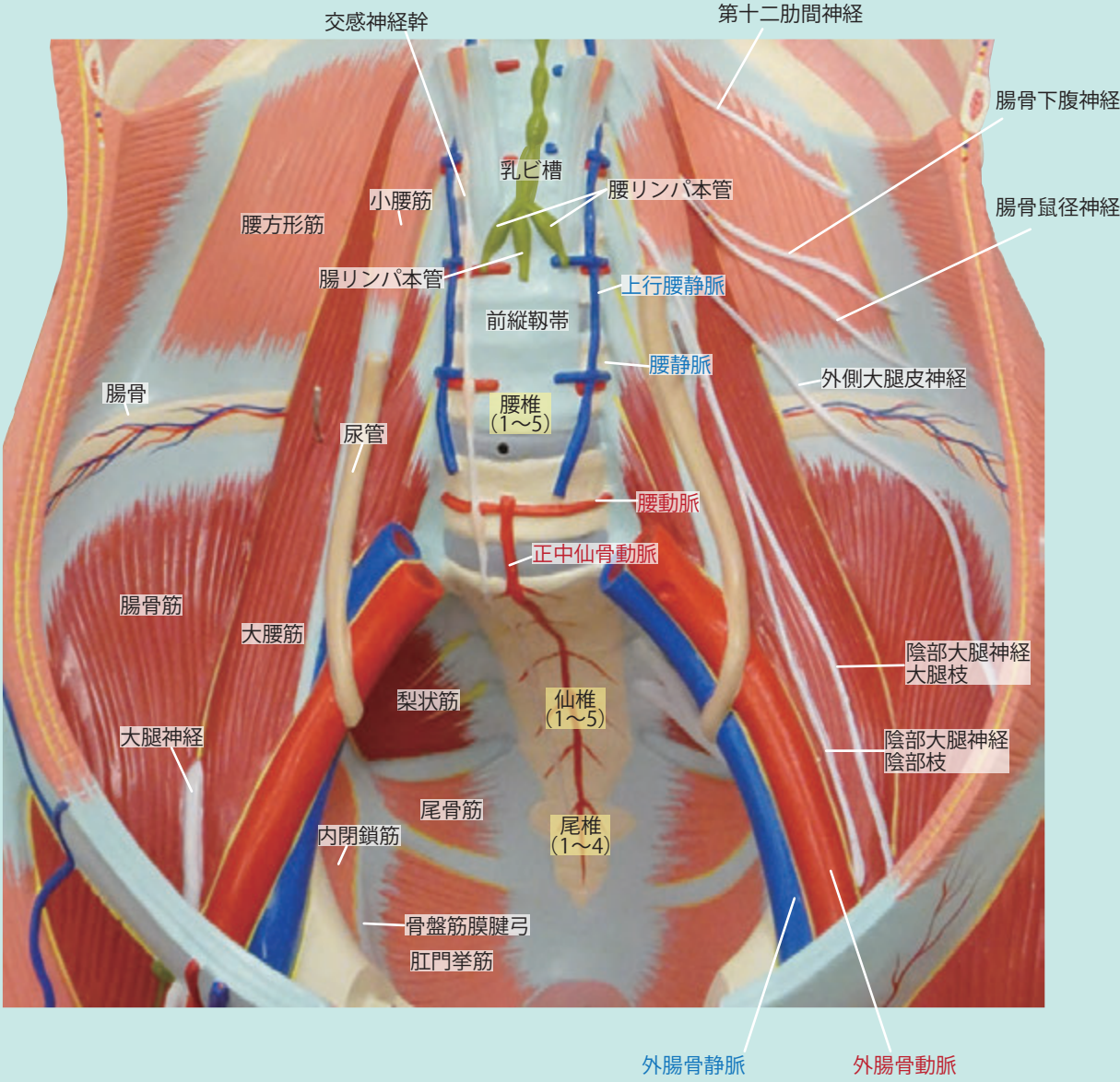
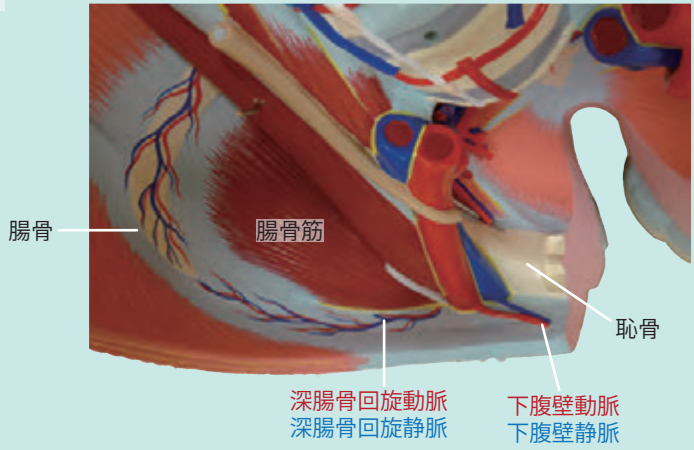
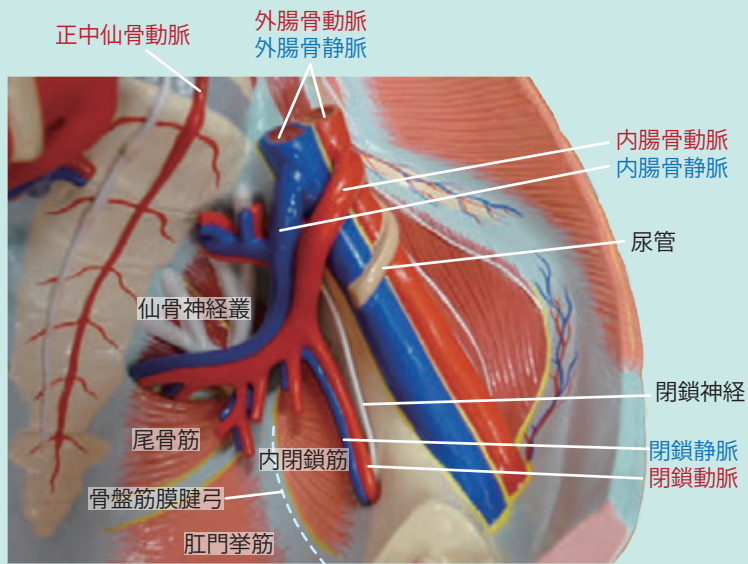
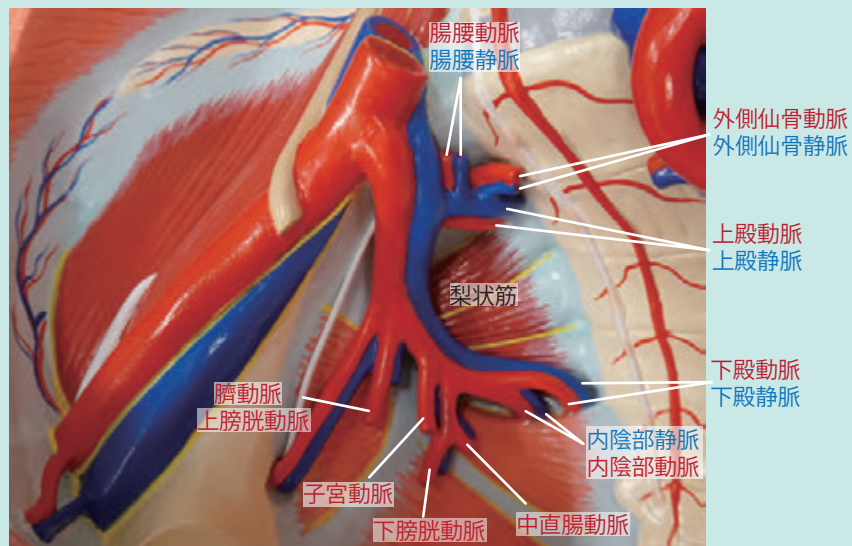
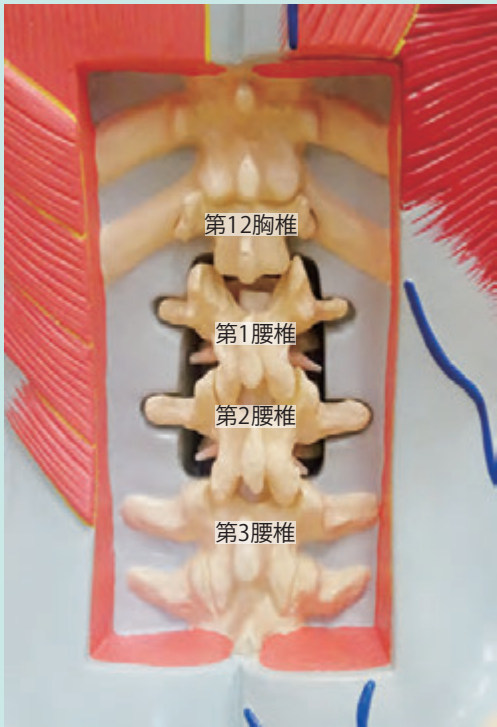
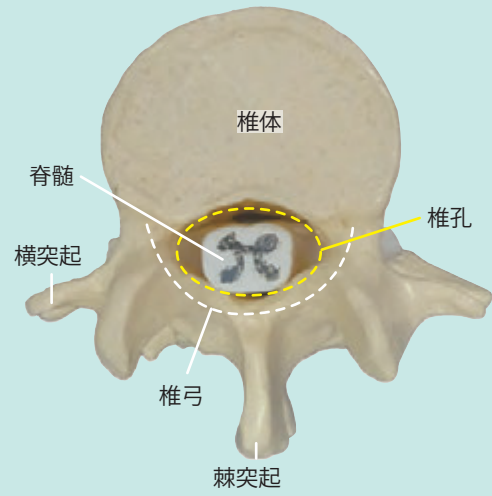
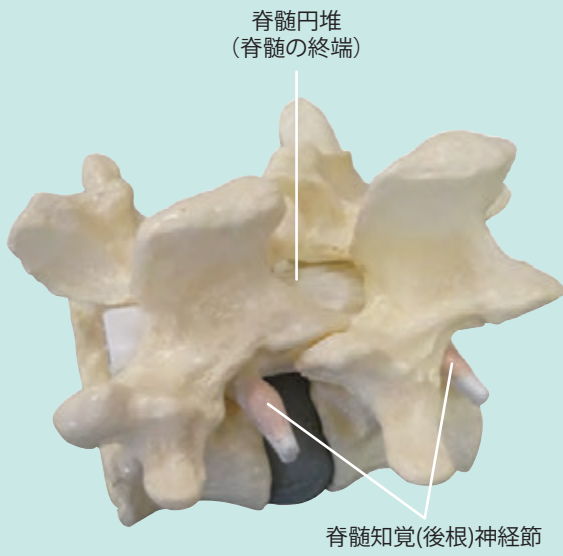
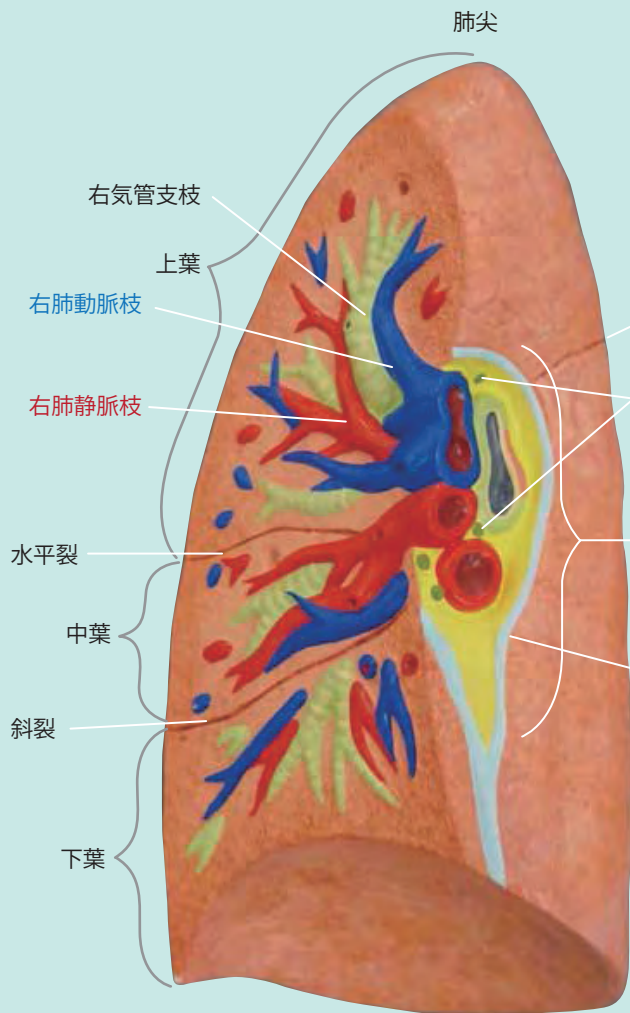
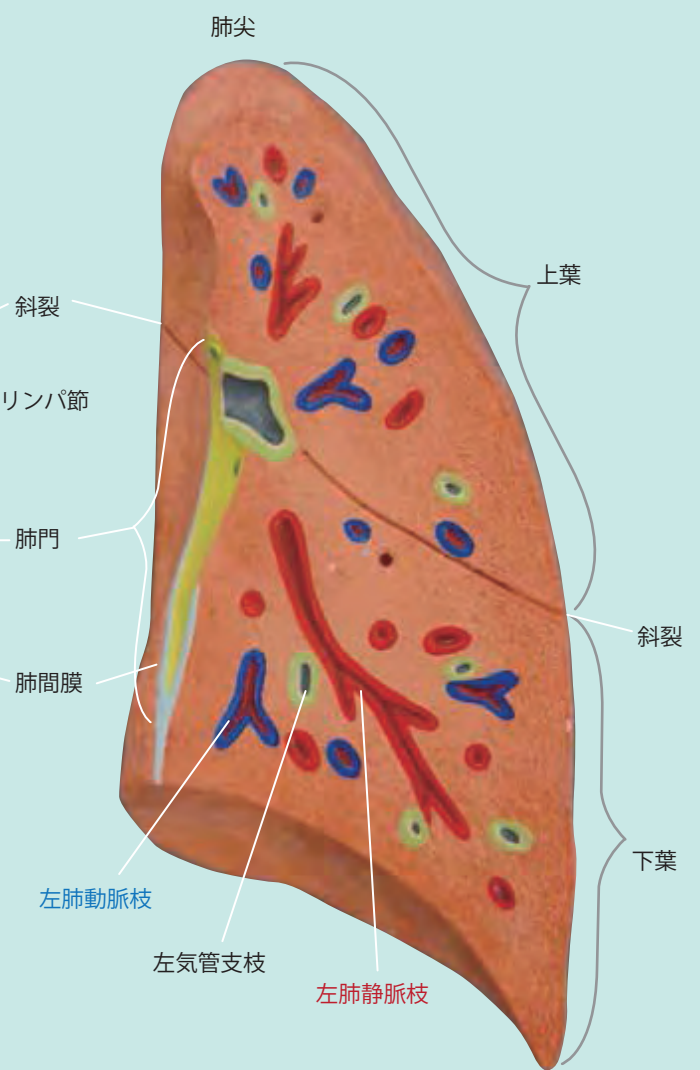
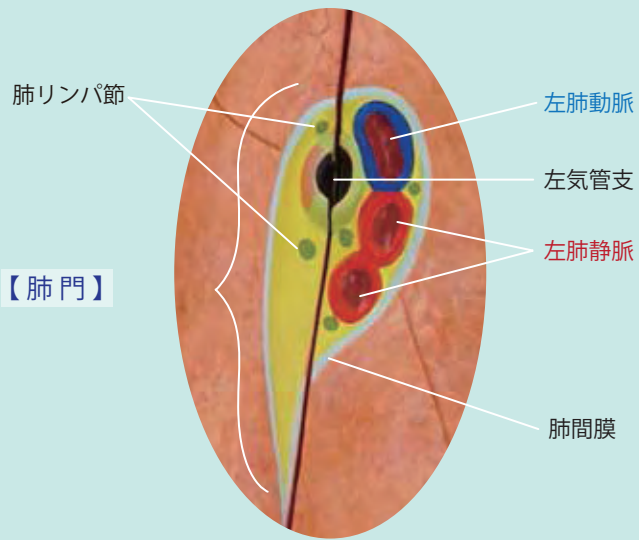
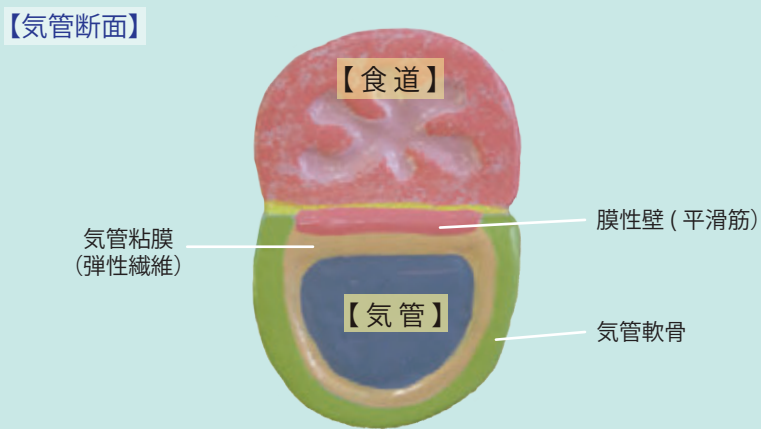
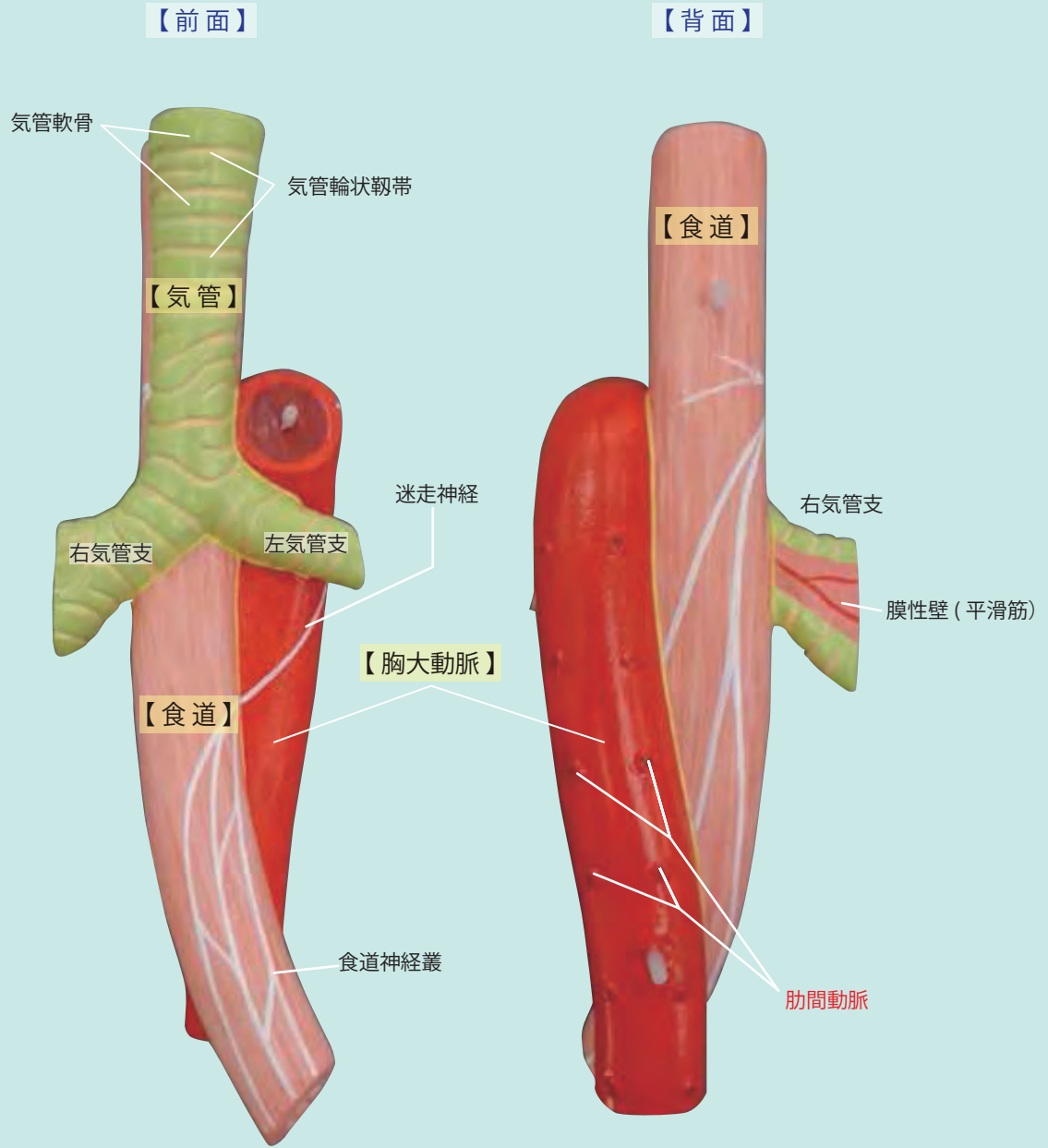
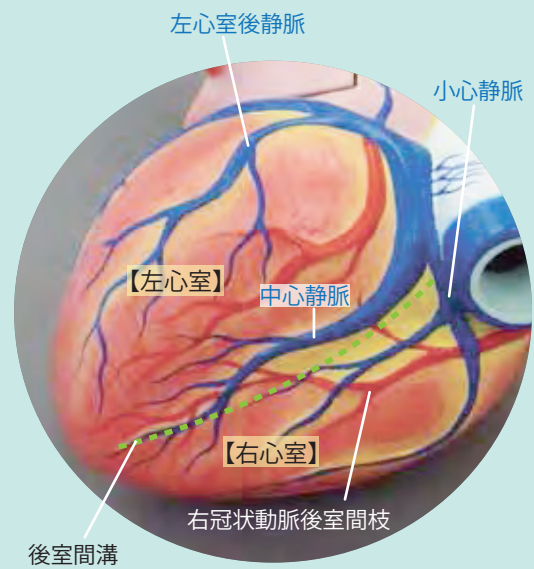
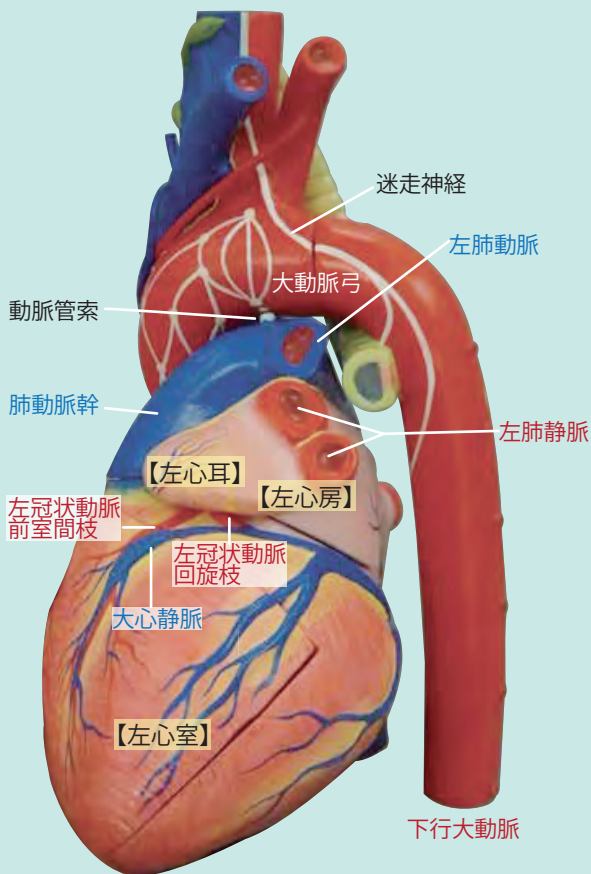
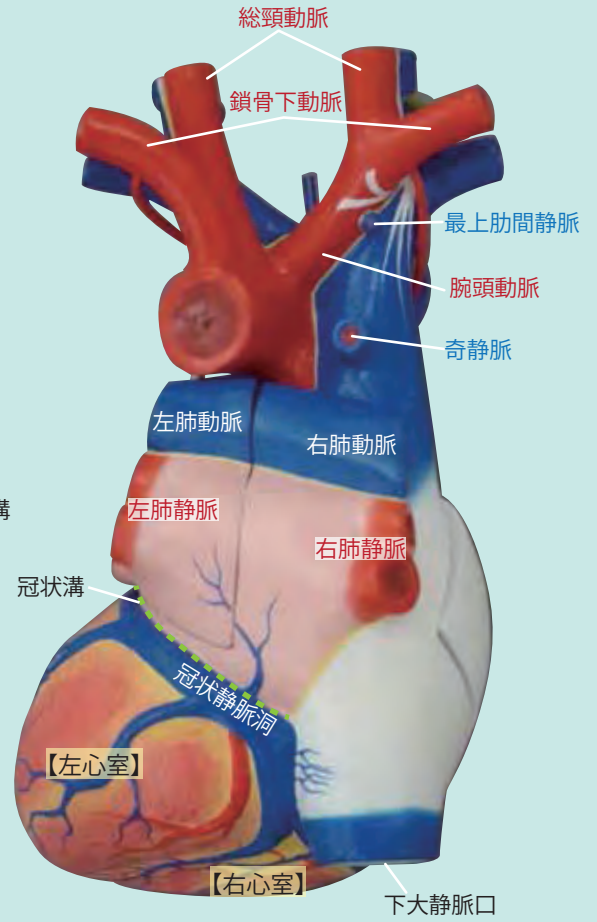
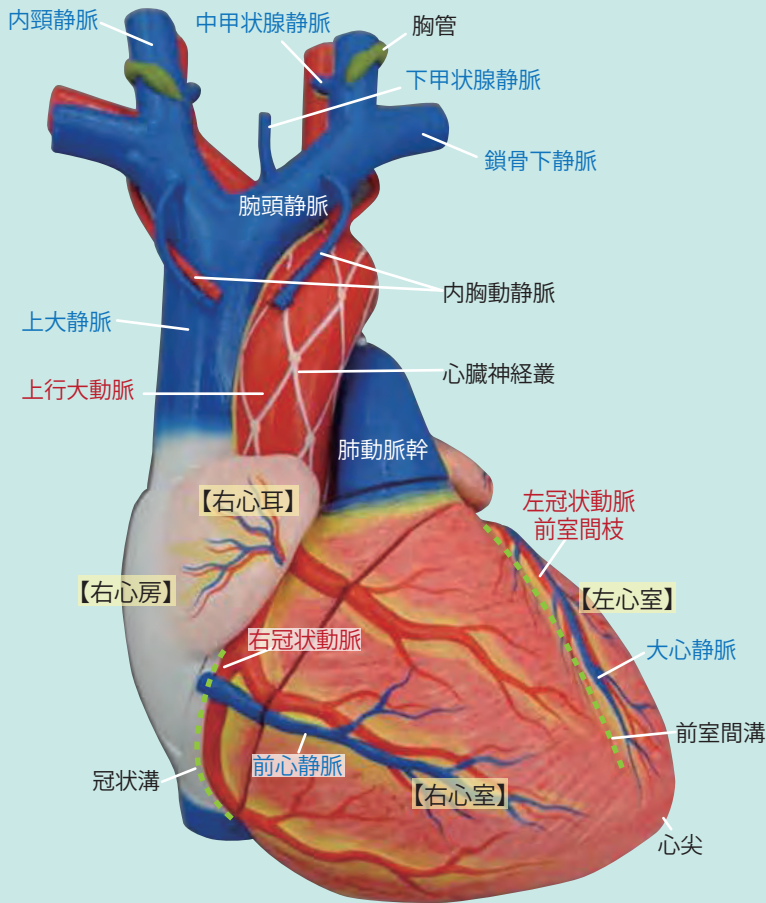
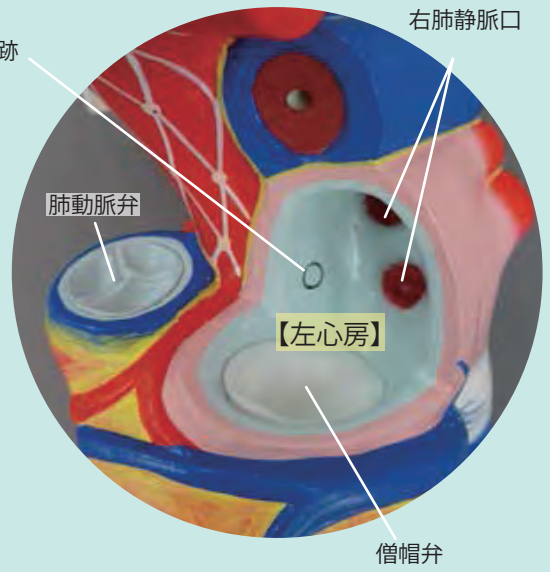
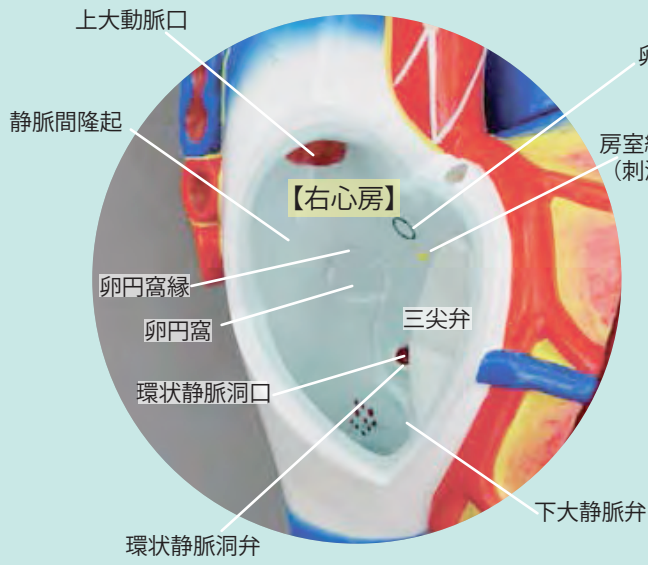
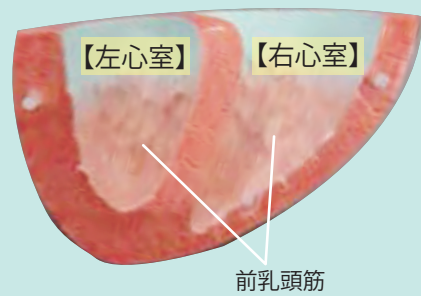
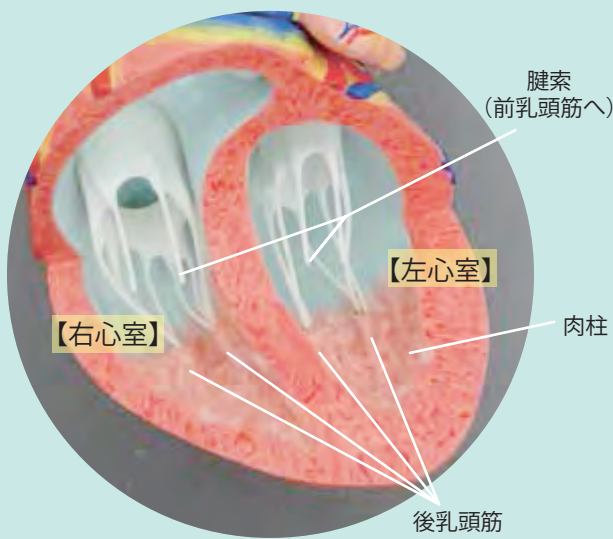
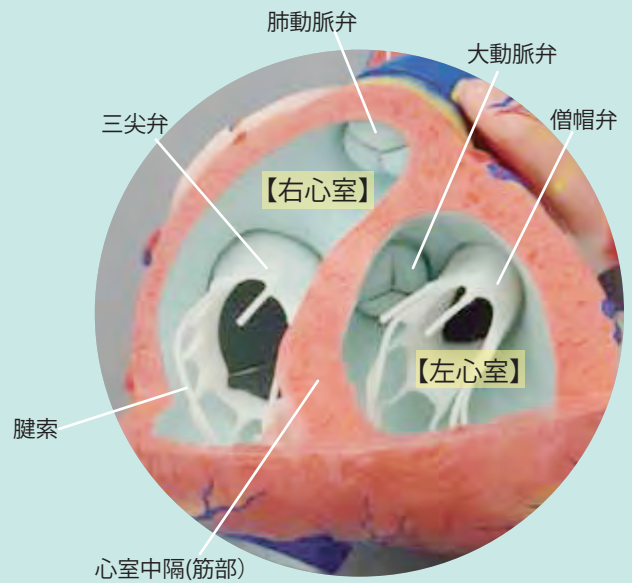
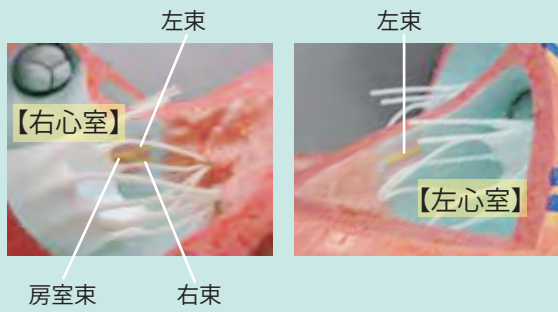
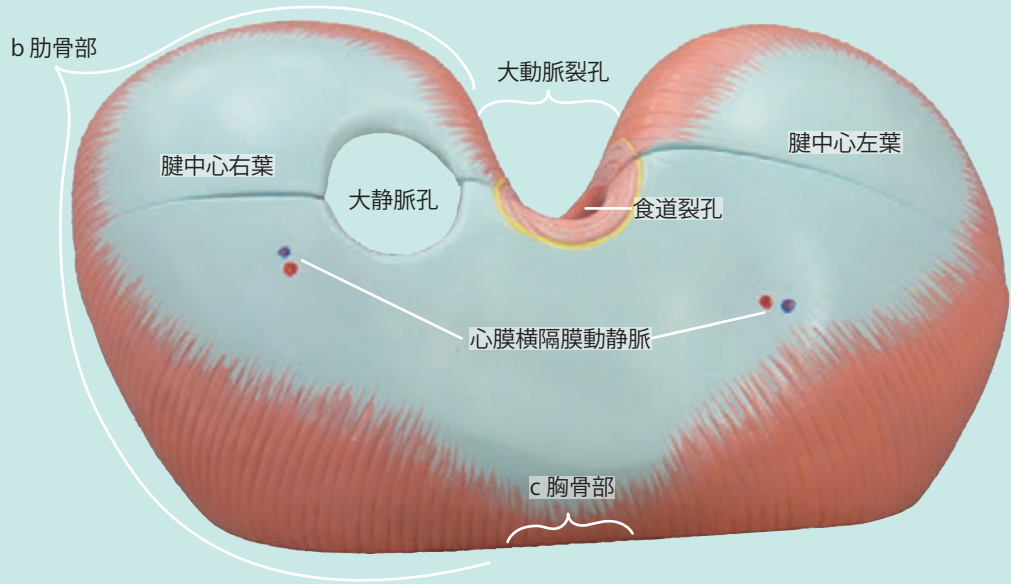
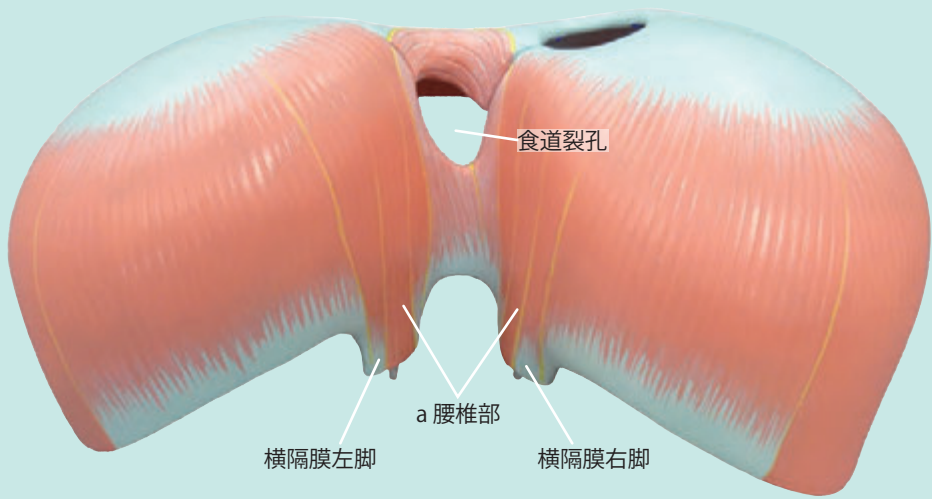
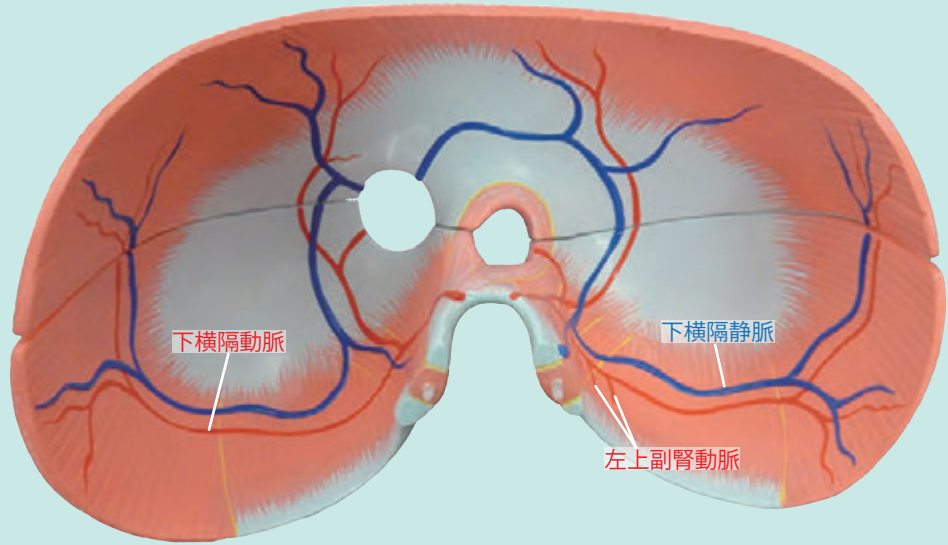
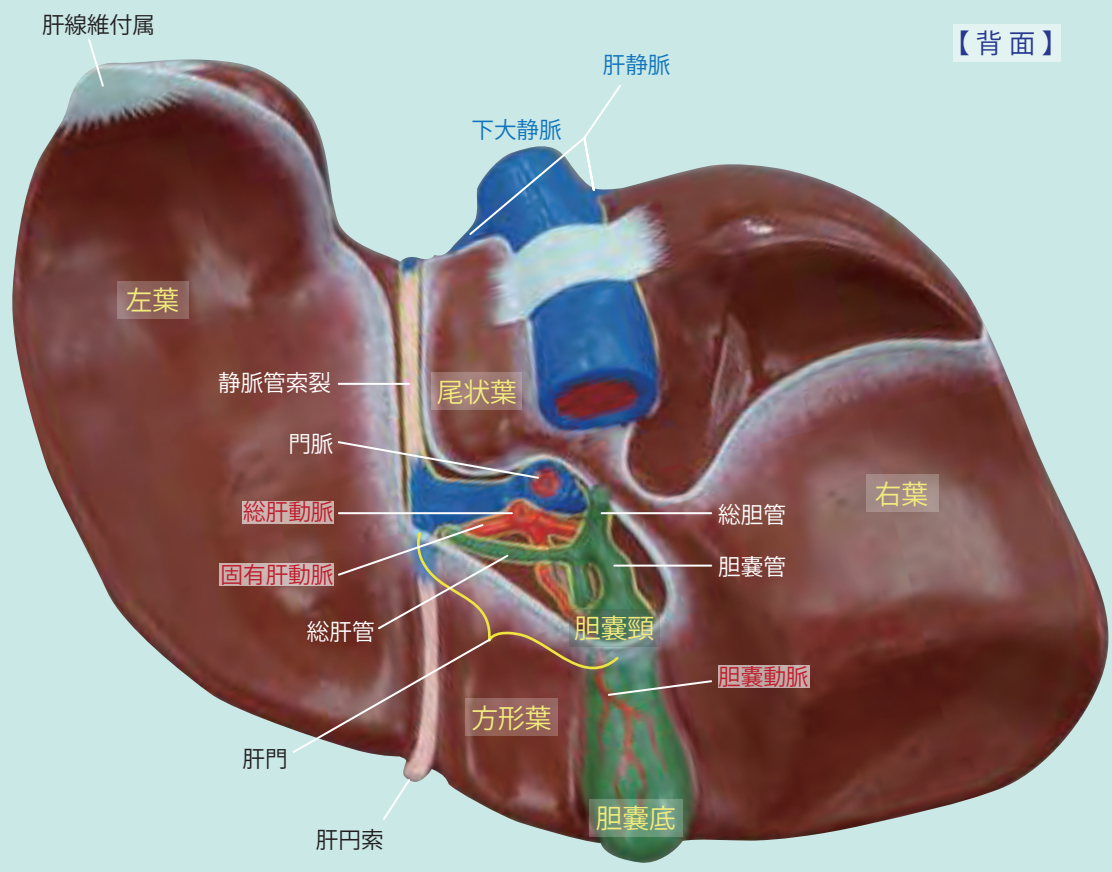
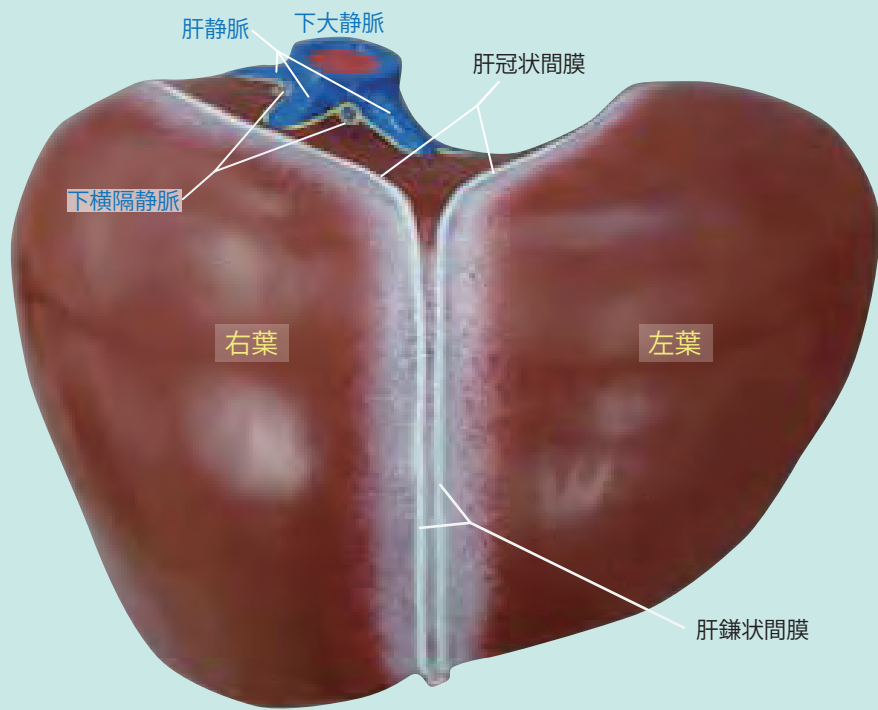
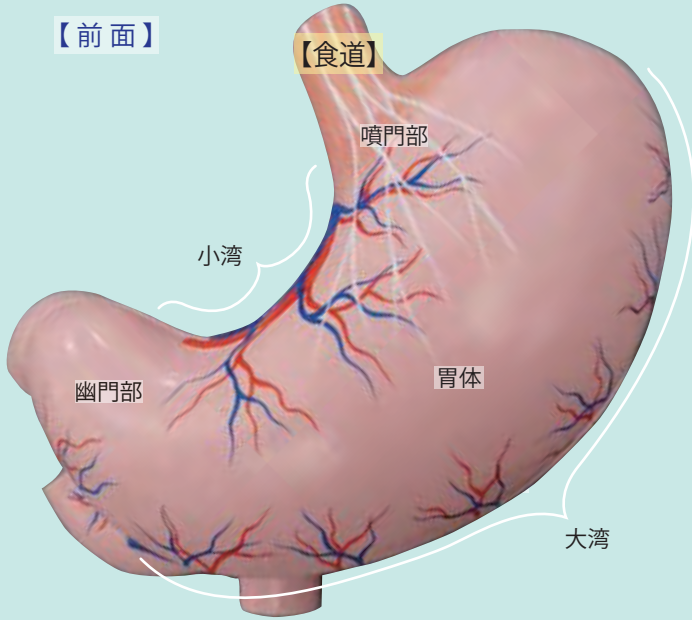
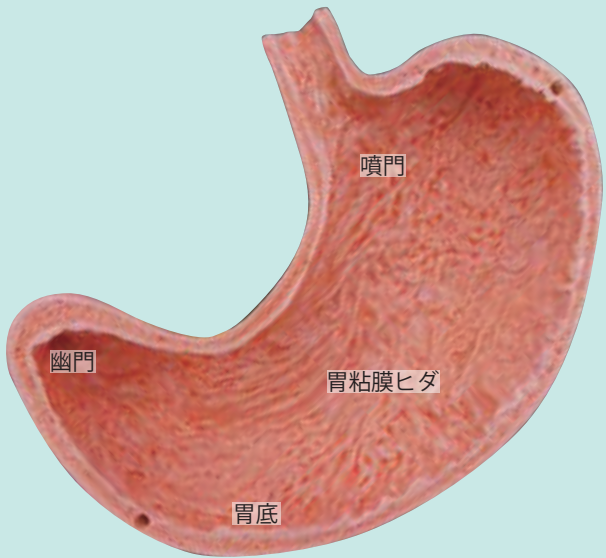
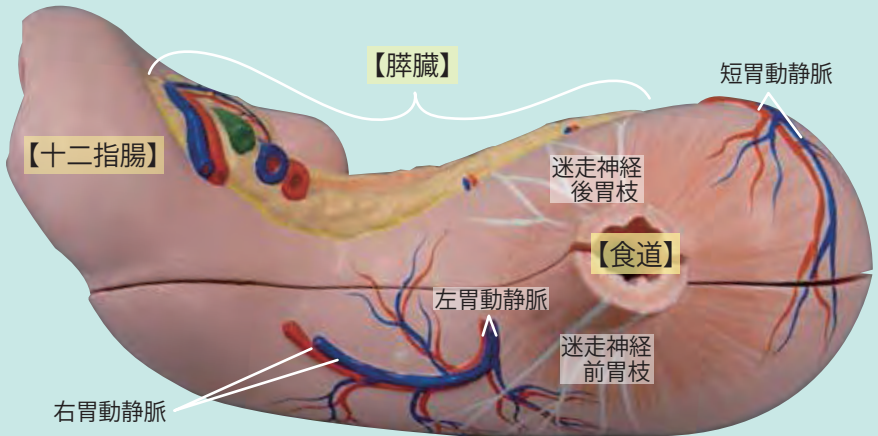
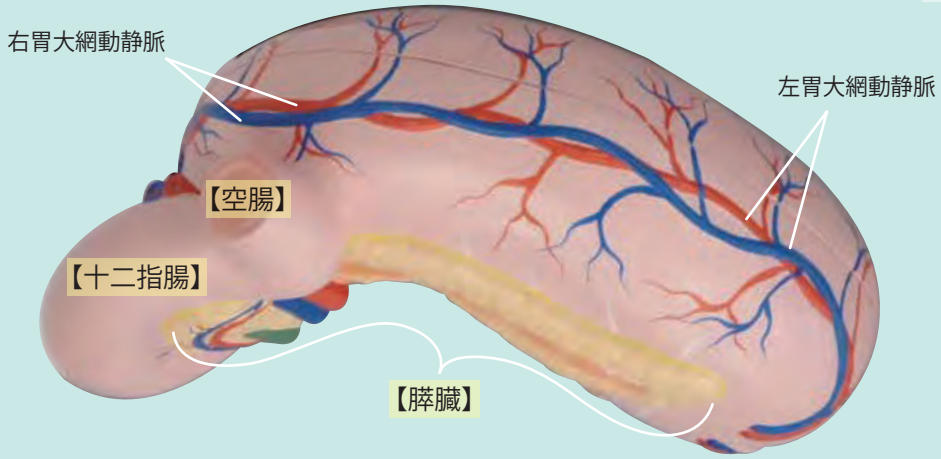
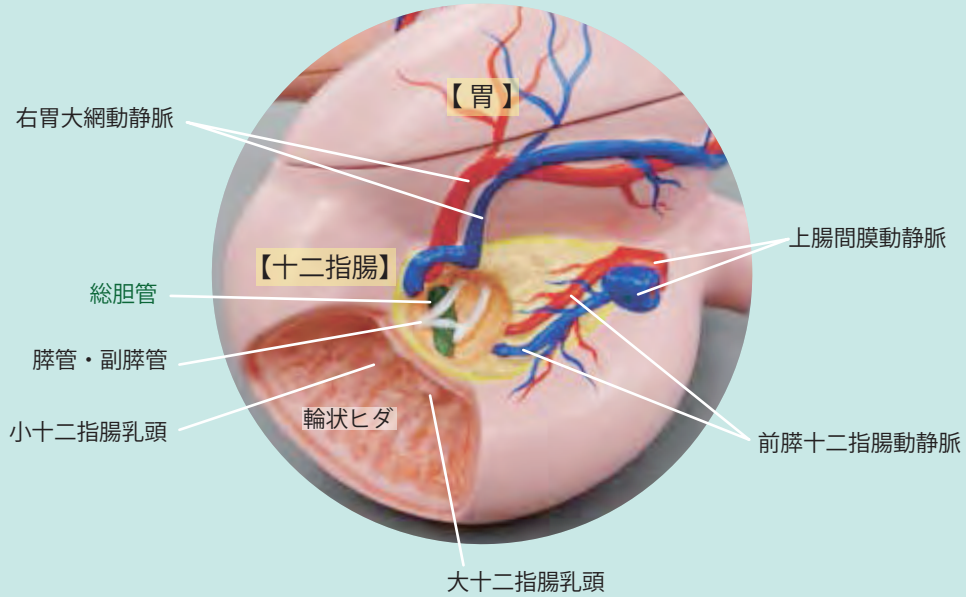
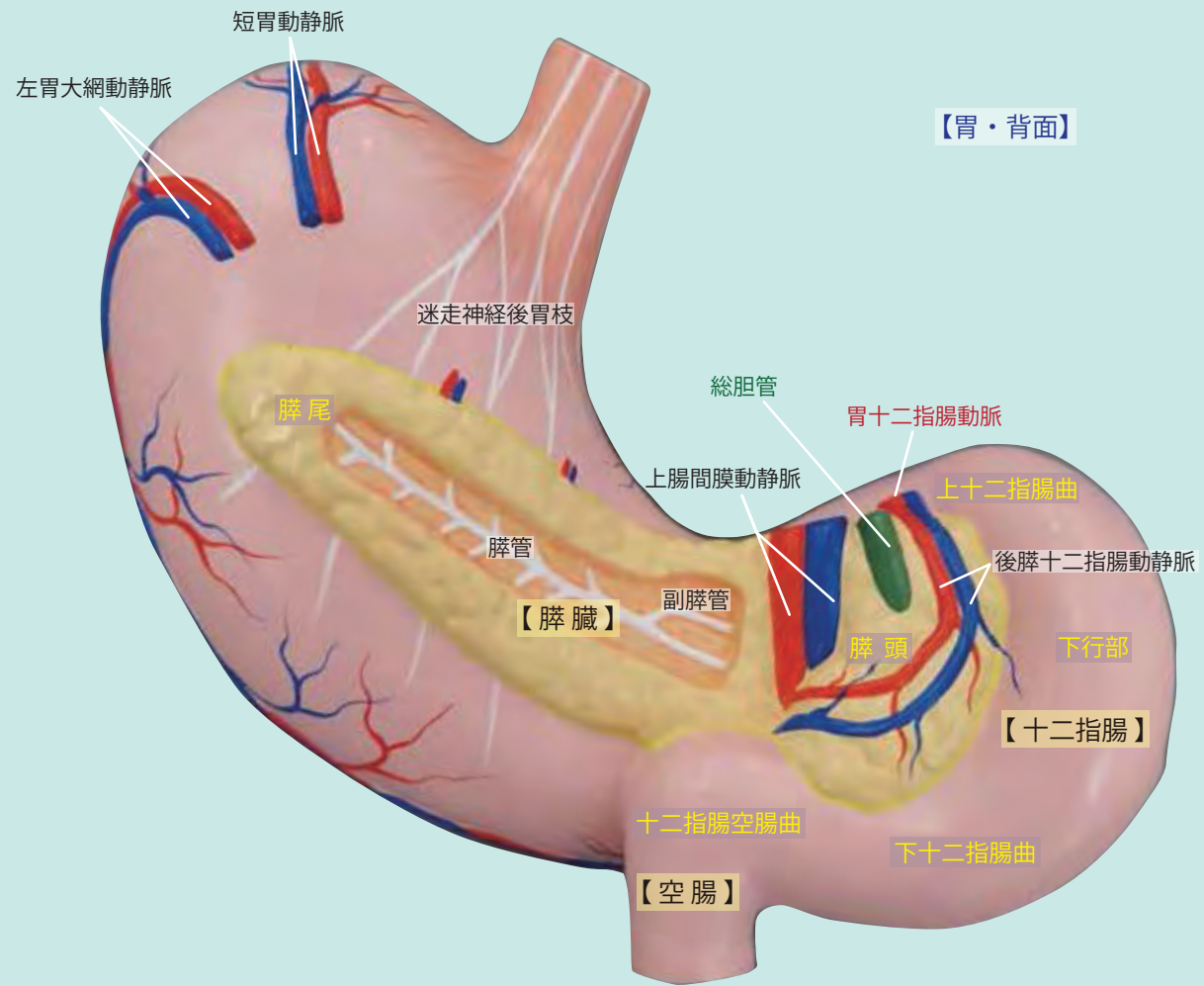
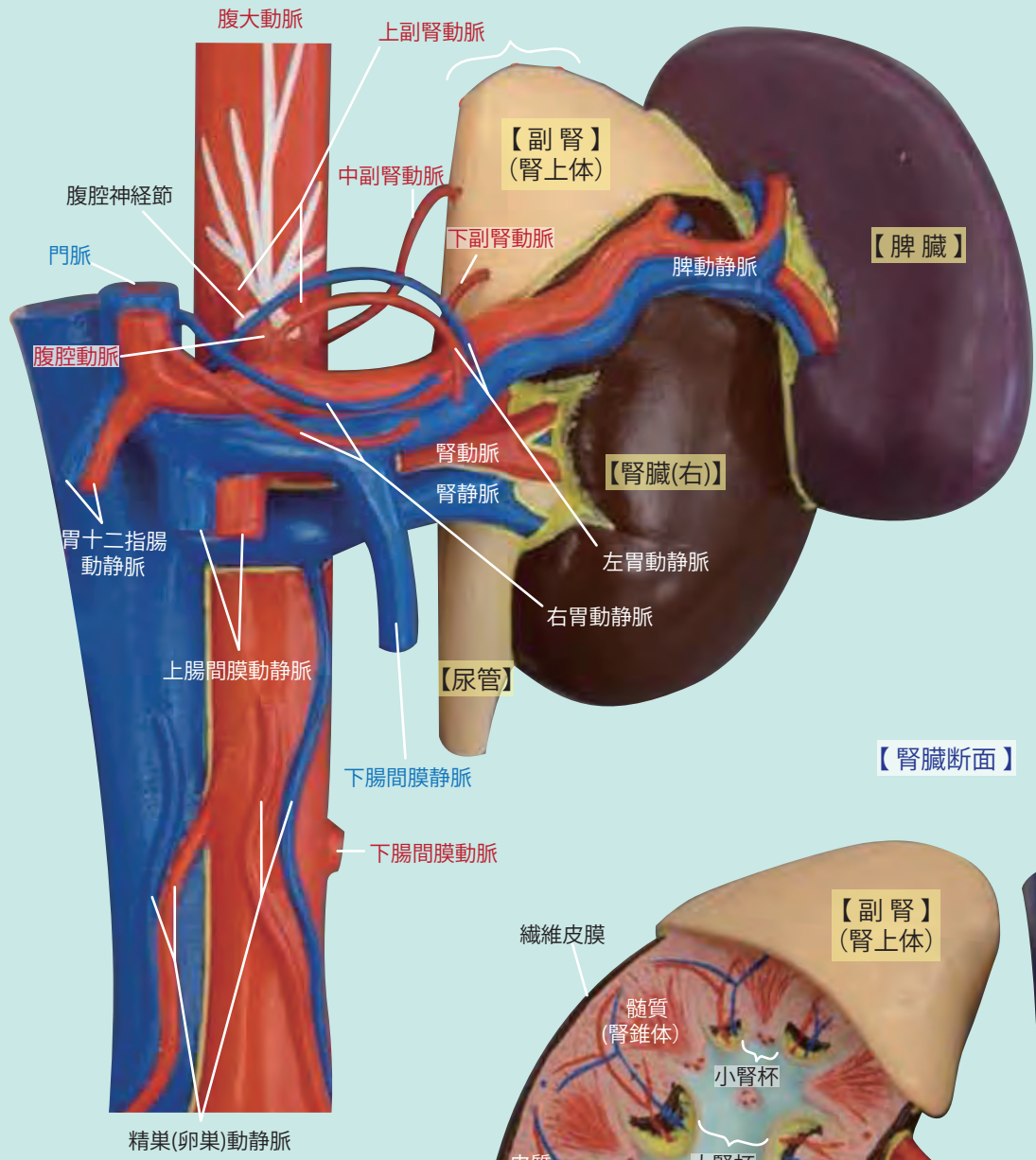
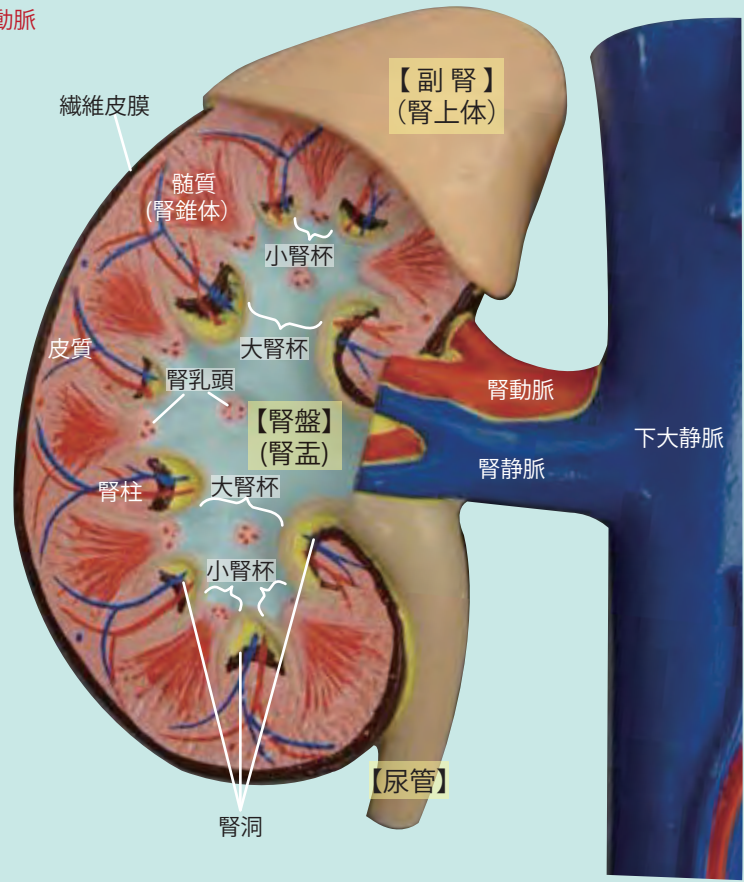
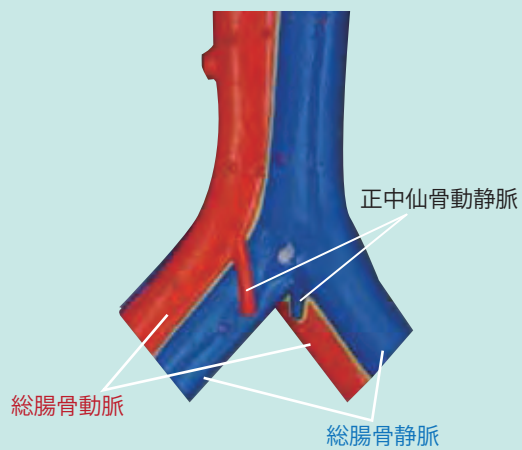
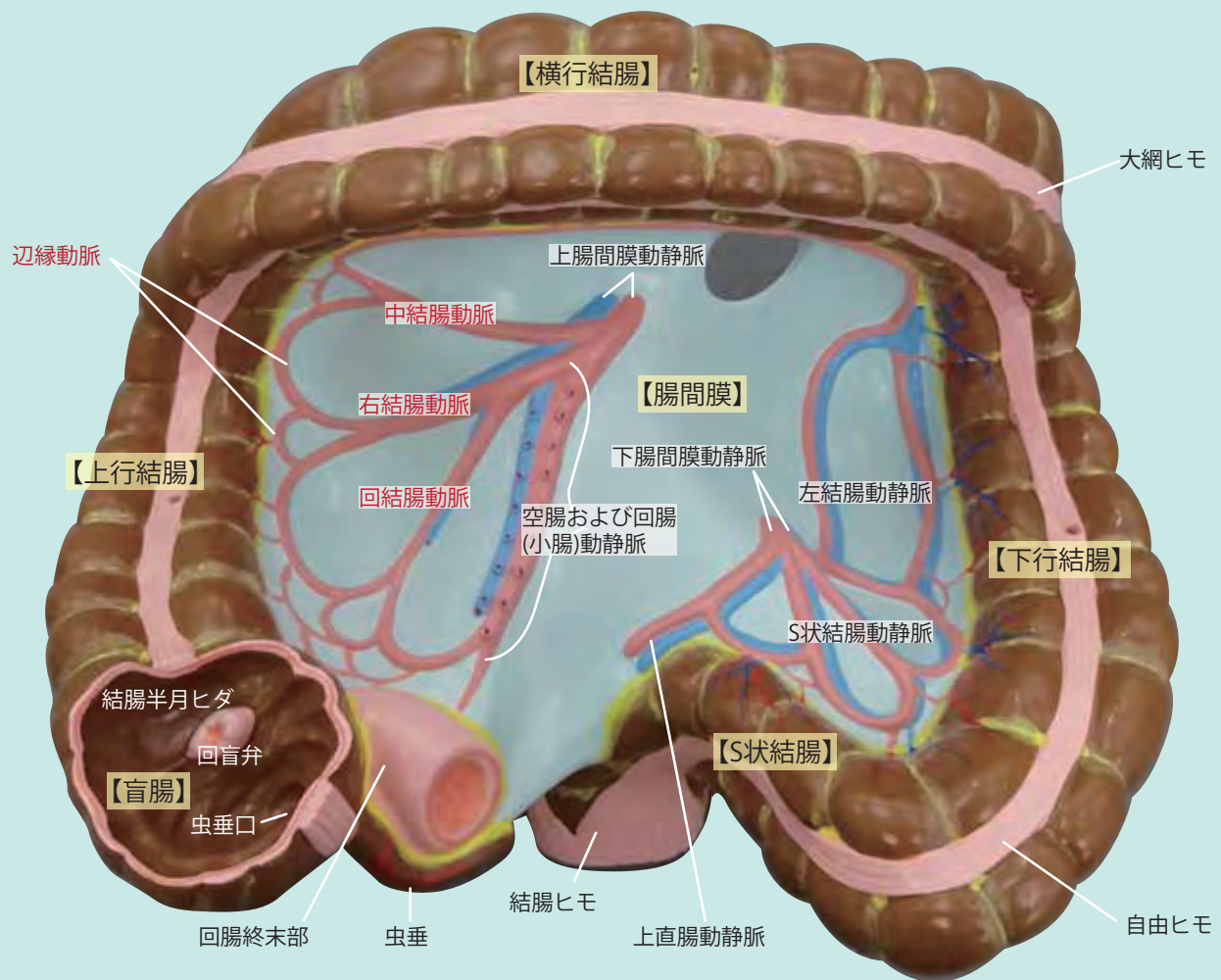
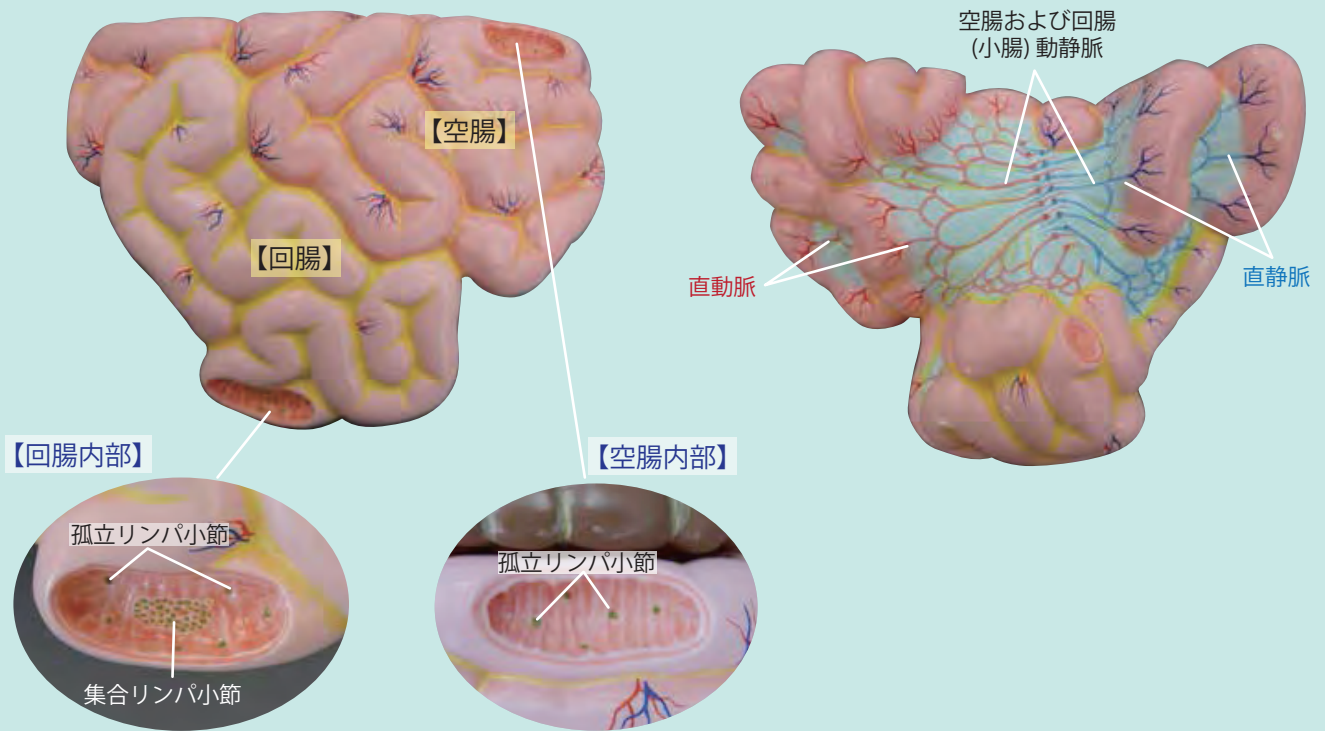
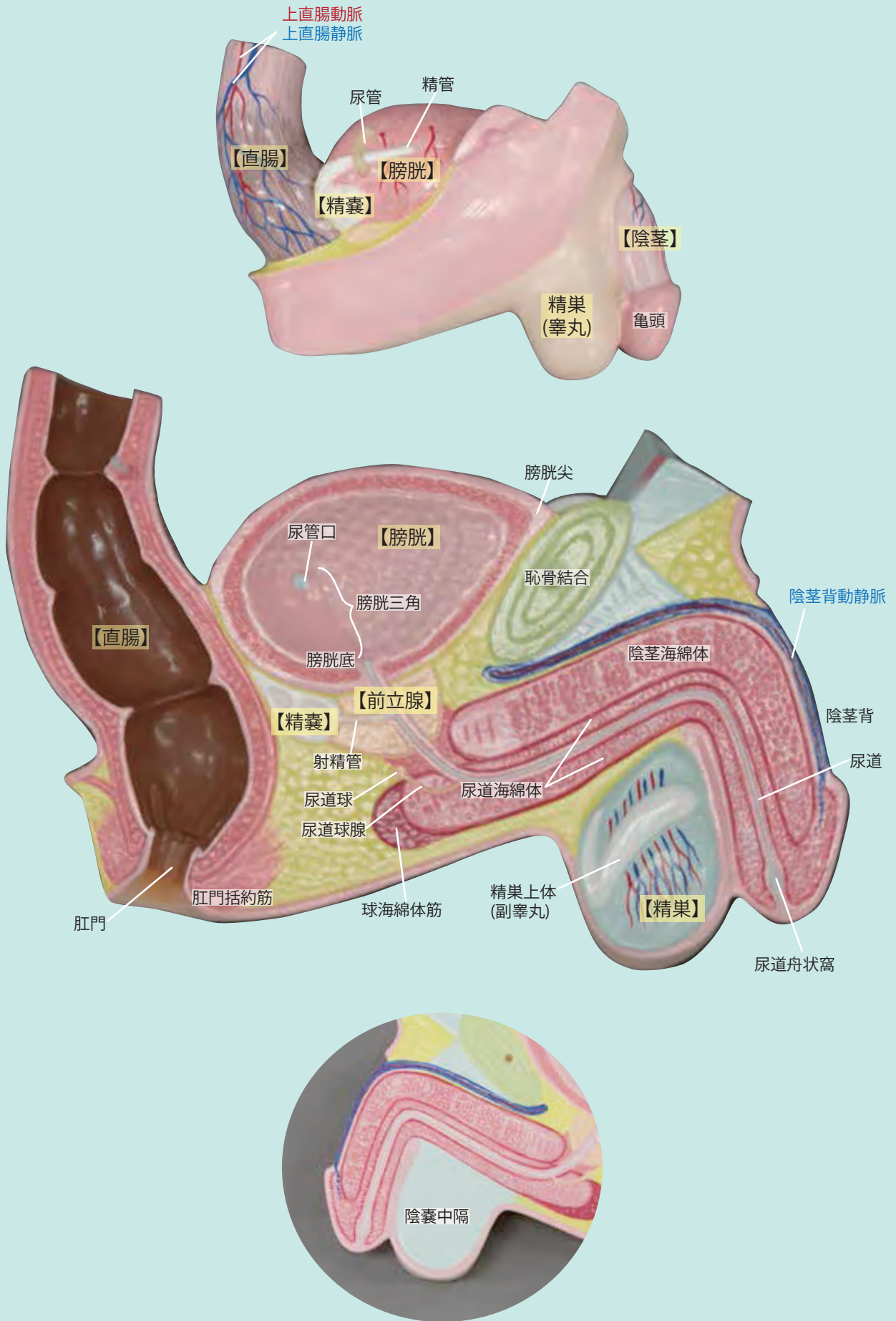
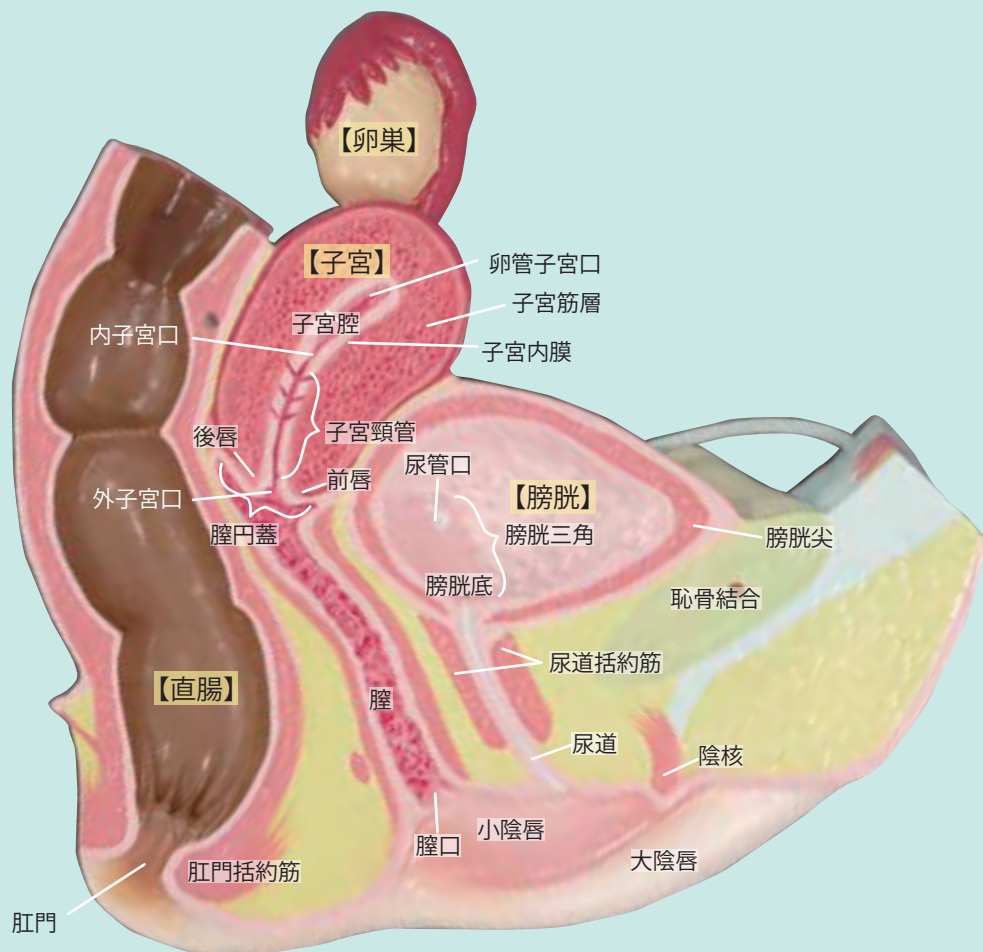
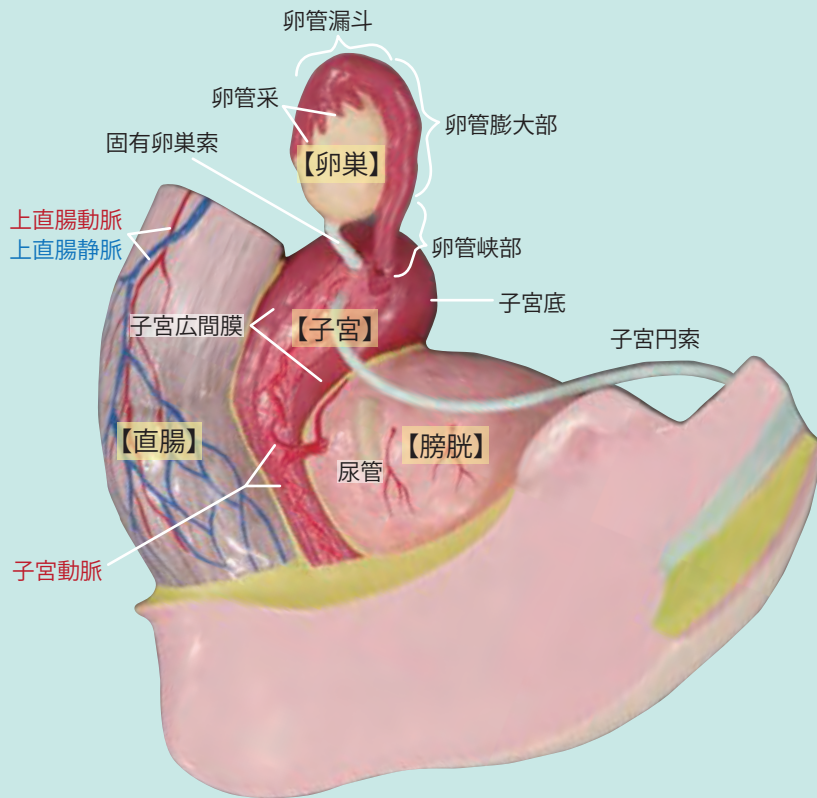
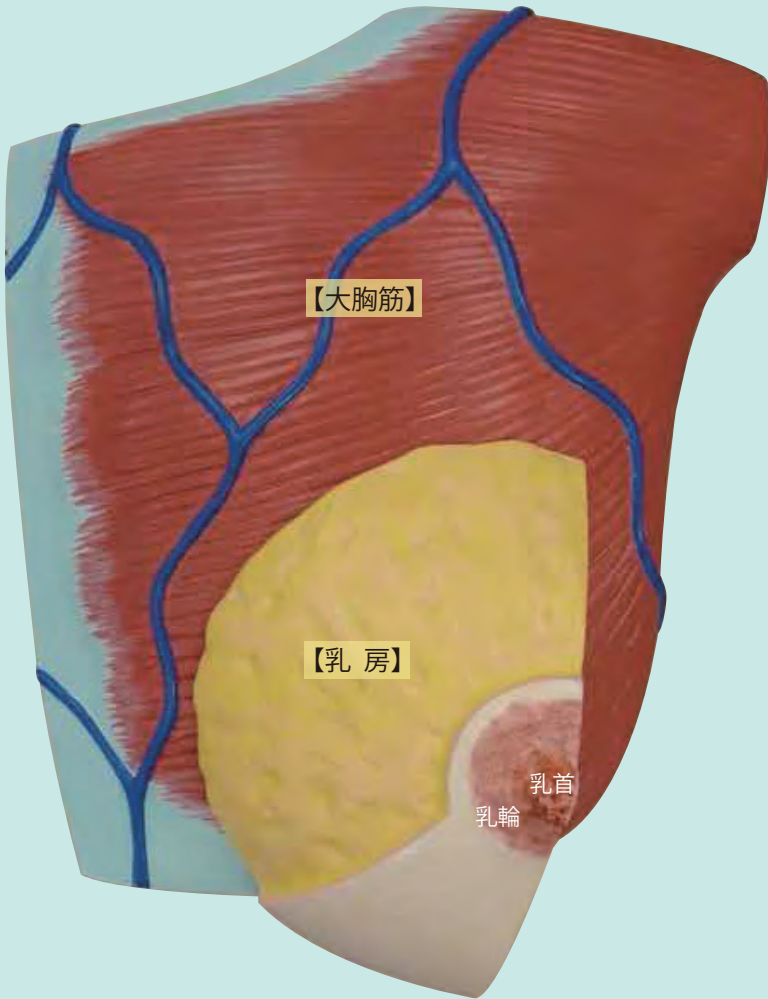
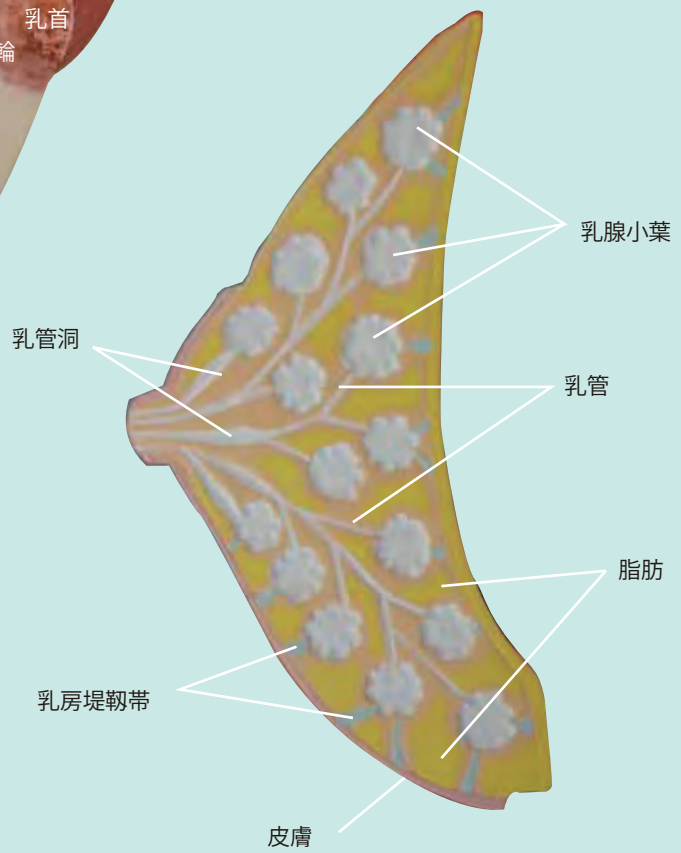
47	Front of head
48	Right side of head
49	Left side of head
50	Back of head / Front of neck
51	Inside skull base
52	Nasopharyngeal larynx
53	Cervical head section
54	Parts of the brain
55	Midsection of the brain
56	Brain cross section / Bottom surface
57	Chest / upper arm
58	Upper body back
59	Abdomen / Thigh
60	Buttocks / Thigh
61	Chest-abdominal wall (front)
62	Chest wall (Posterior)
63	Abdominal wall (Posterior)
64	Pelvic cavity
65	Lumbar
66	Lung
67	Trachea / esophagus
68	Heart
69	Atria / Ventricles
70	Diaphragm
71	Liver
72	Stomach
73	Pancreas / Duodenum
74	Kidney
75	Intestines
76	Male reproductive organs
77	Female reproductive organs
78	Breast

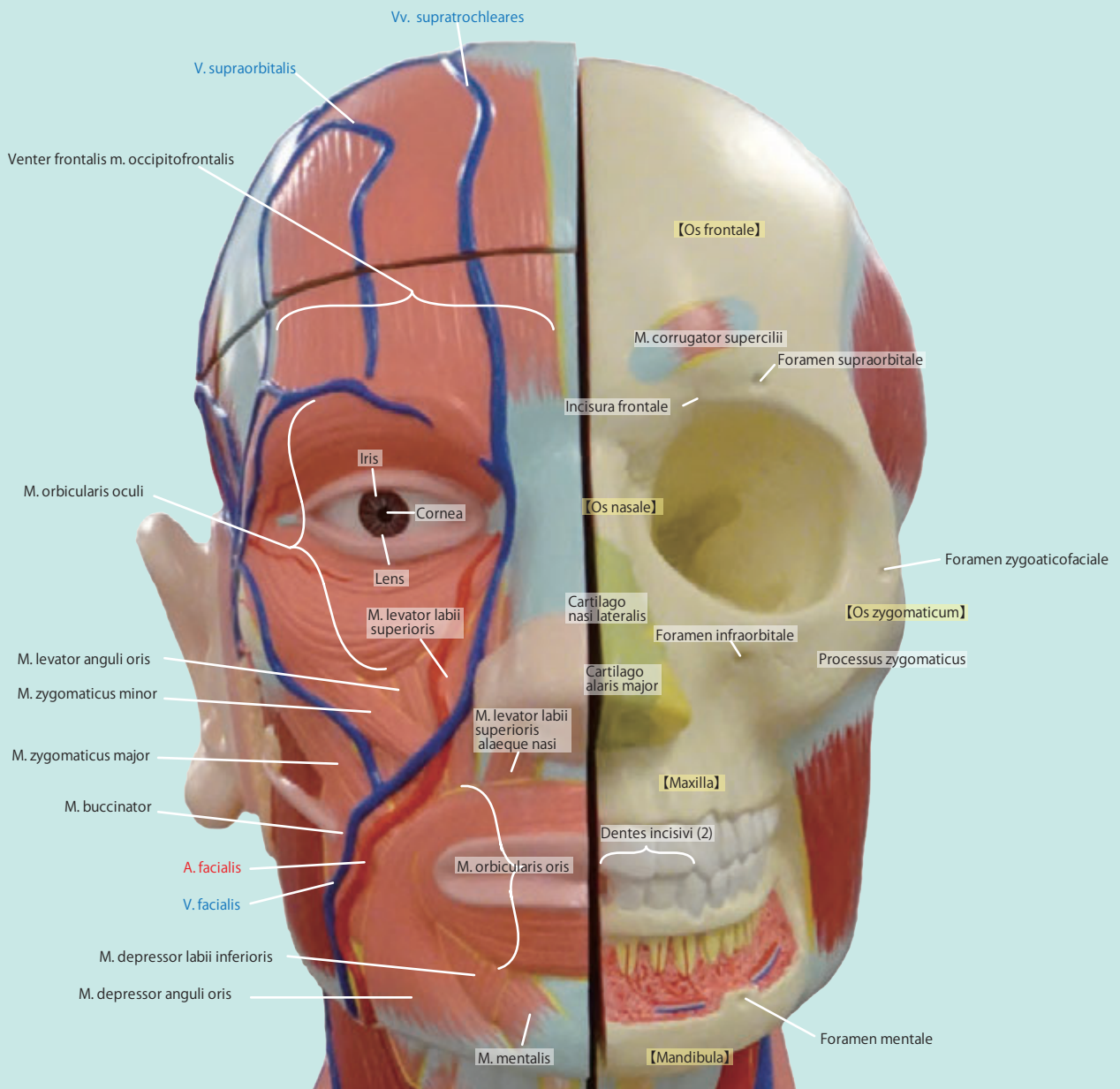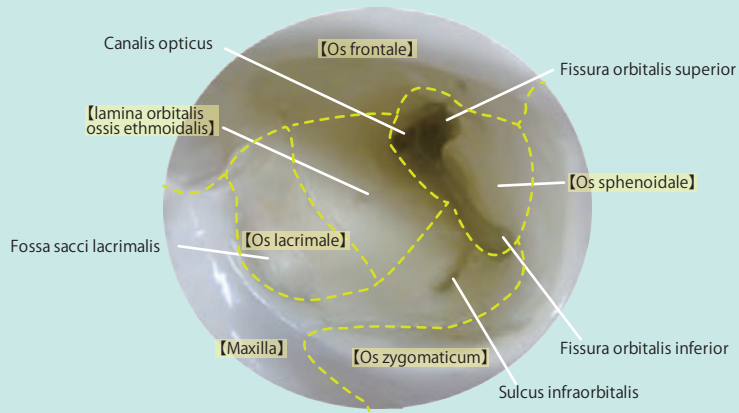
Front of head



Right side of head

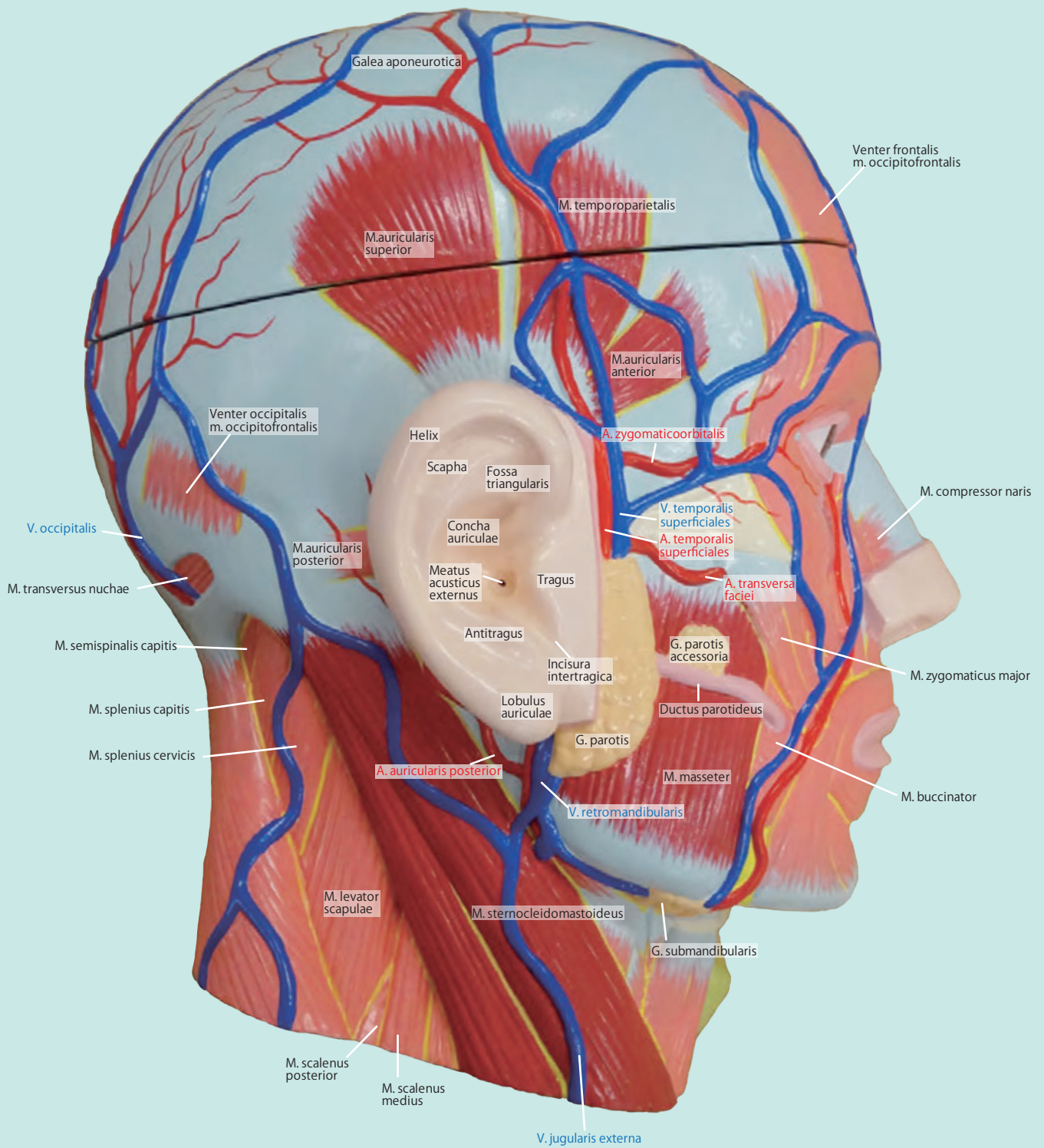
Left side of head

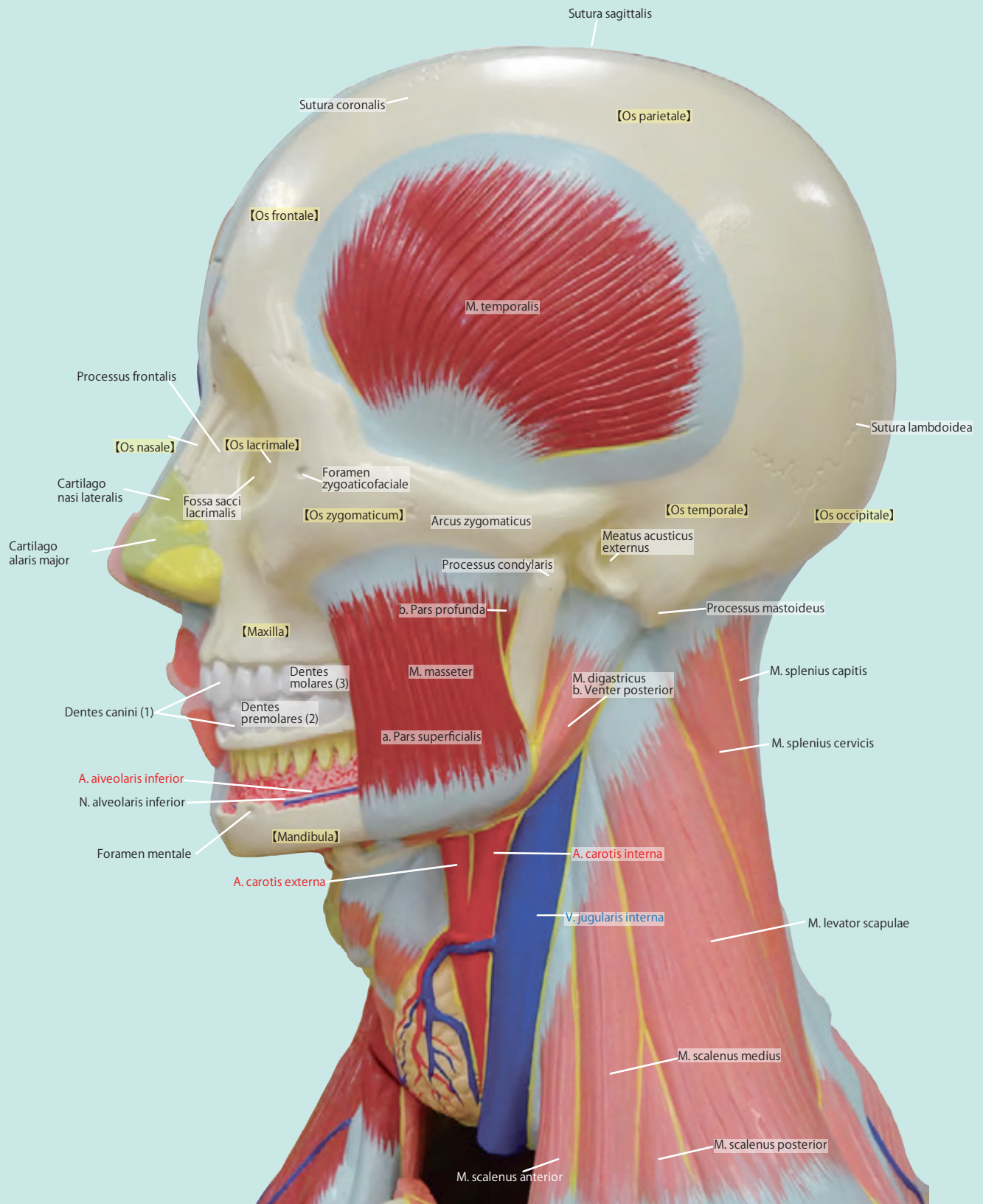
Back of head / Front of neck

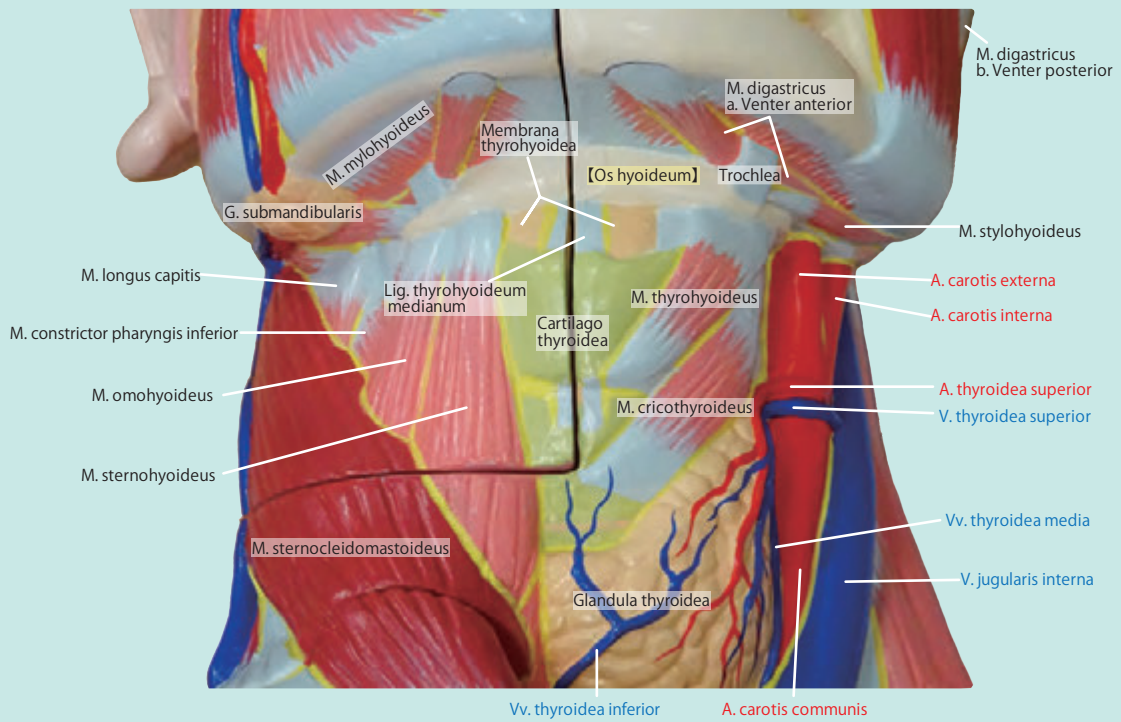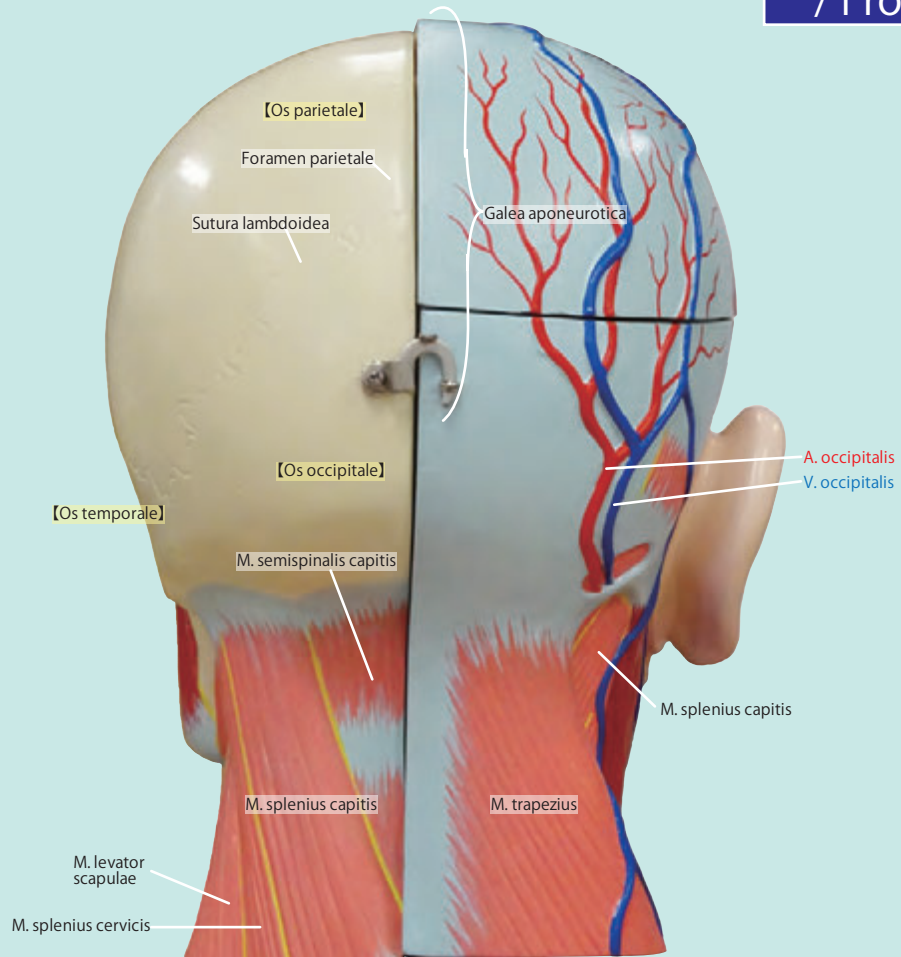
Inside skull base




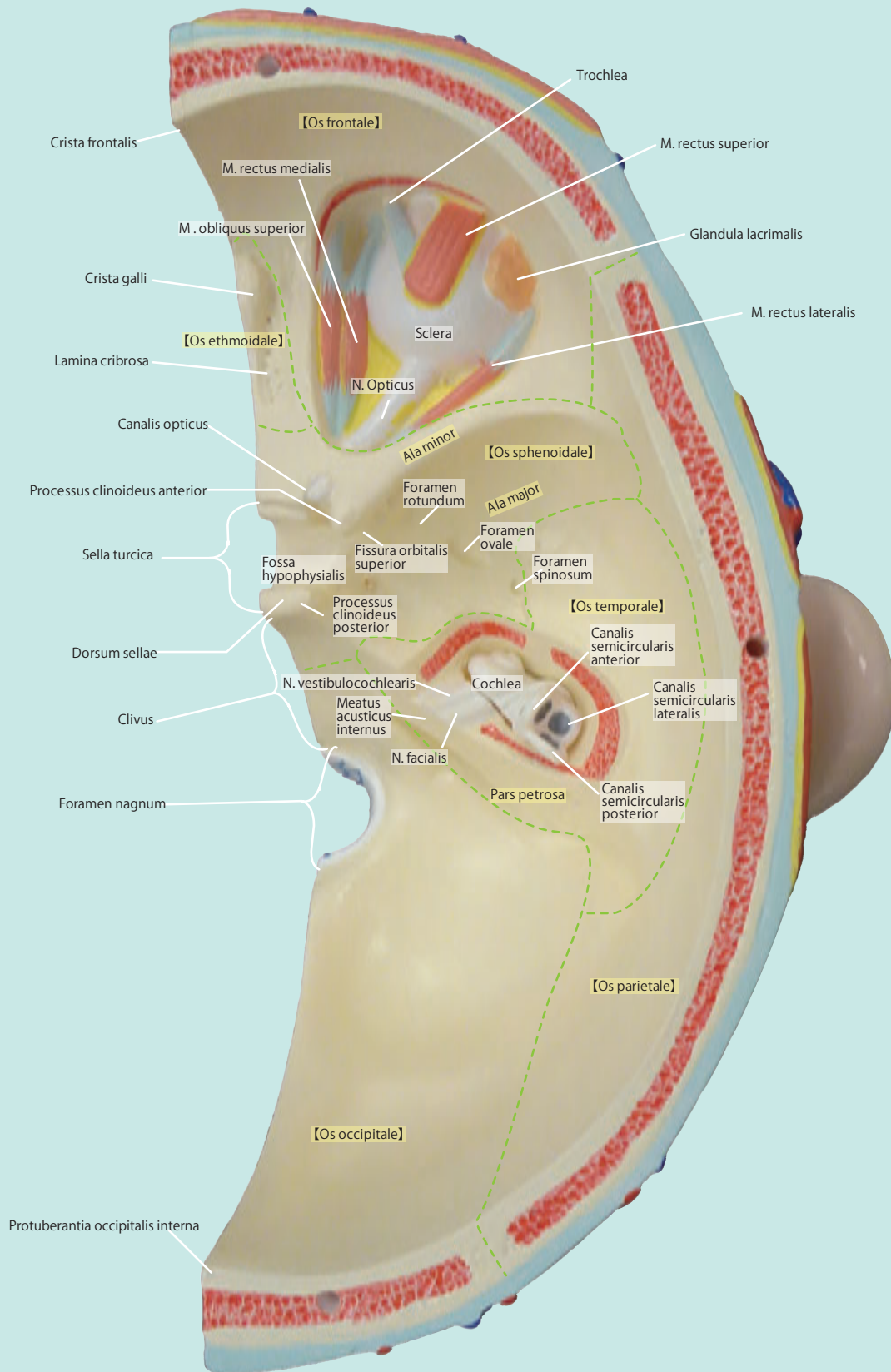
Nasopharyngeal larynx

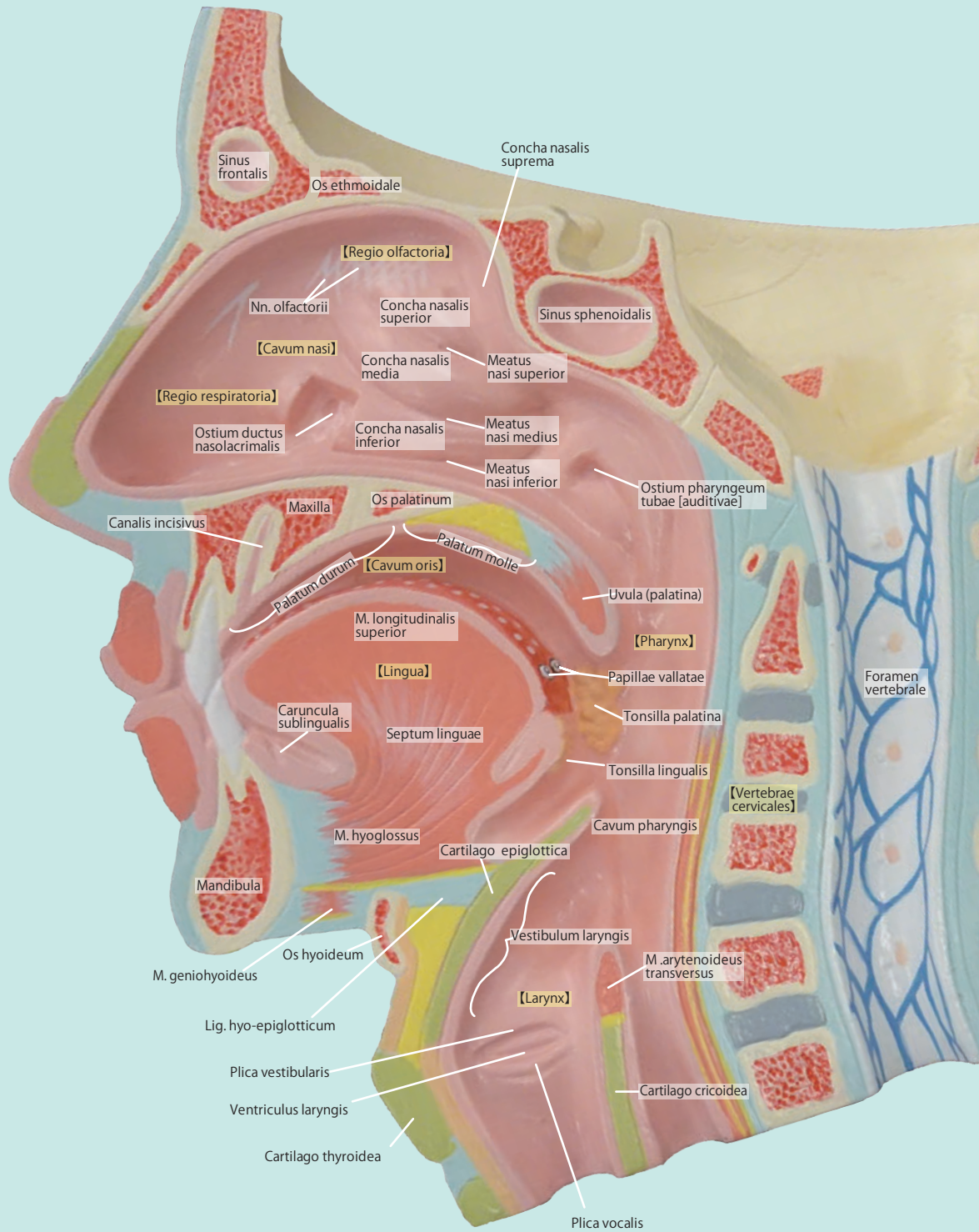
Cervical head section

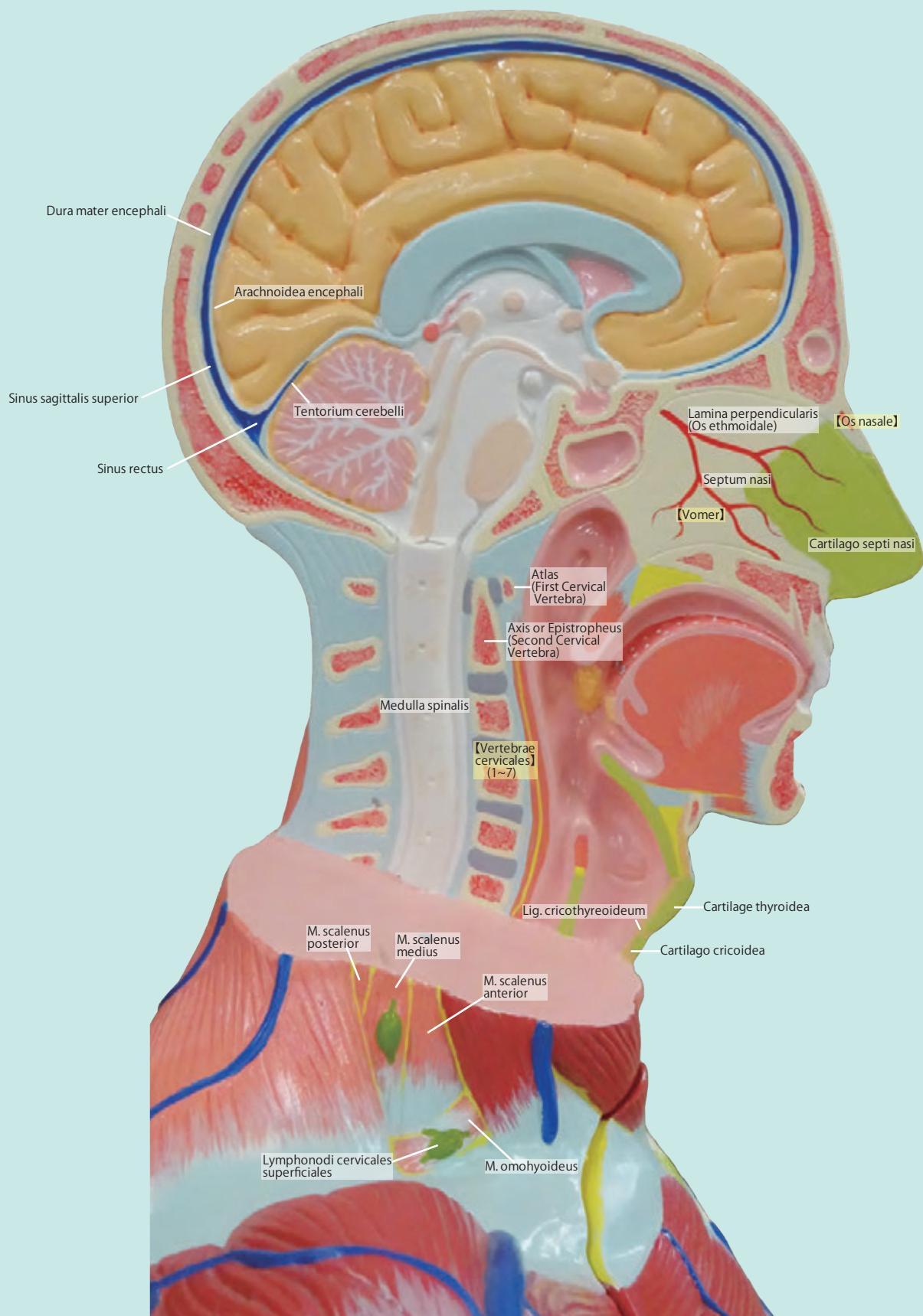Parts of the brain

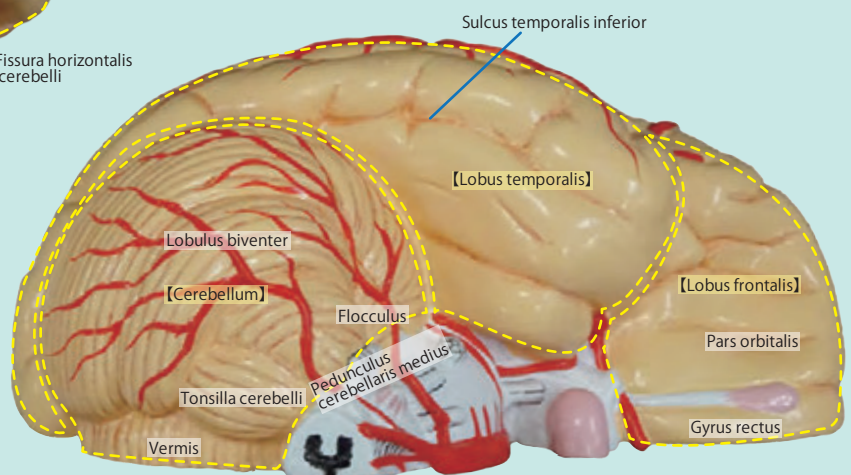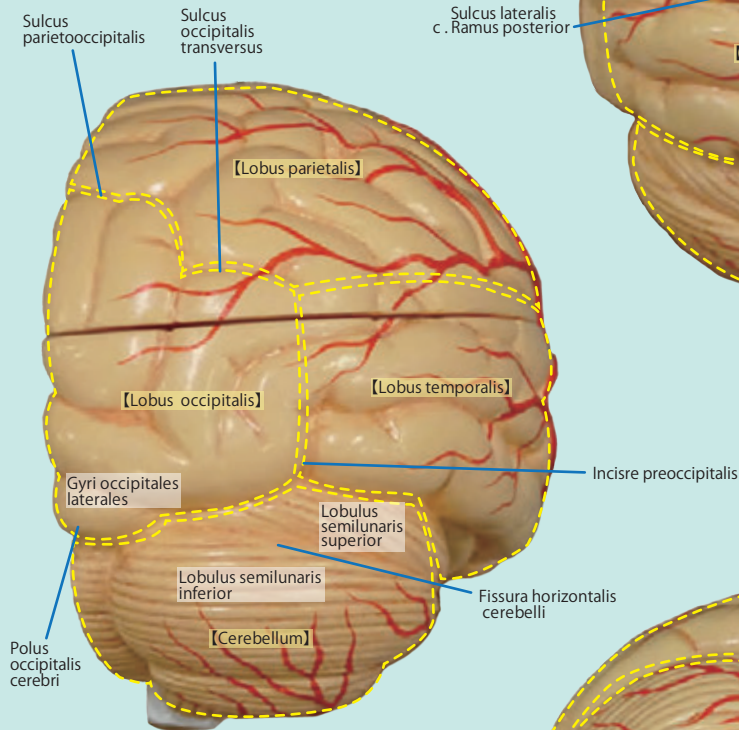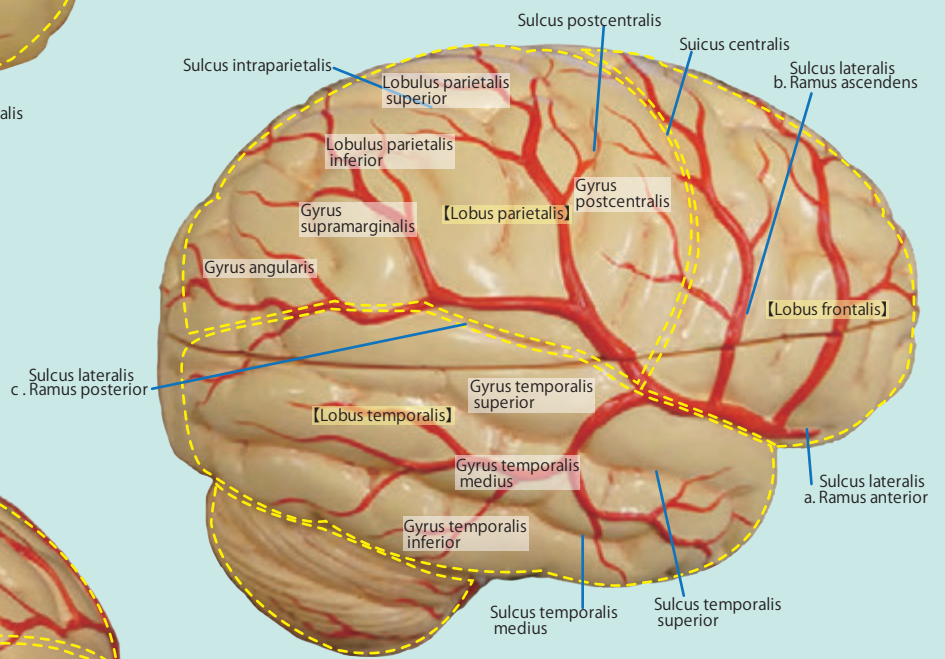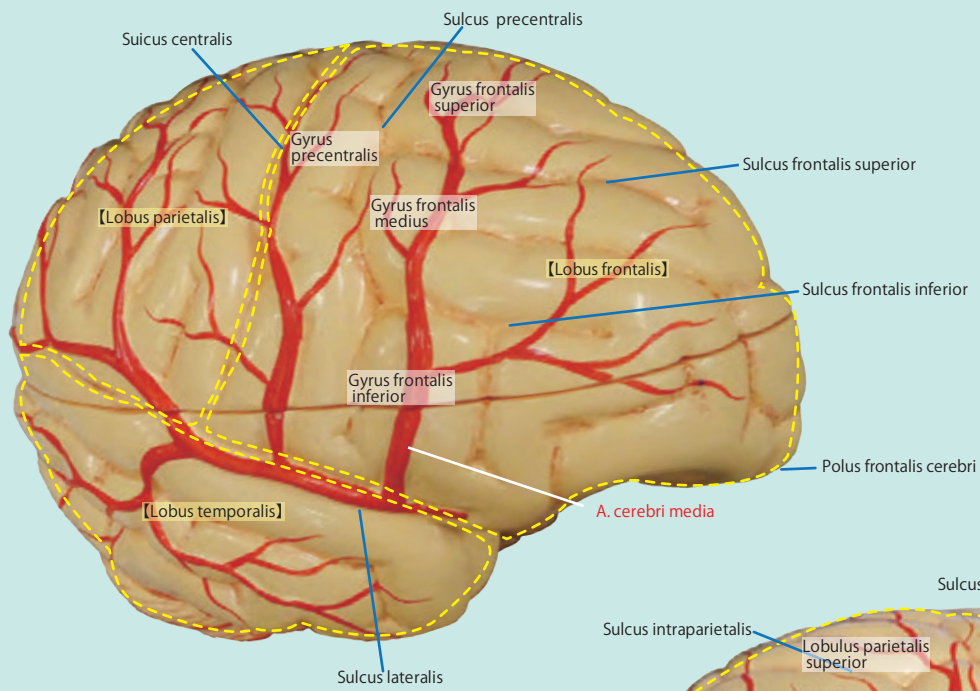
Midsection of the brain

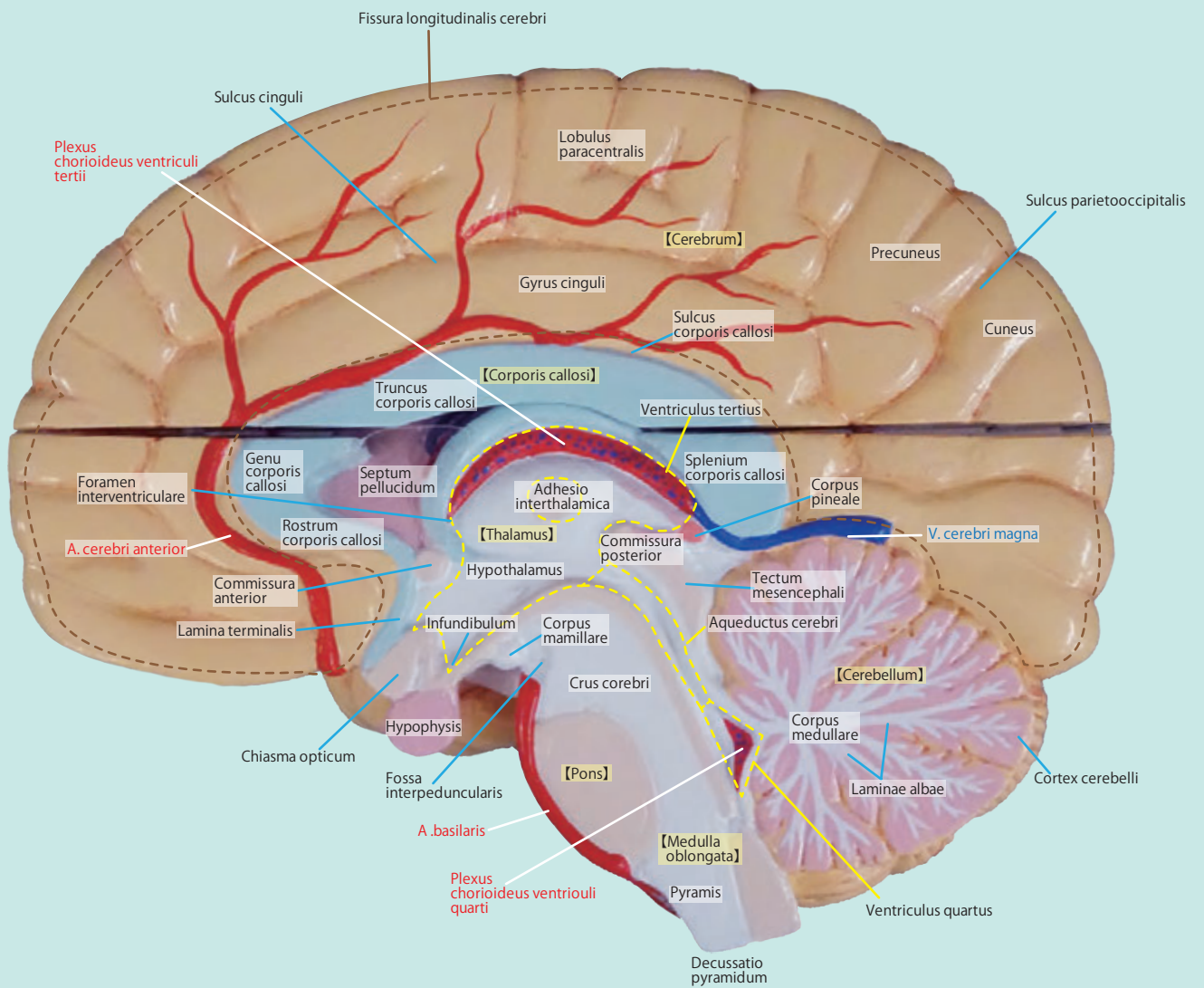
Brain cross section /Bottom surface

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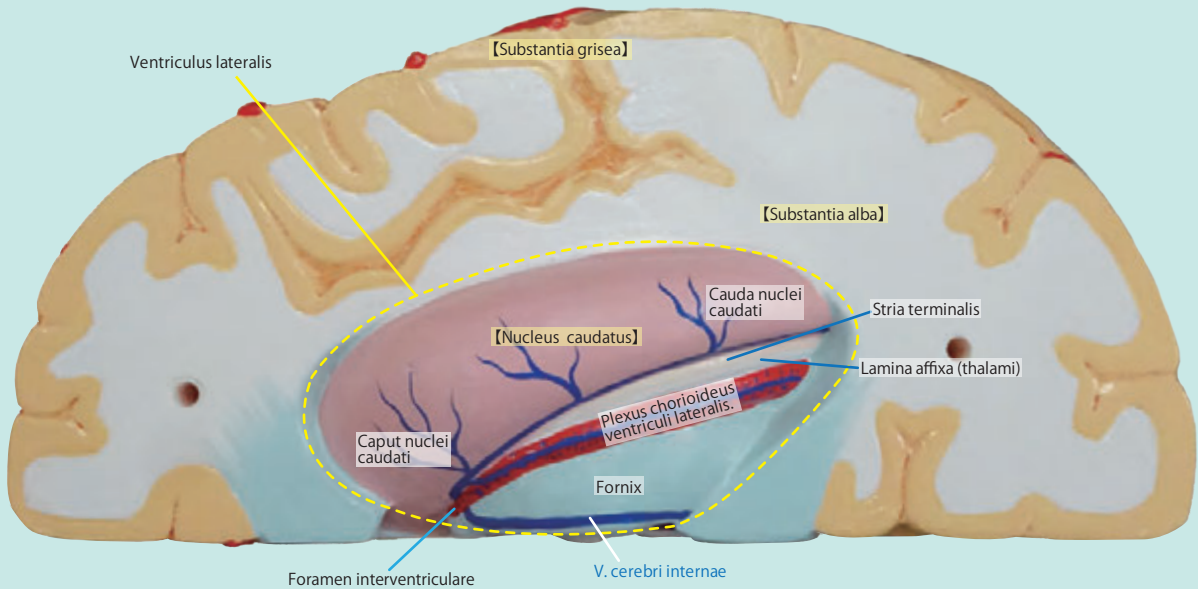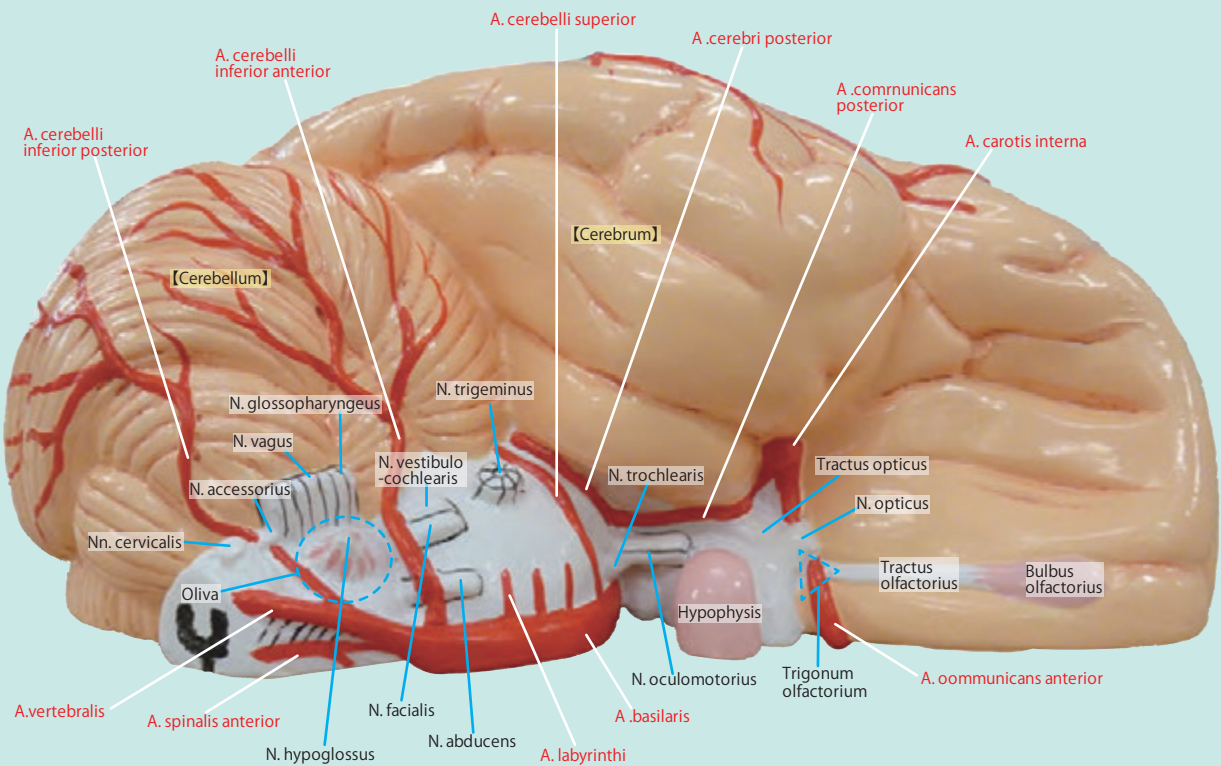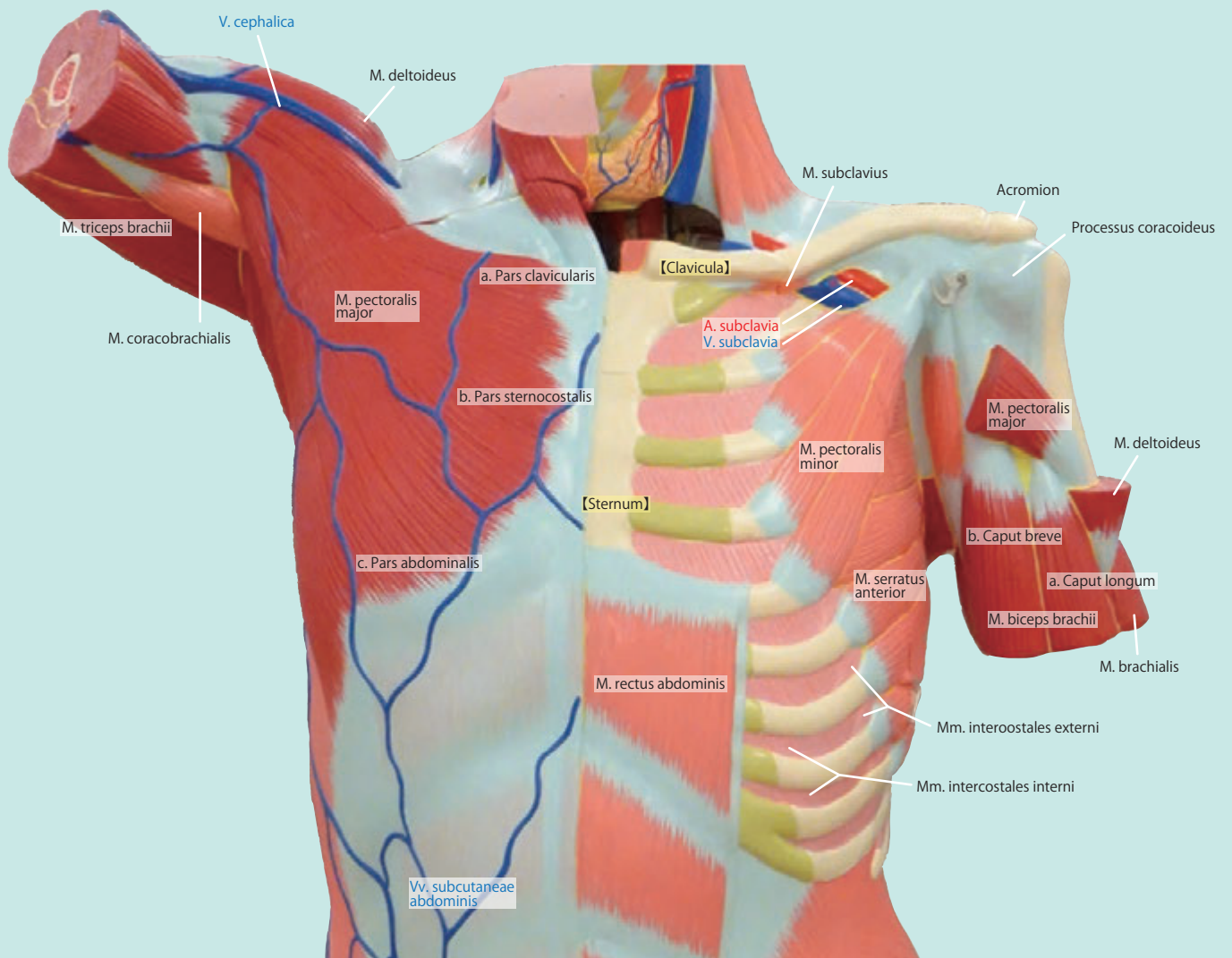
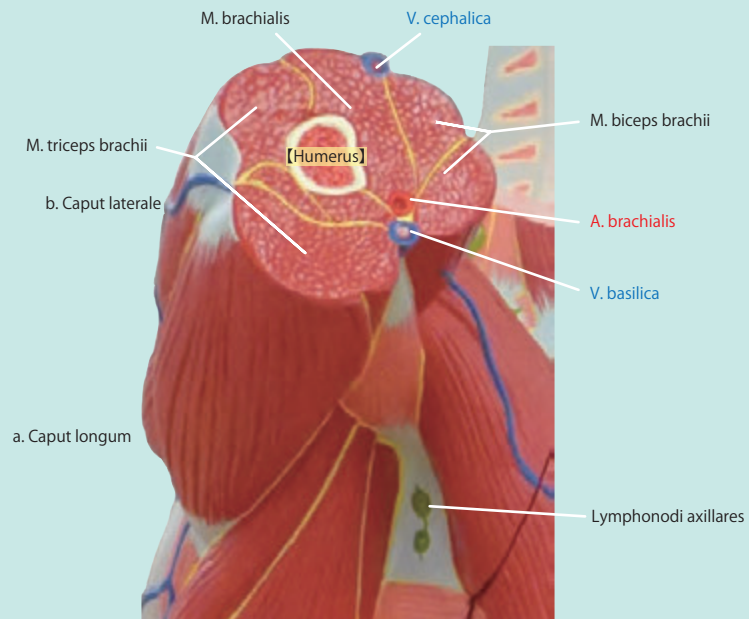
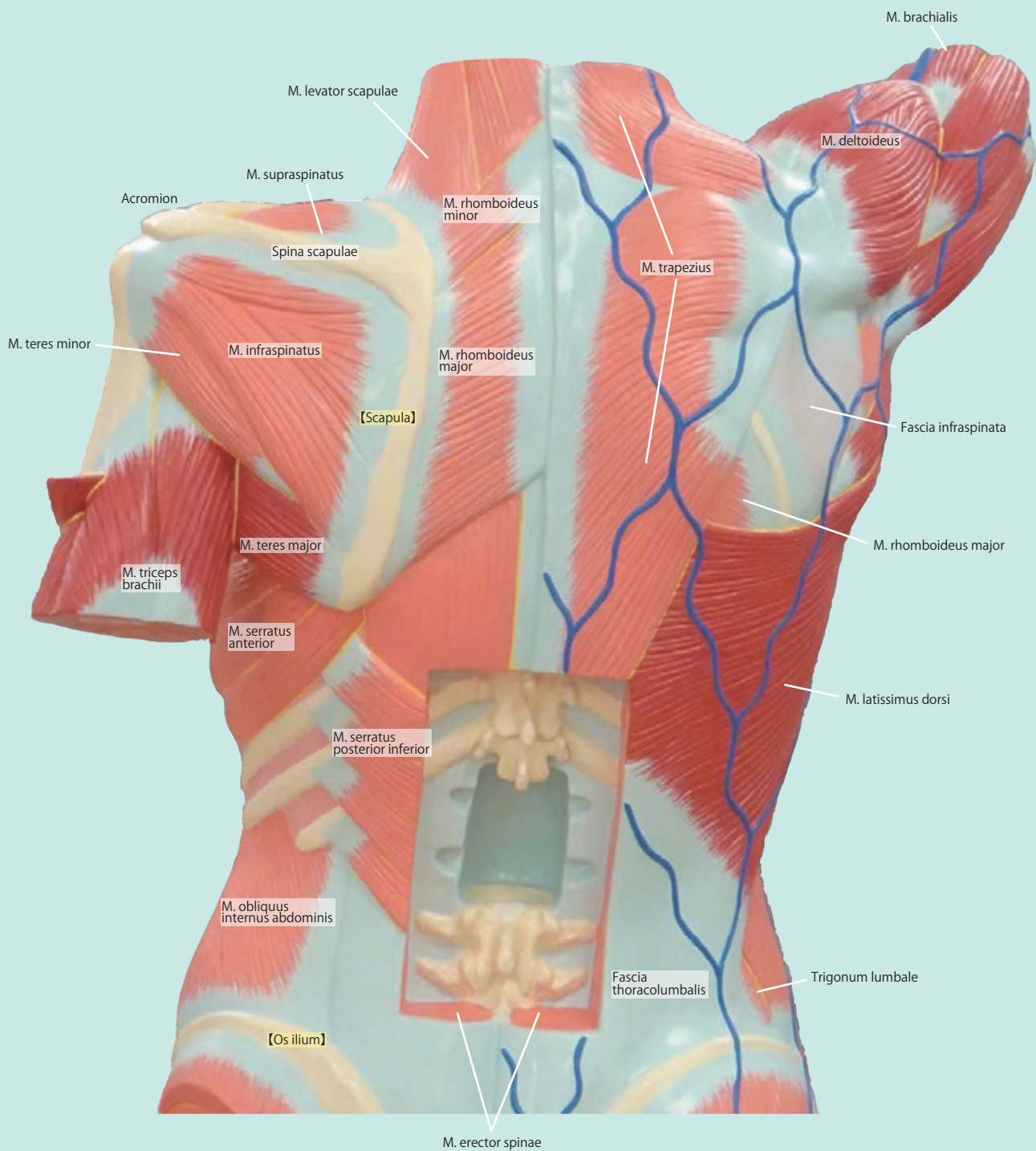
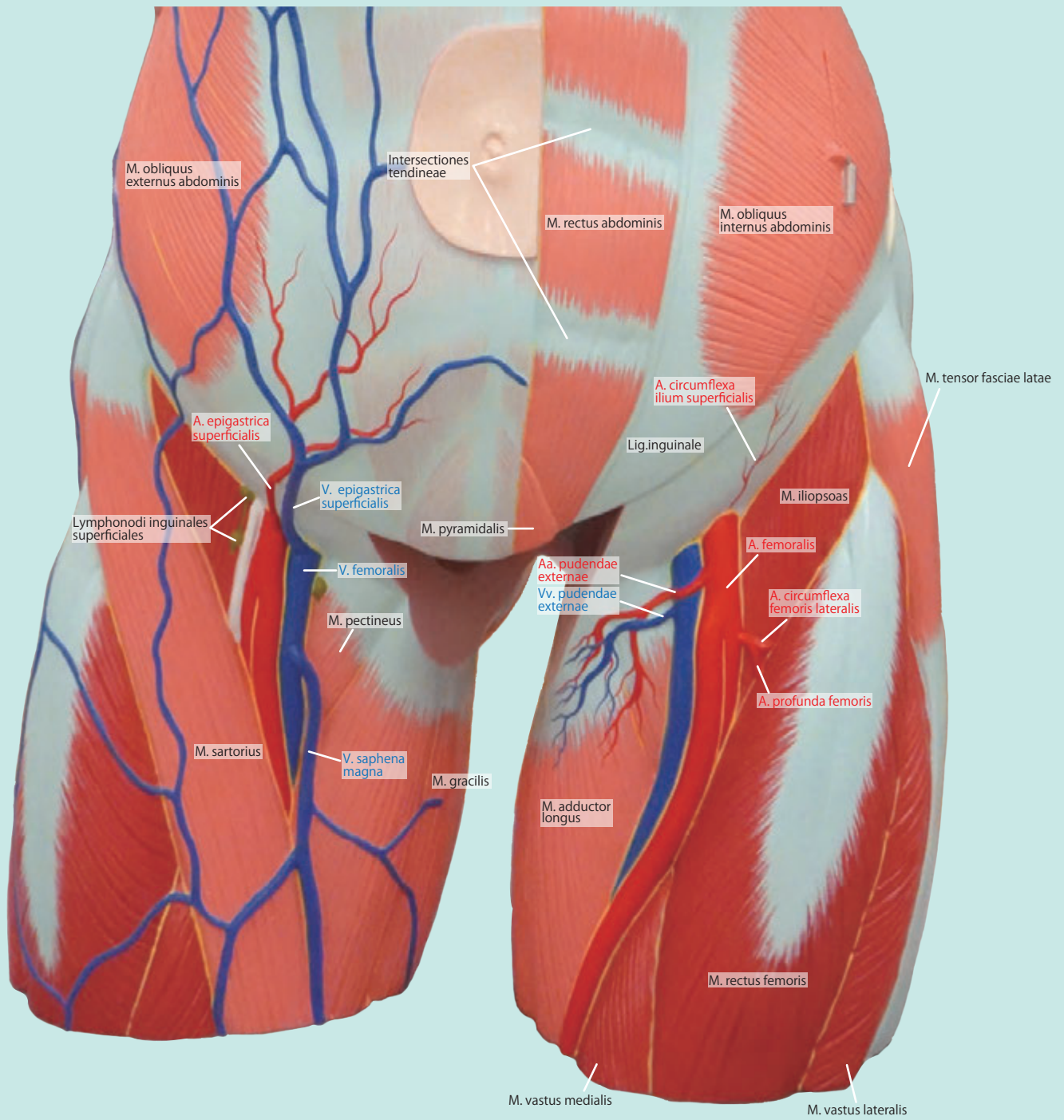
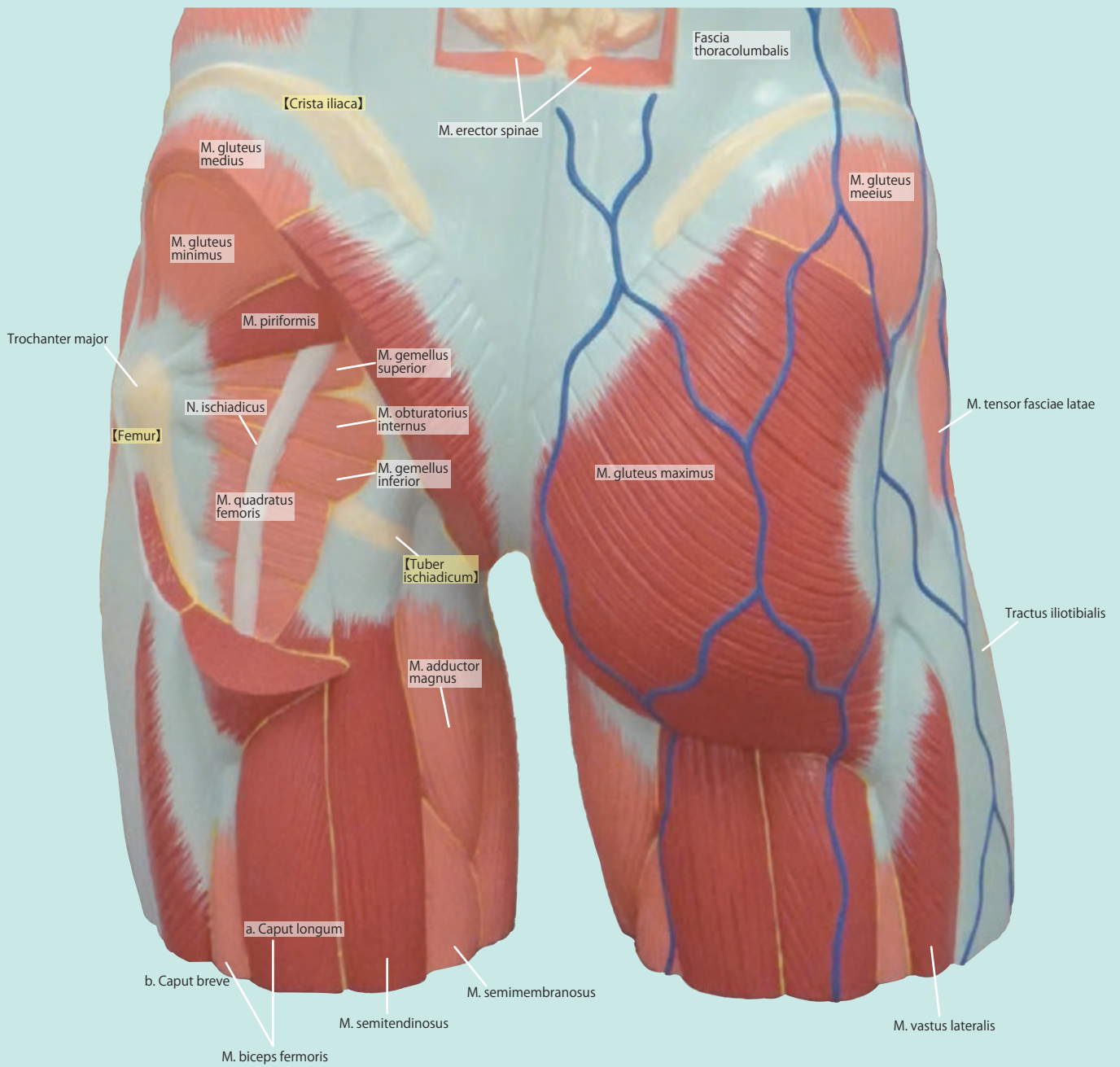
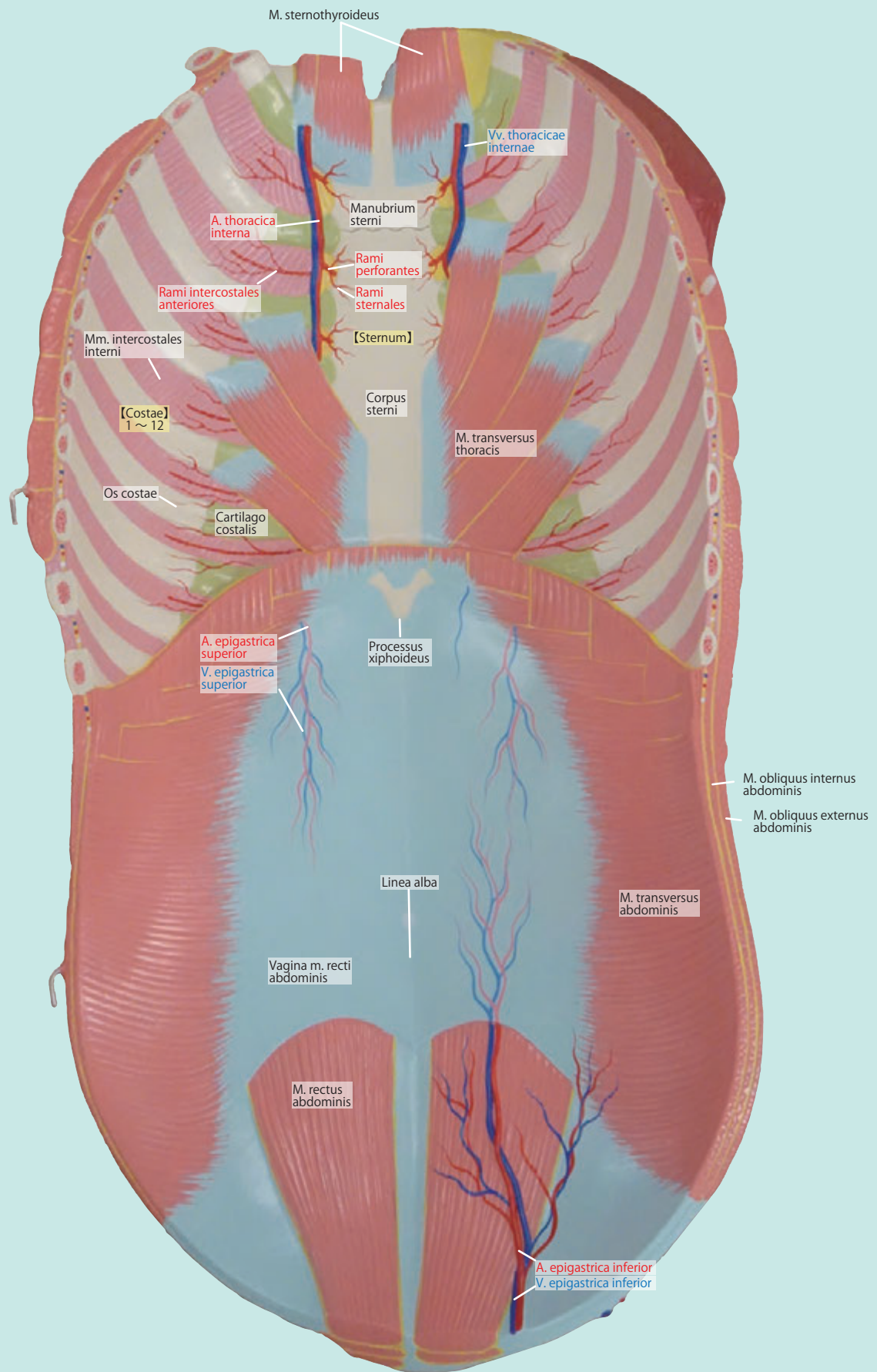
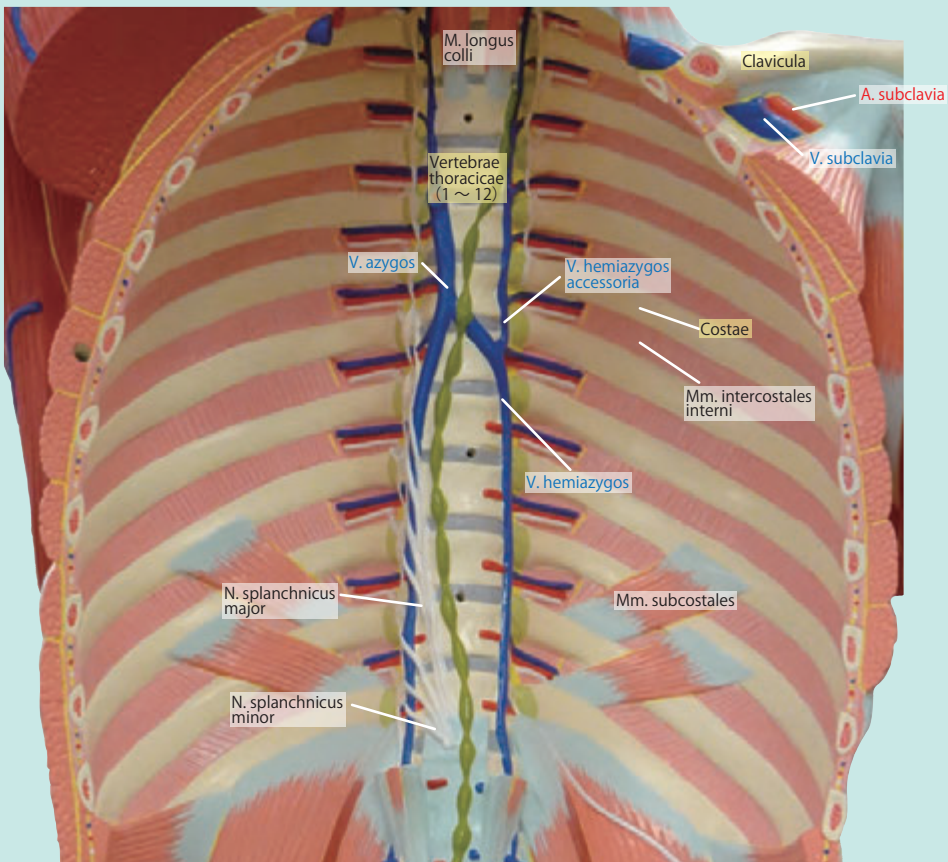
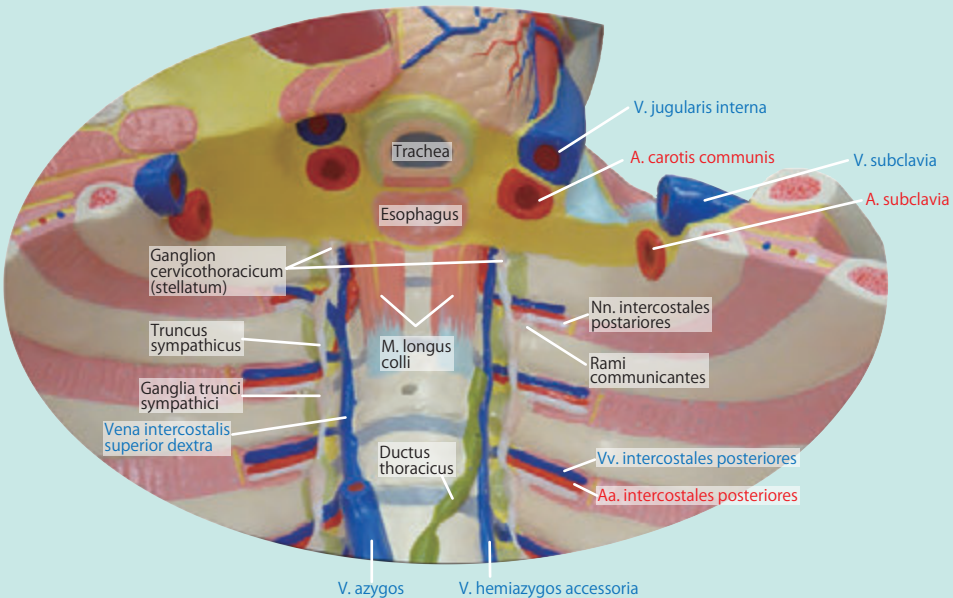
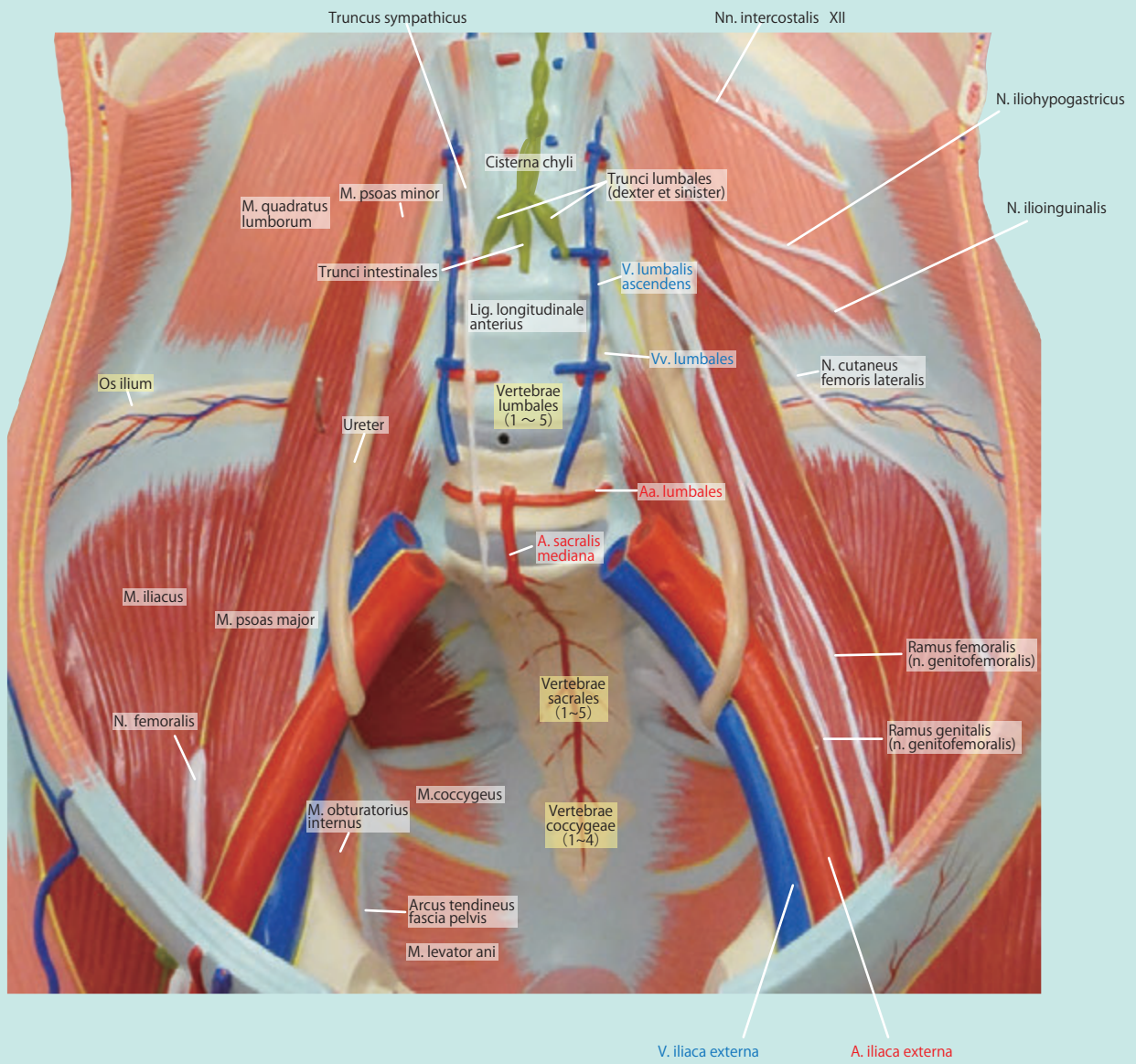
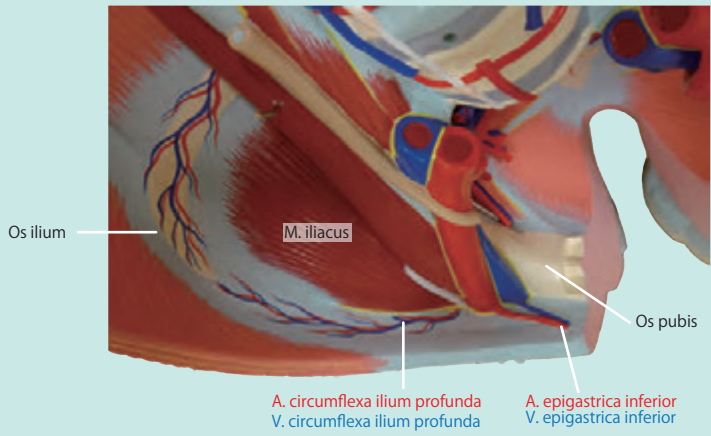
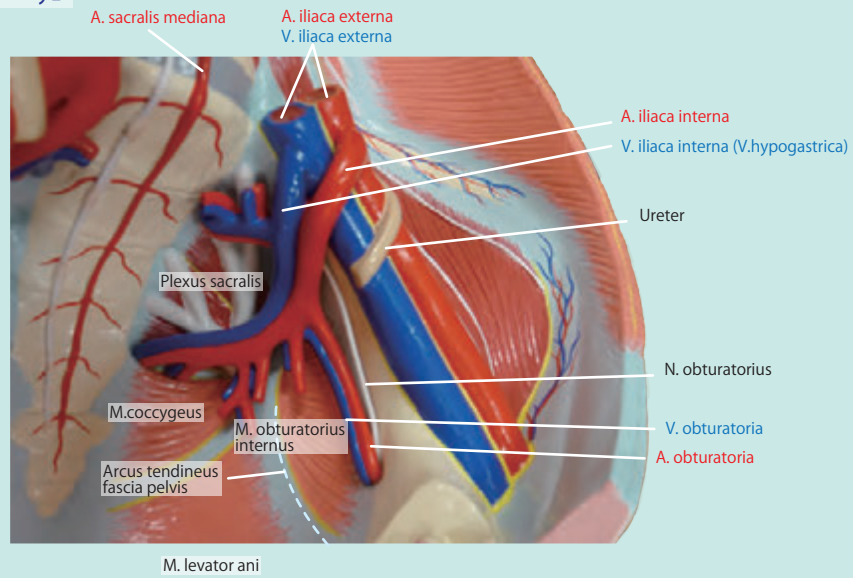
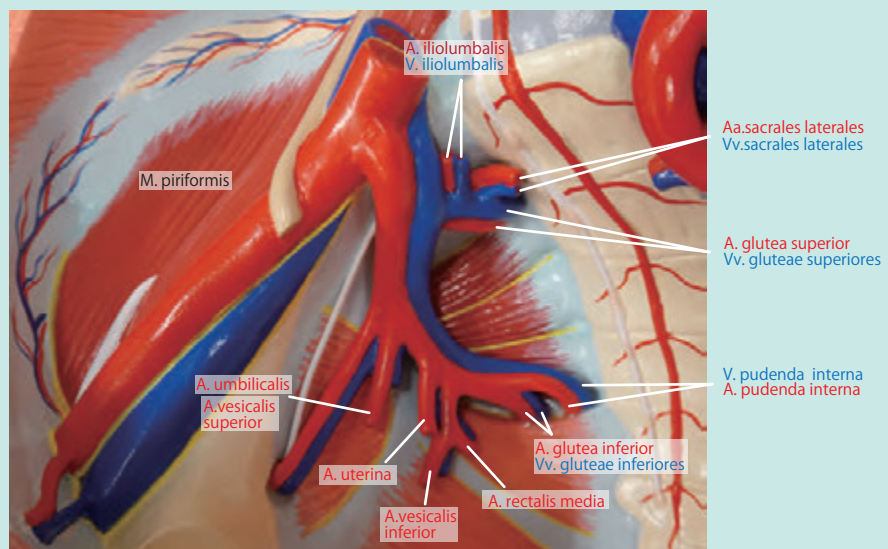
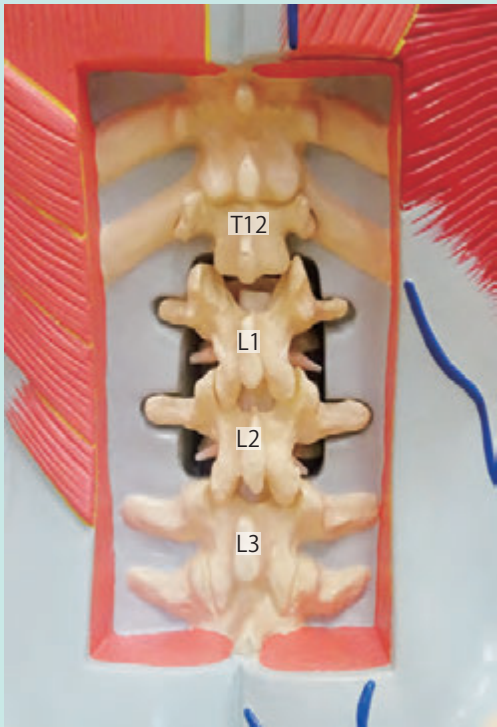
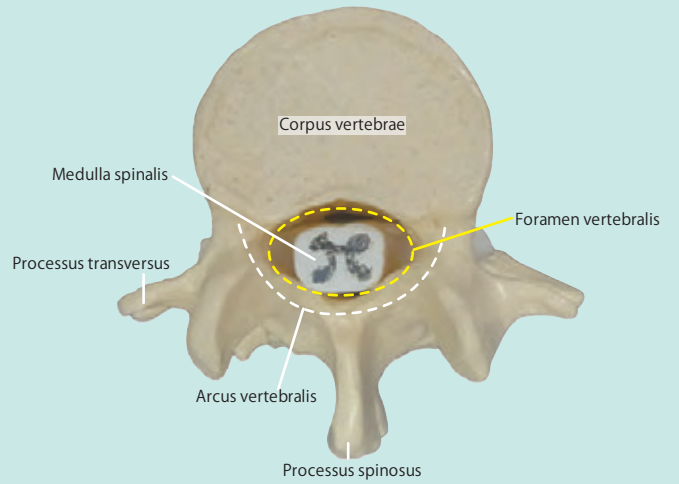
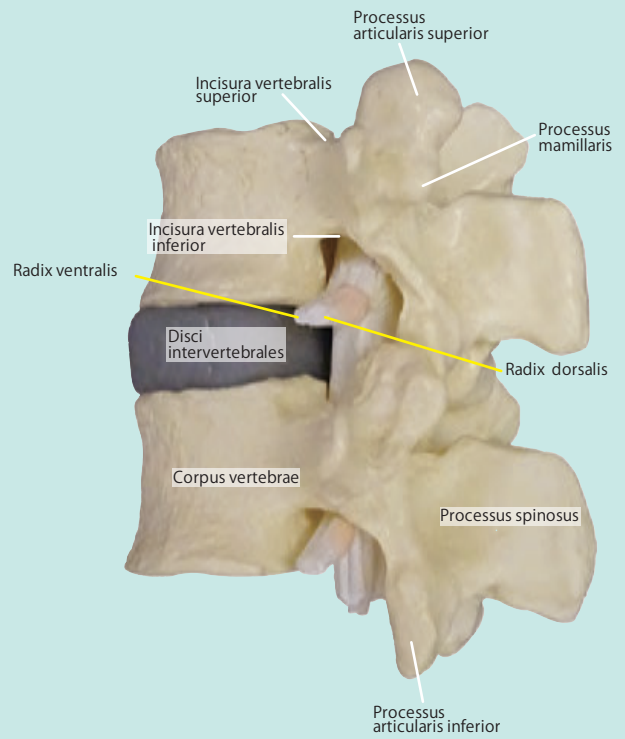
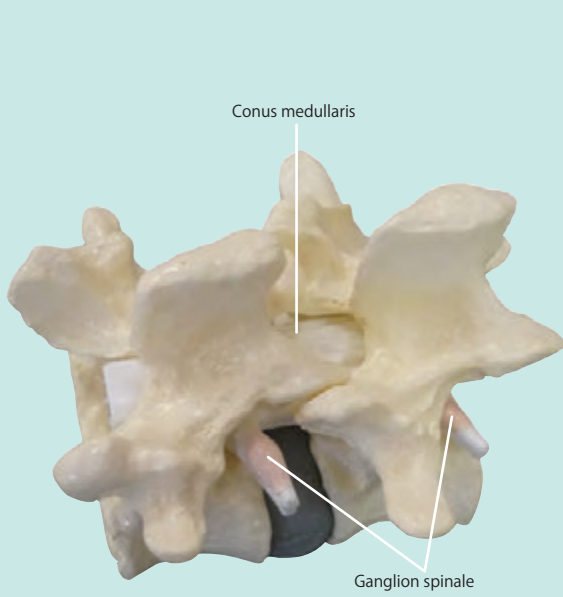
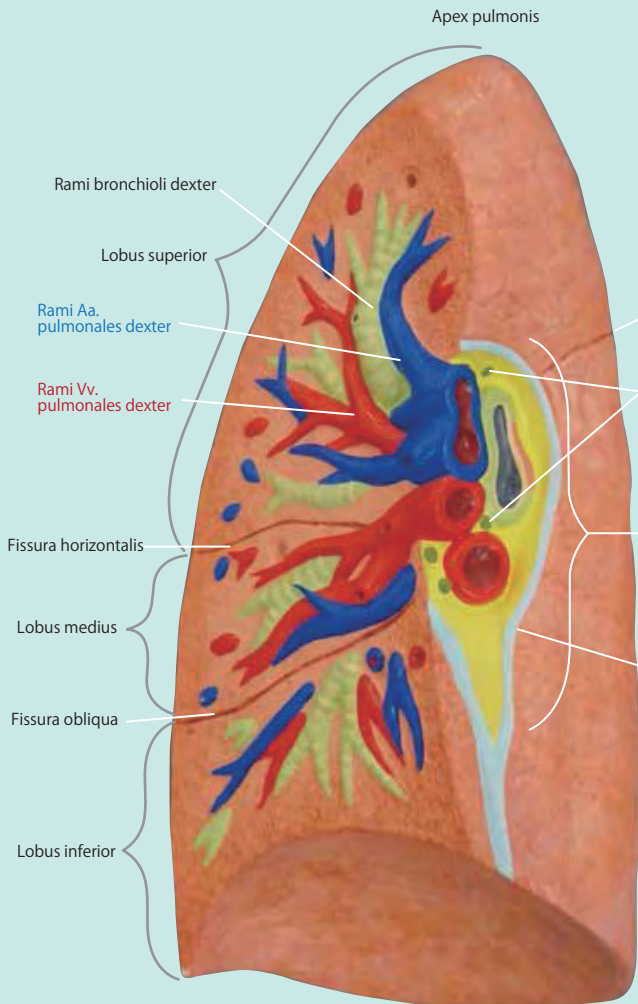
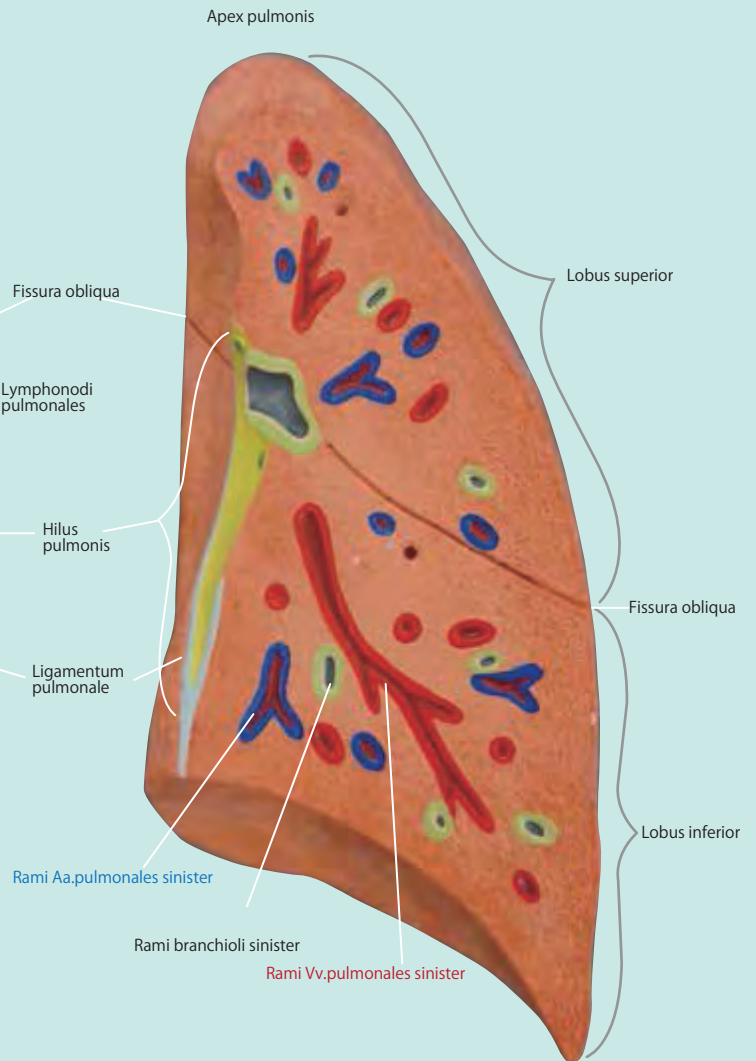
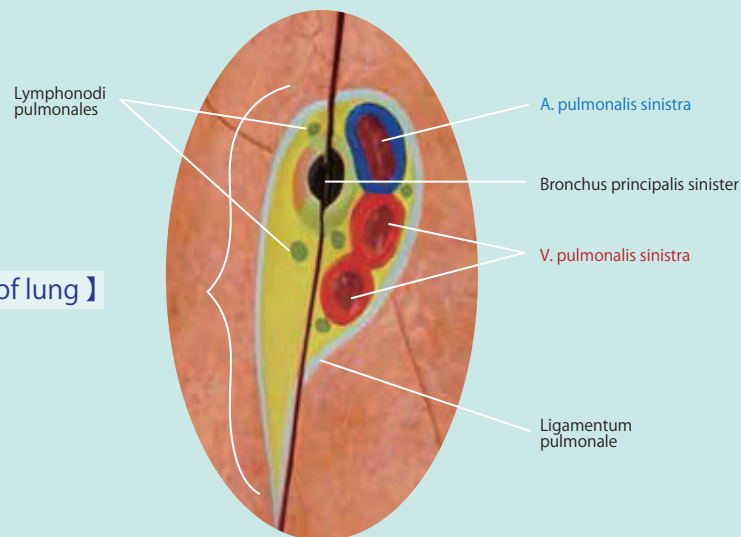
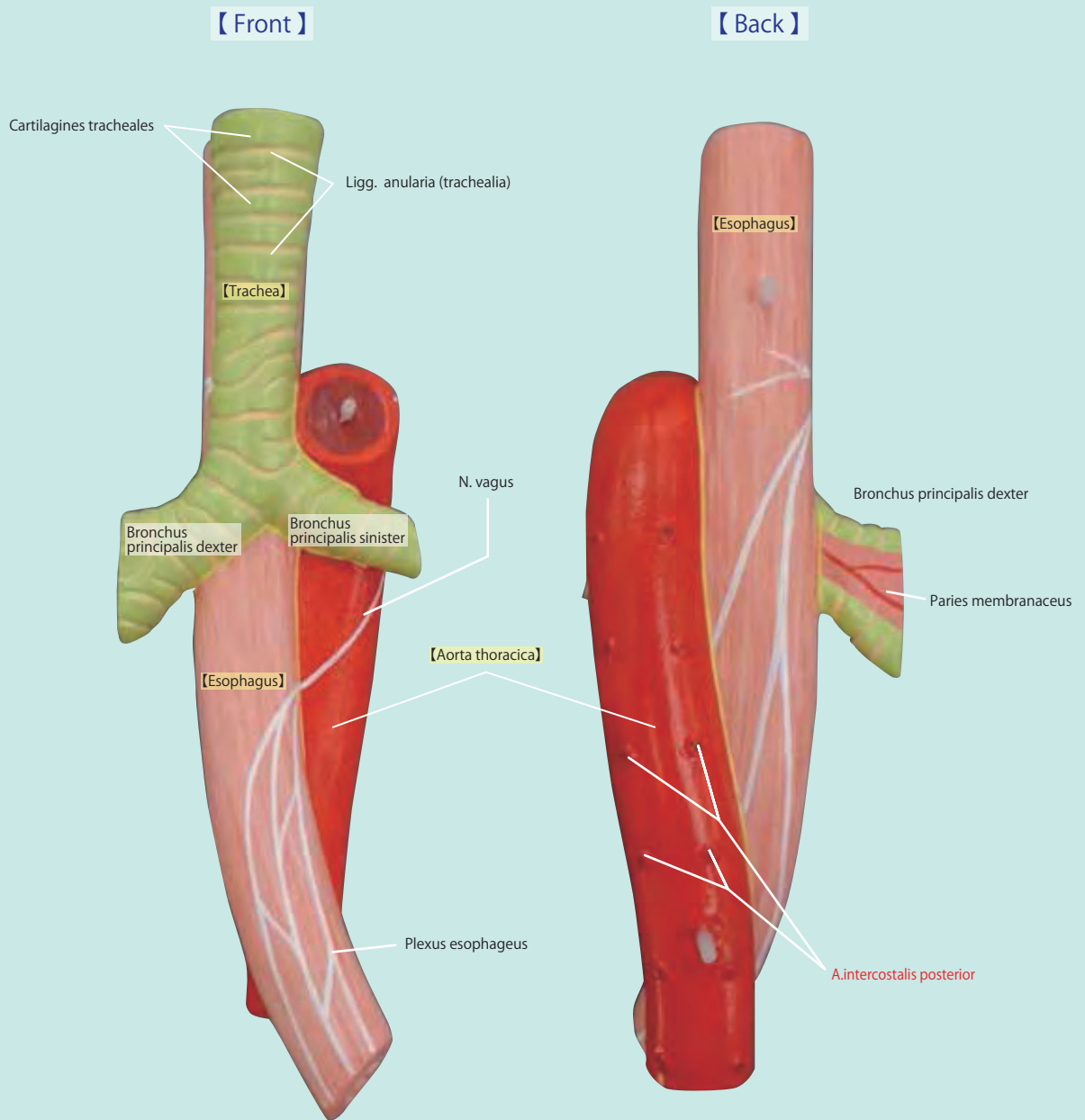
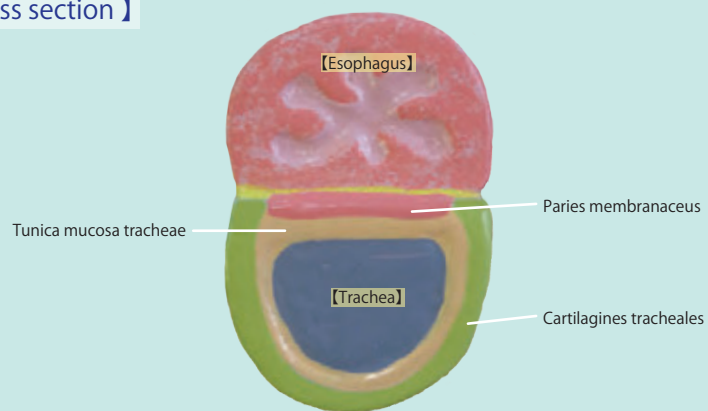
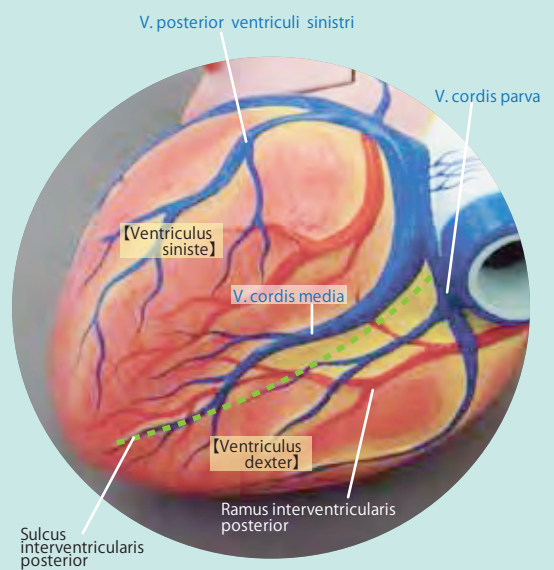
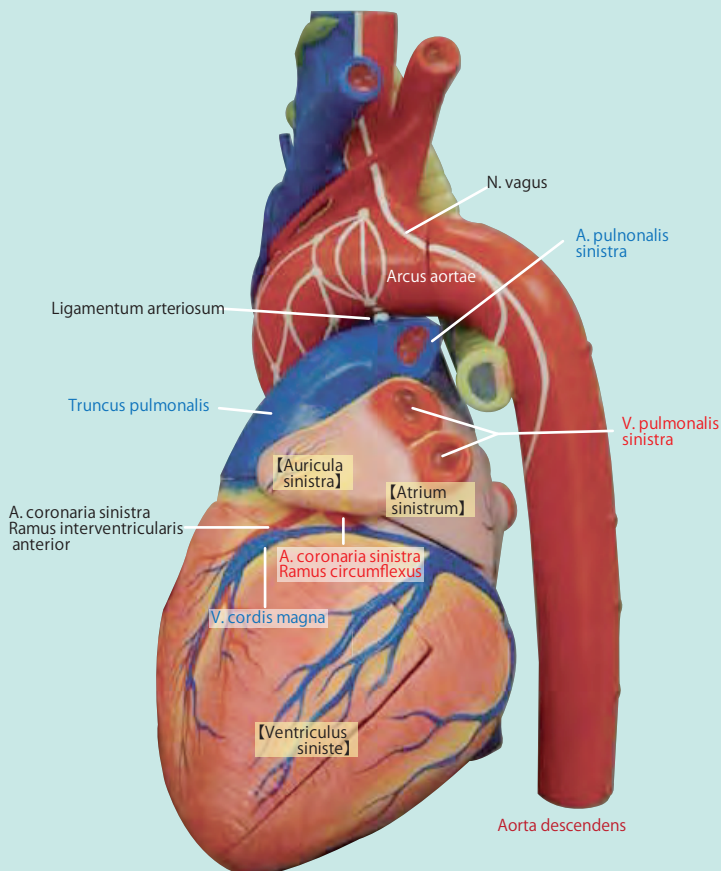
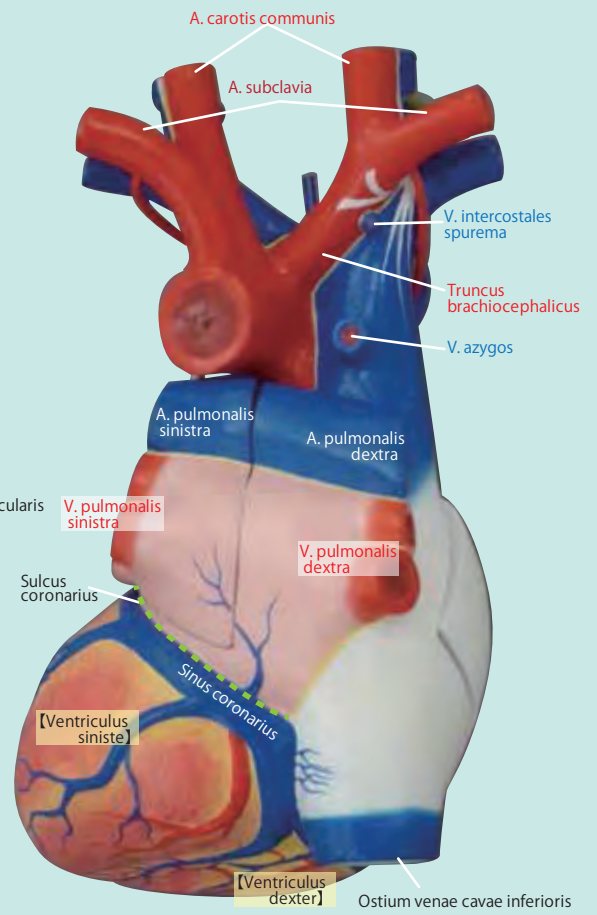
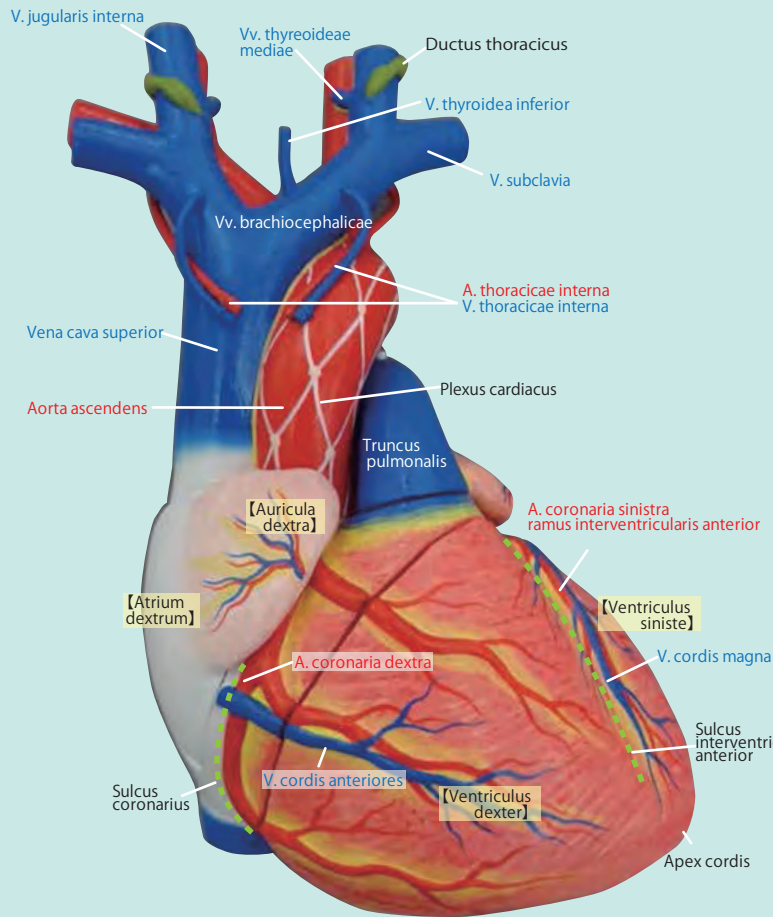
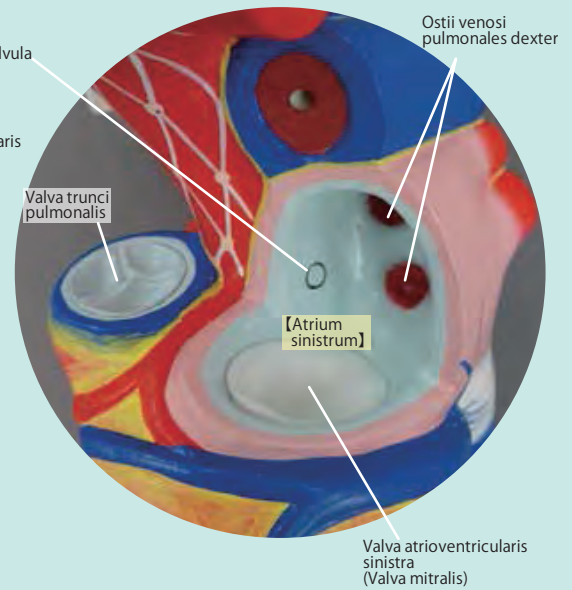
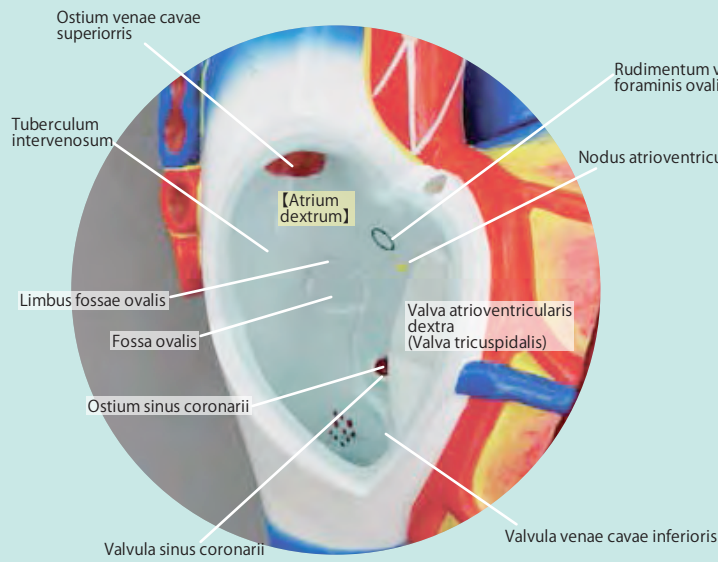
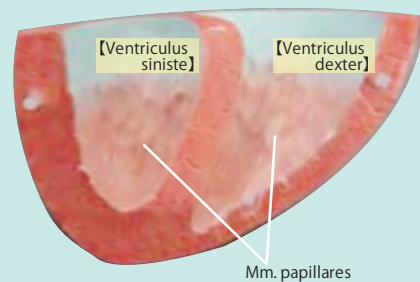
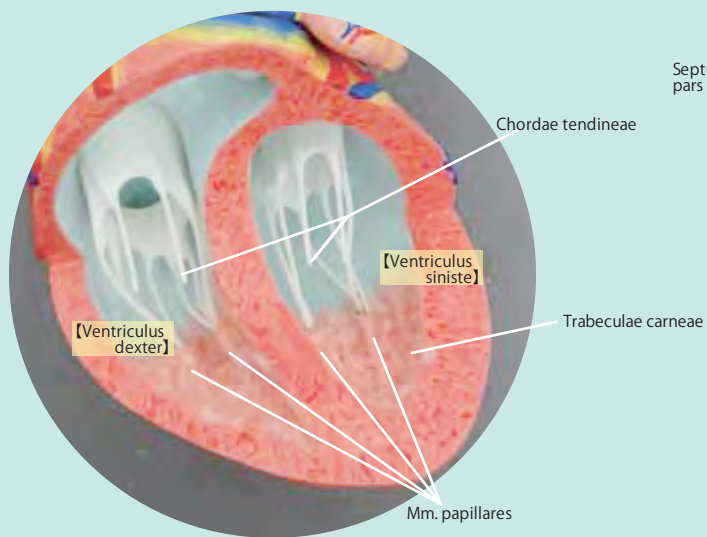
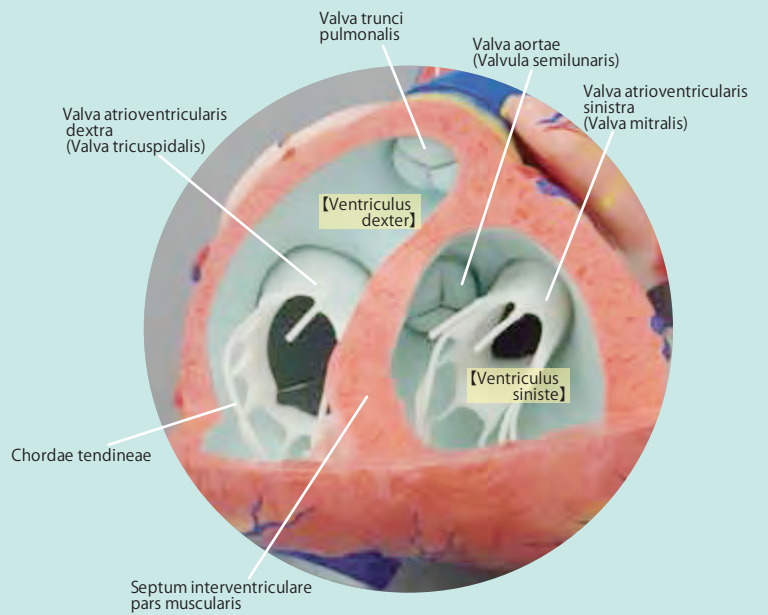
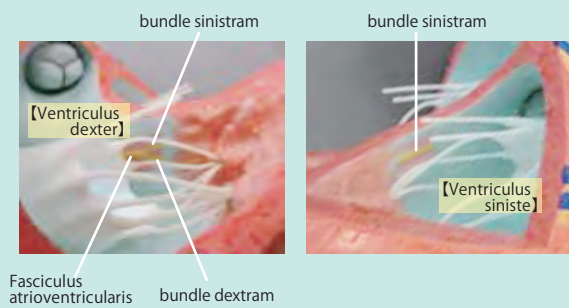
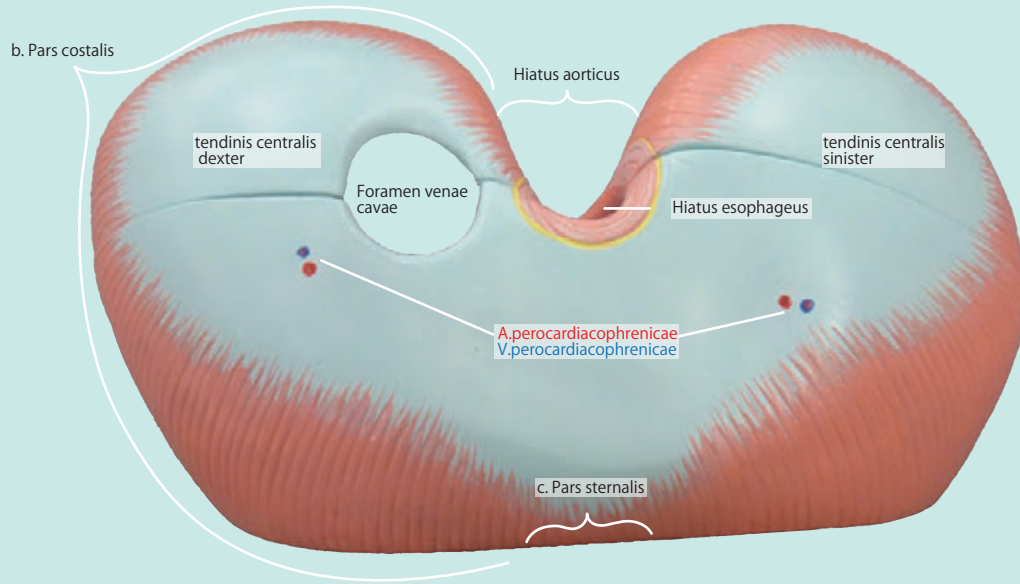
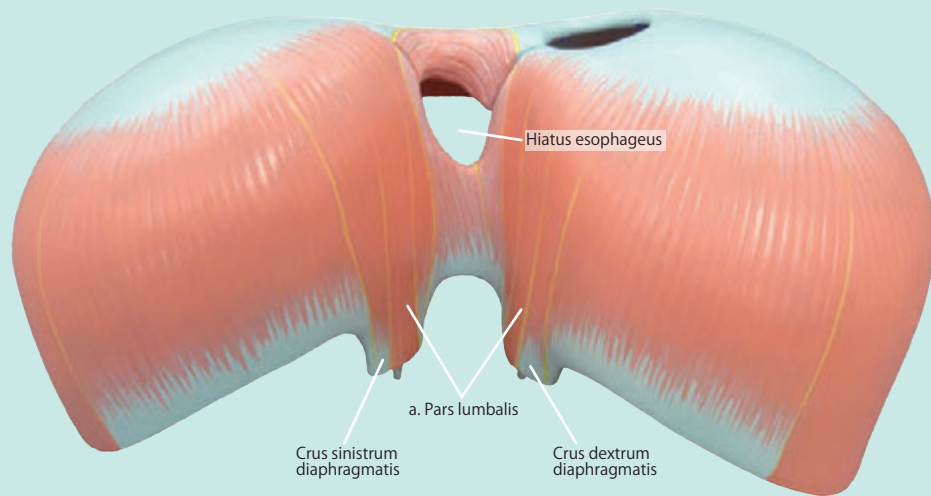
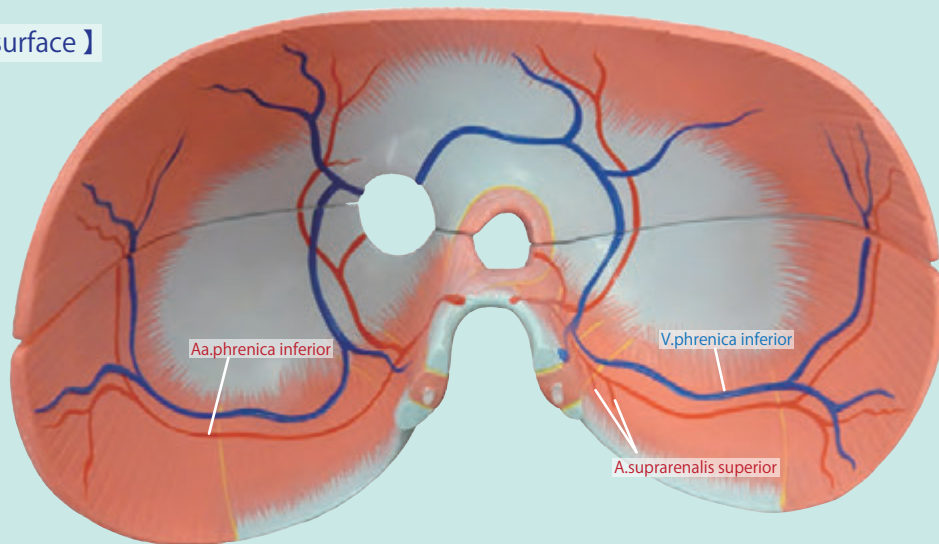
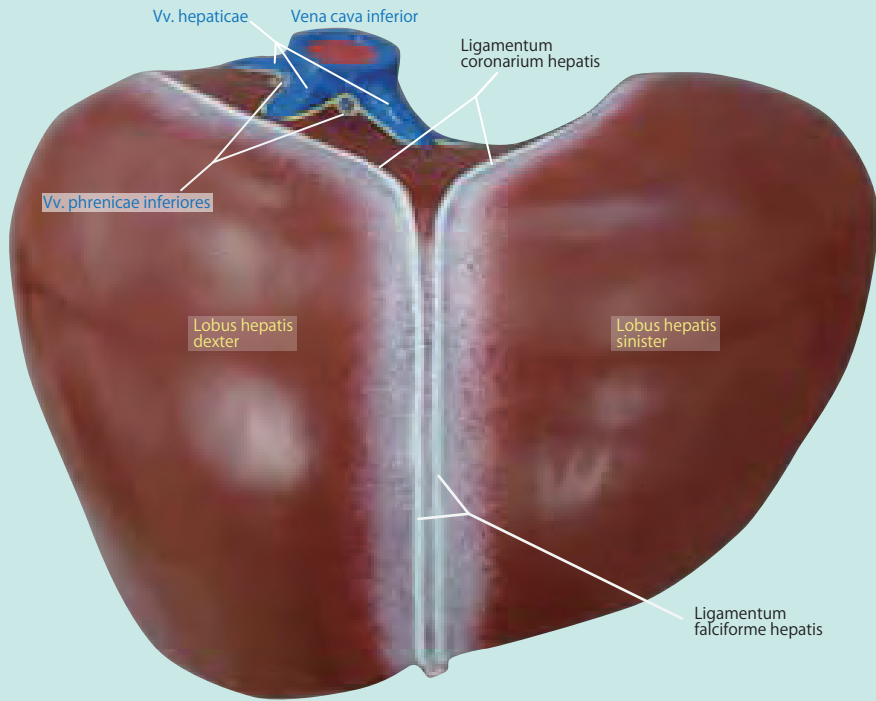
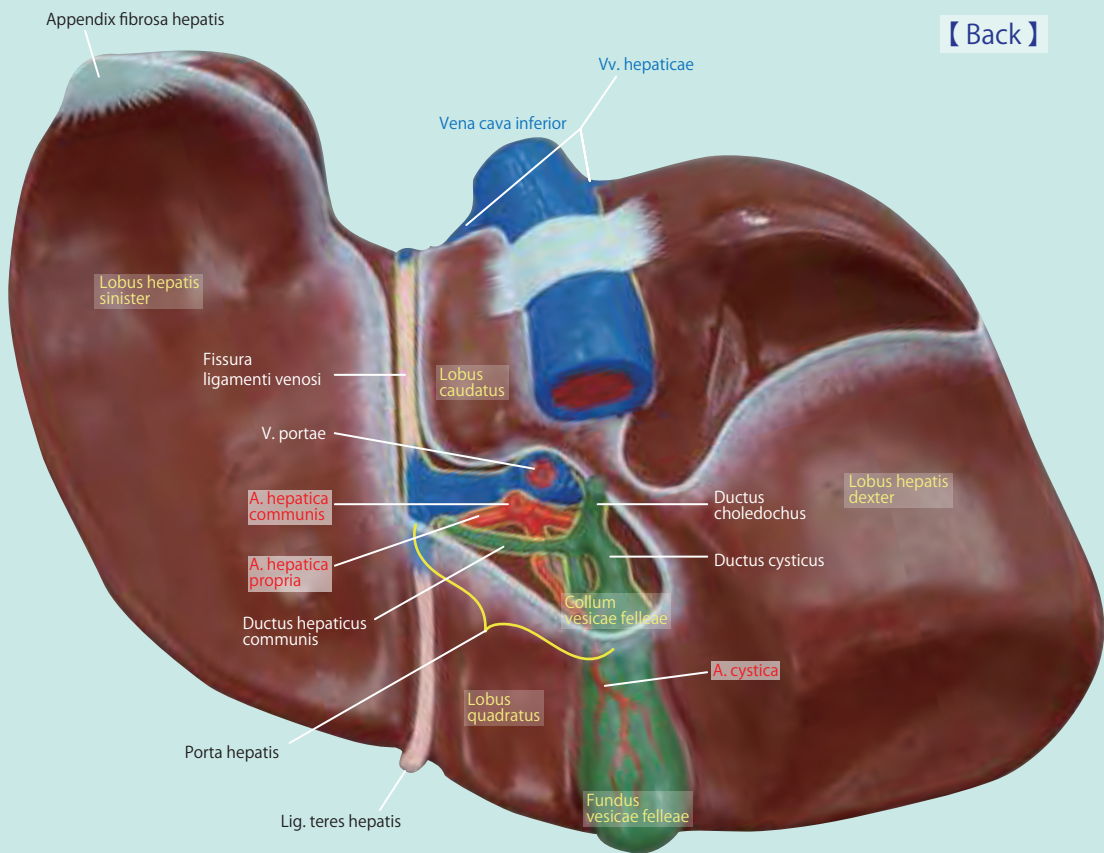
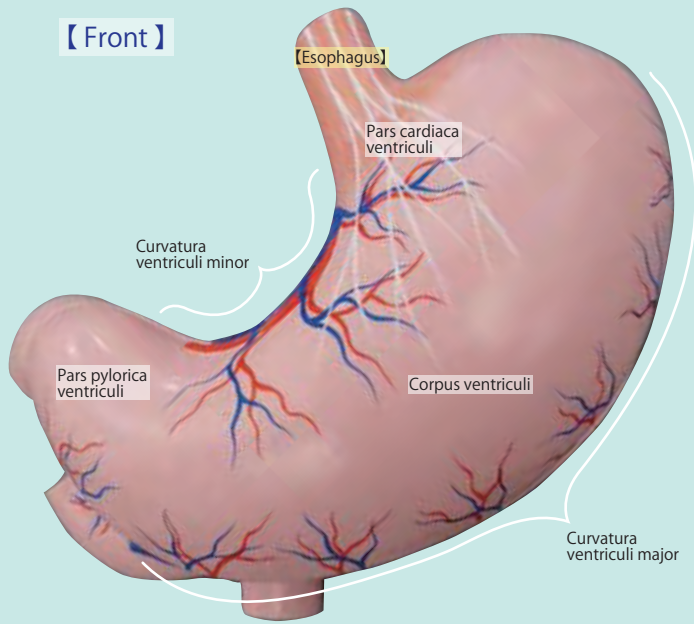
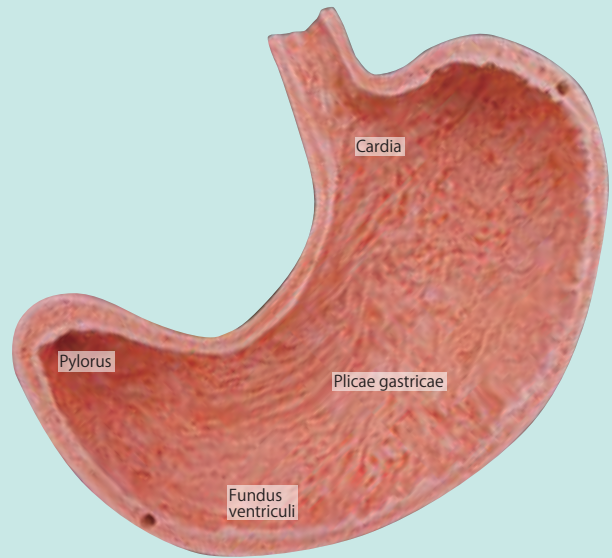
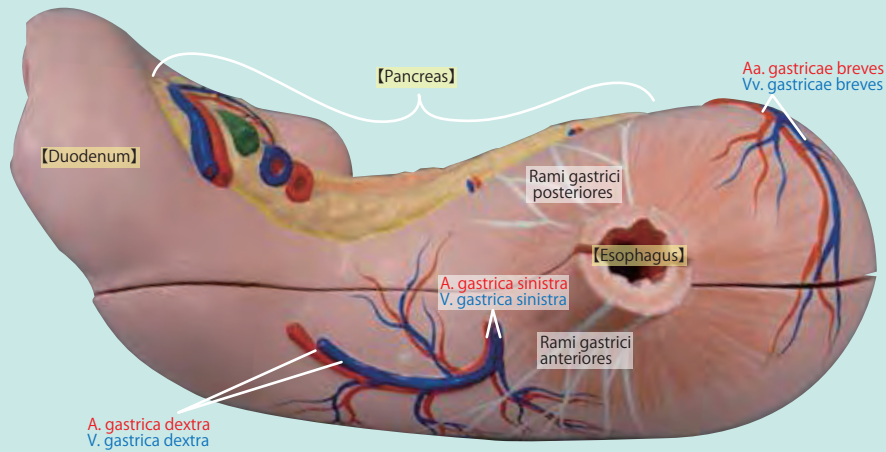
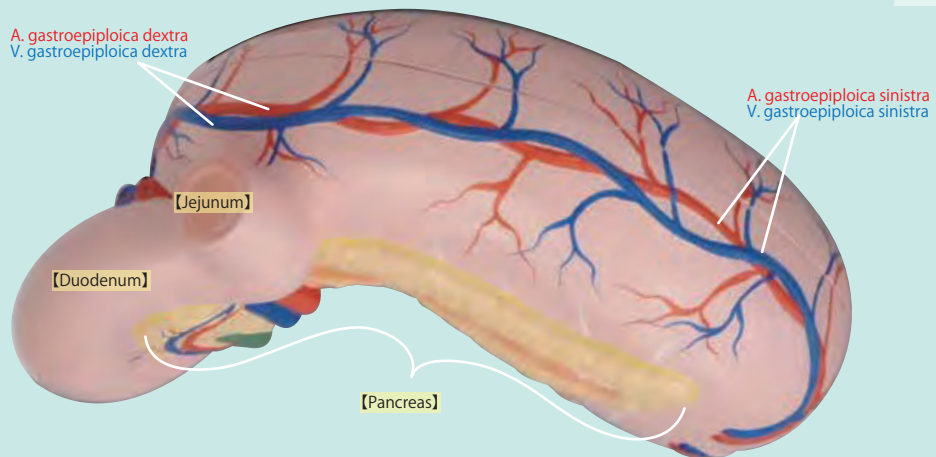
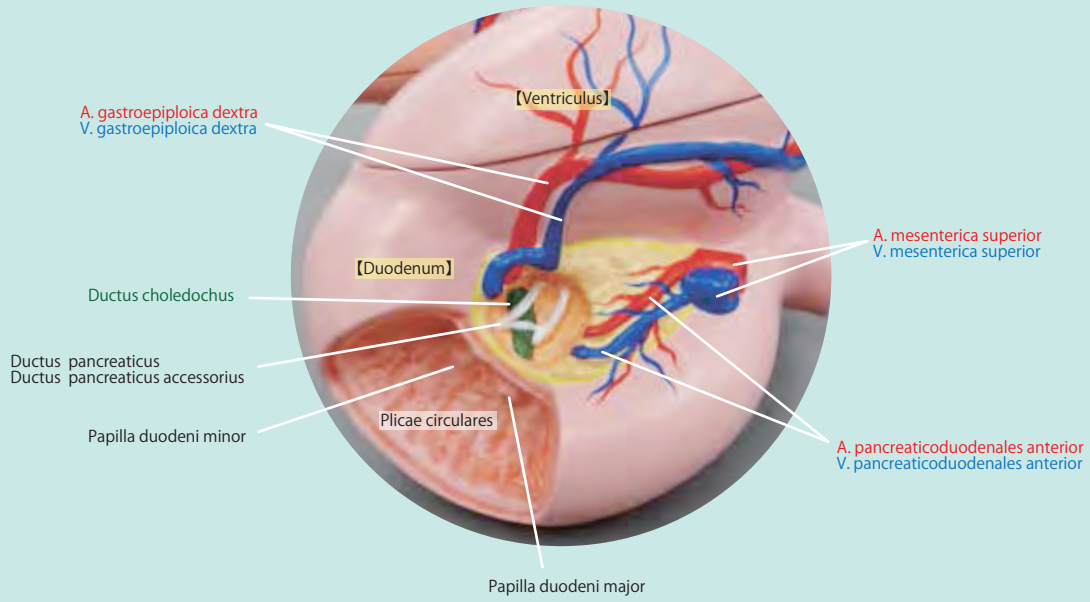
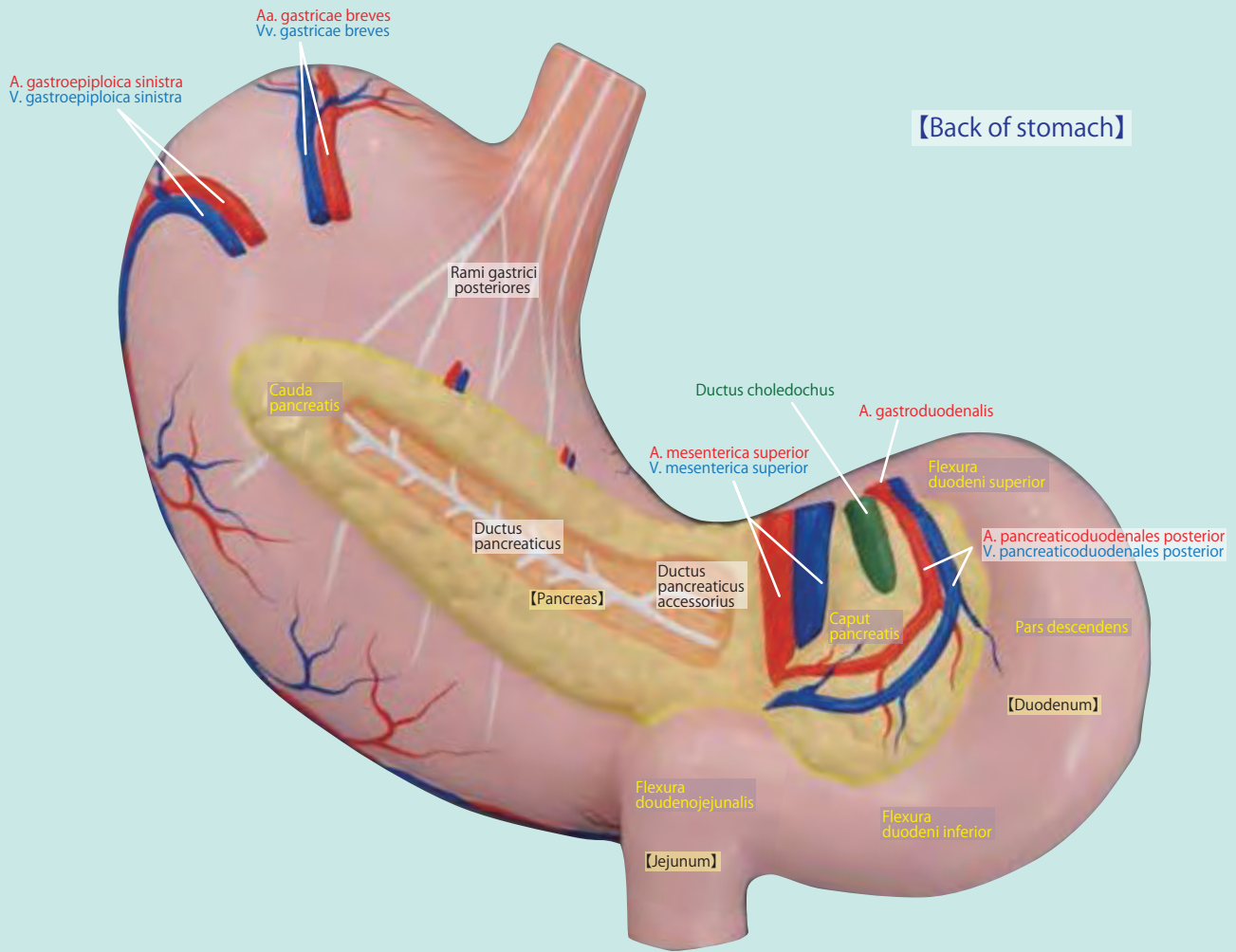
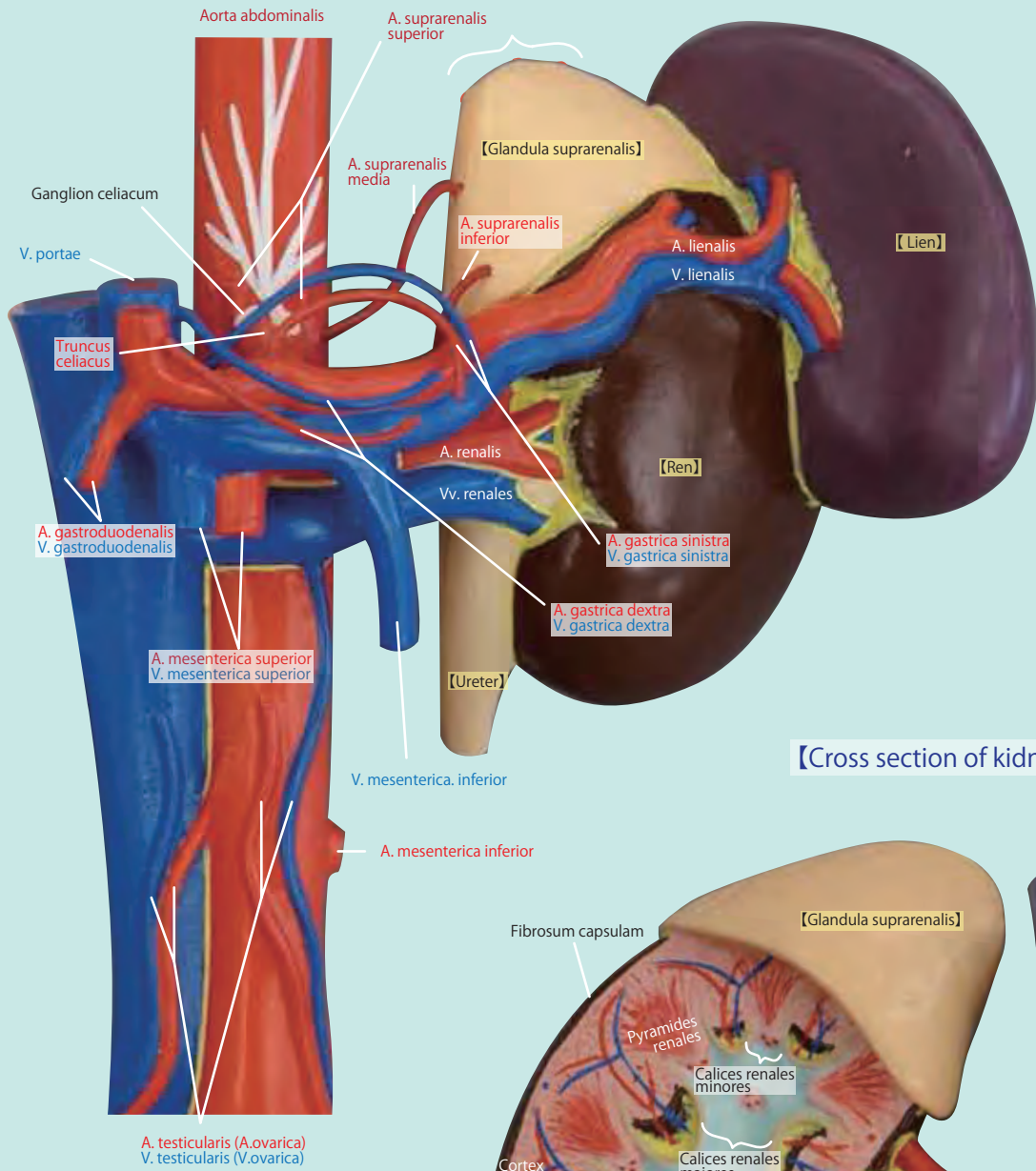
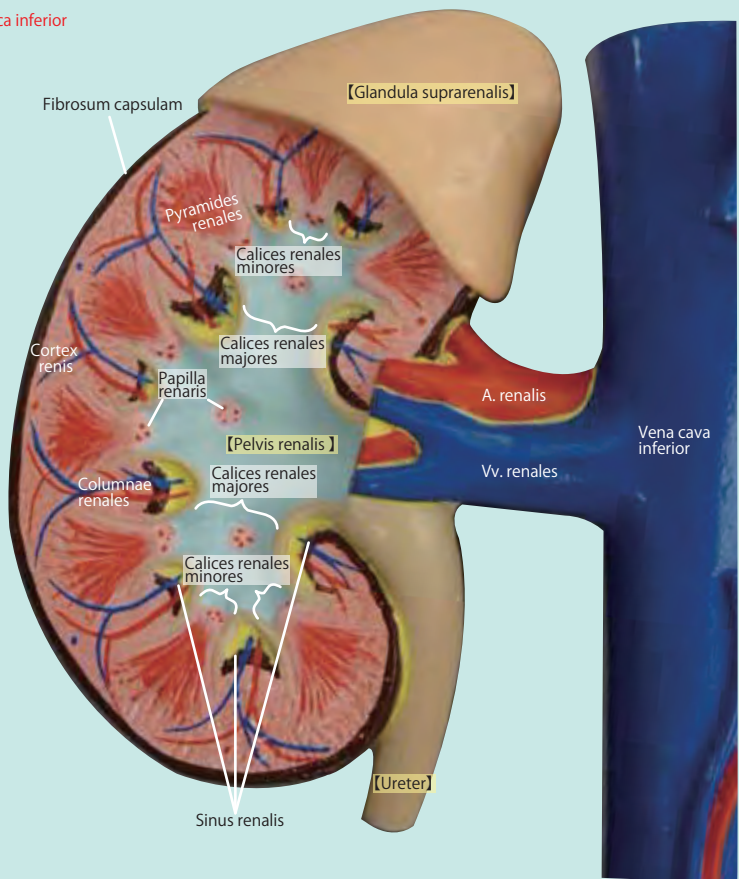
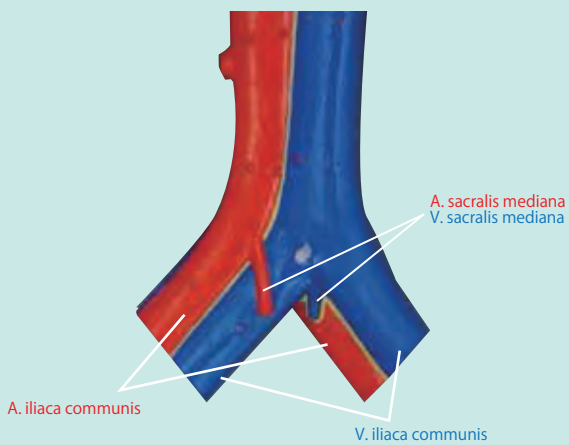
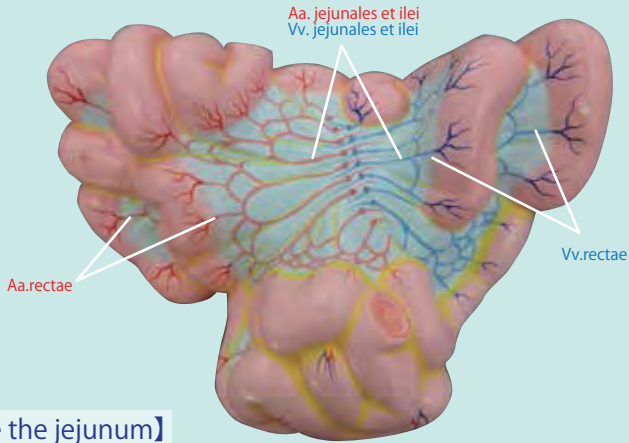
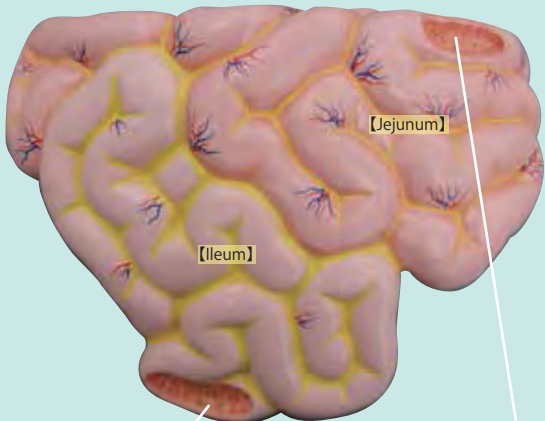
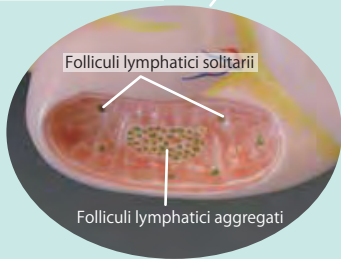
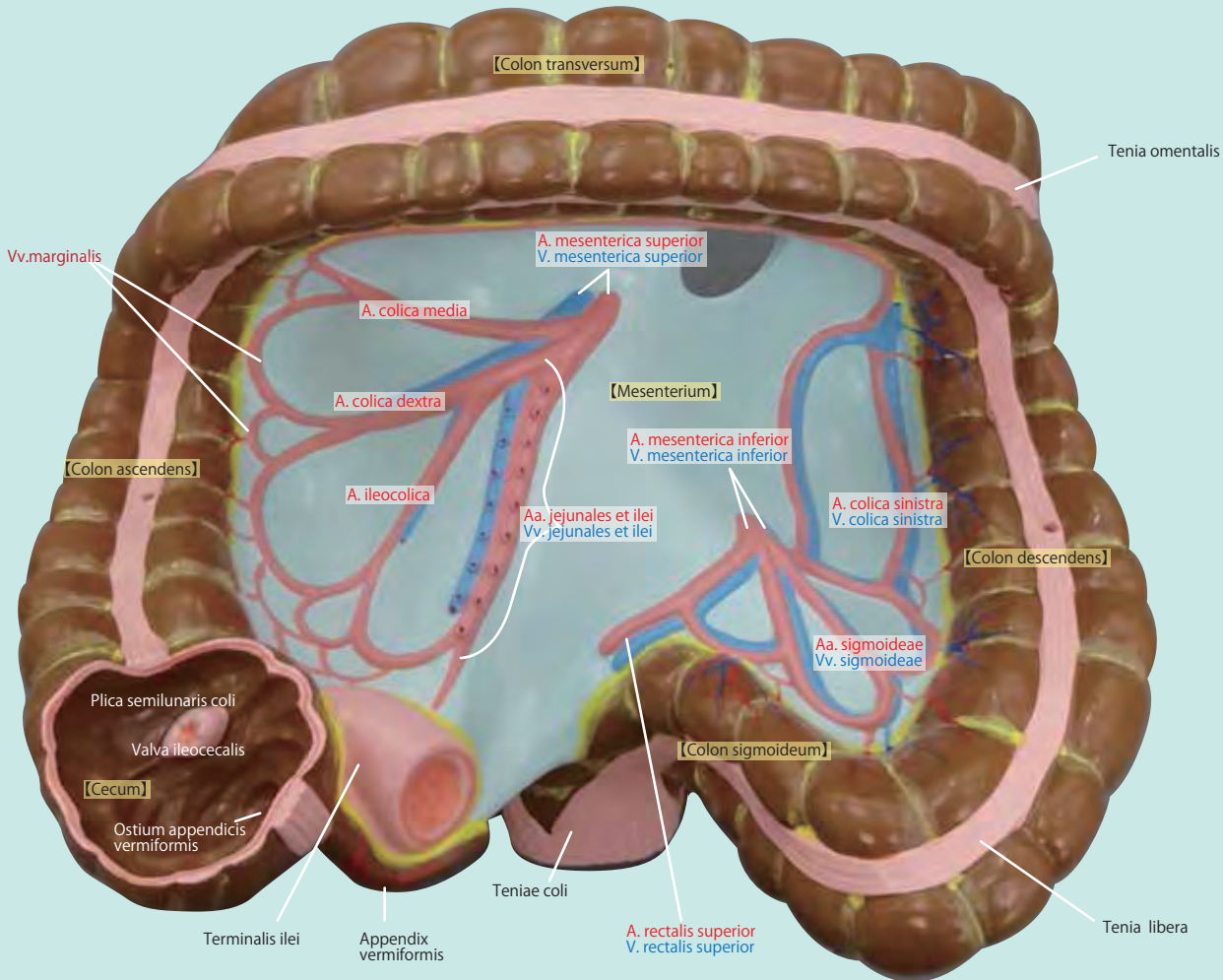
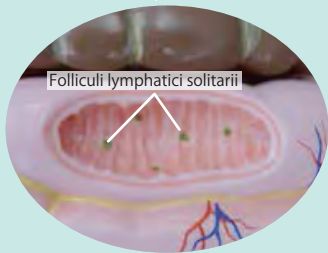
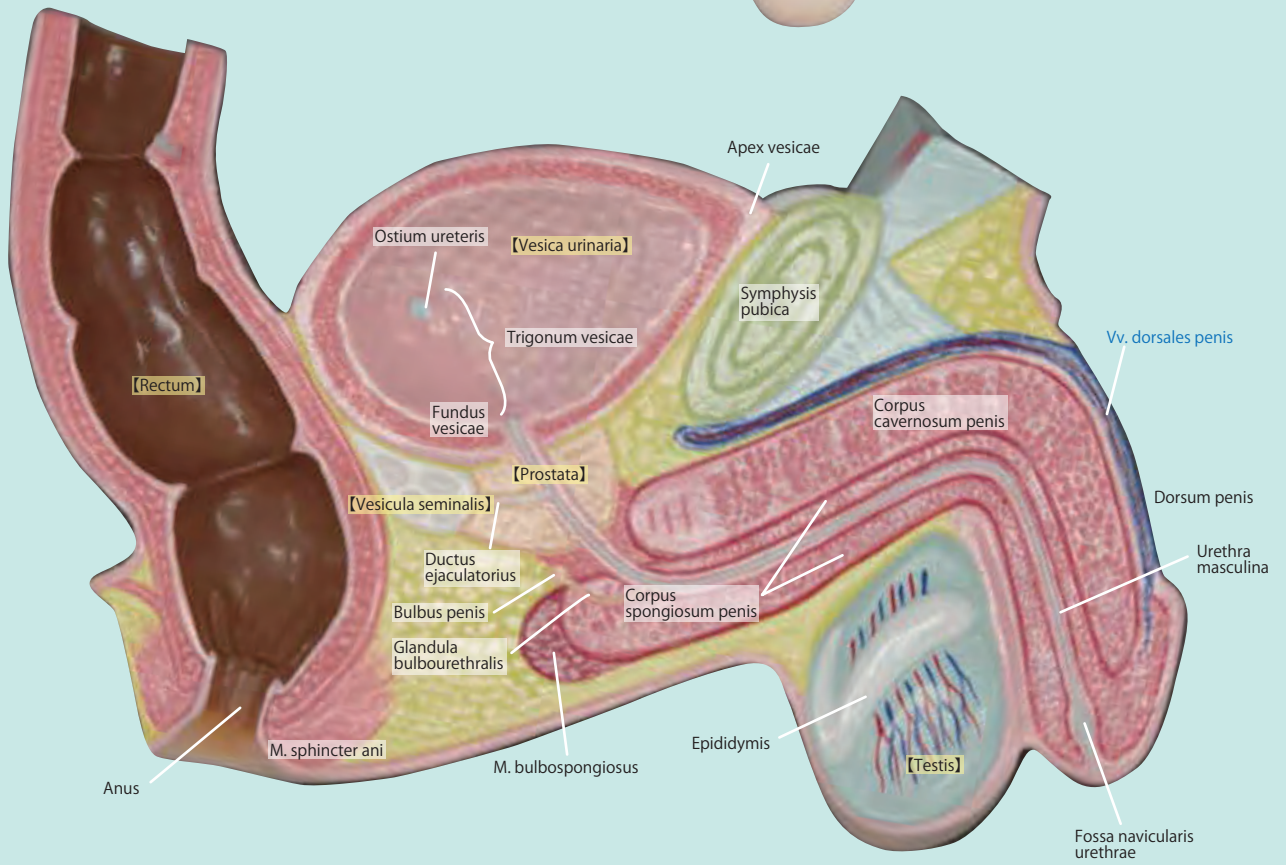
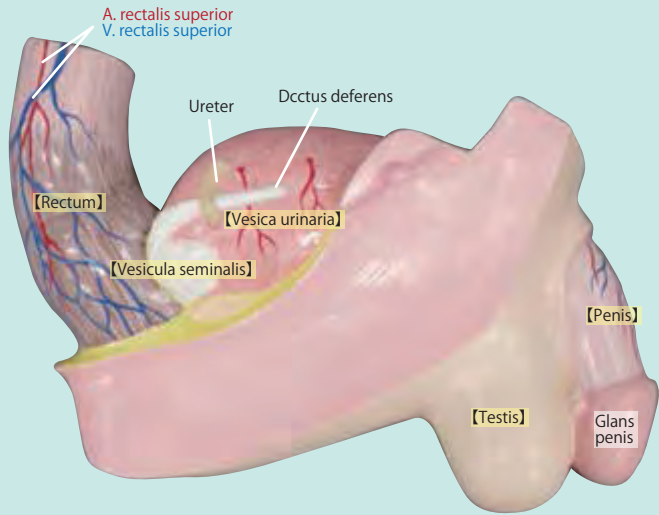
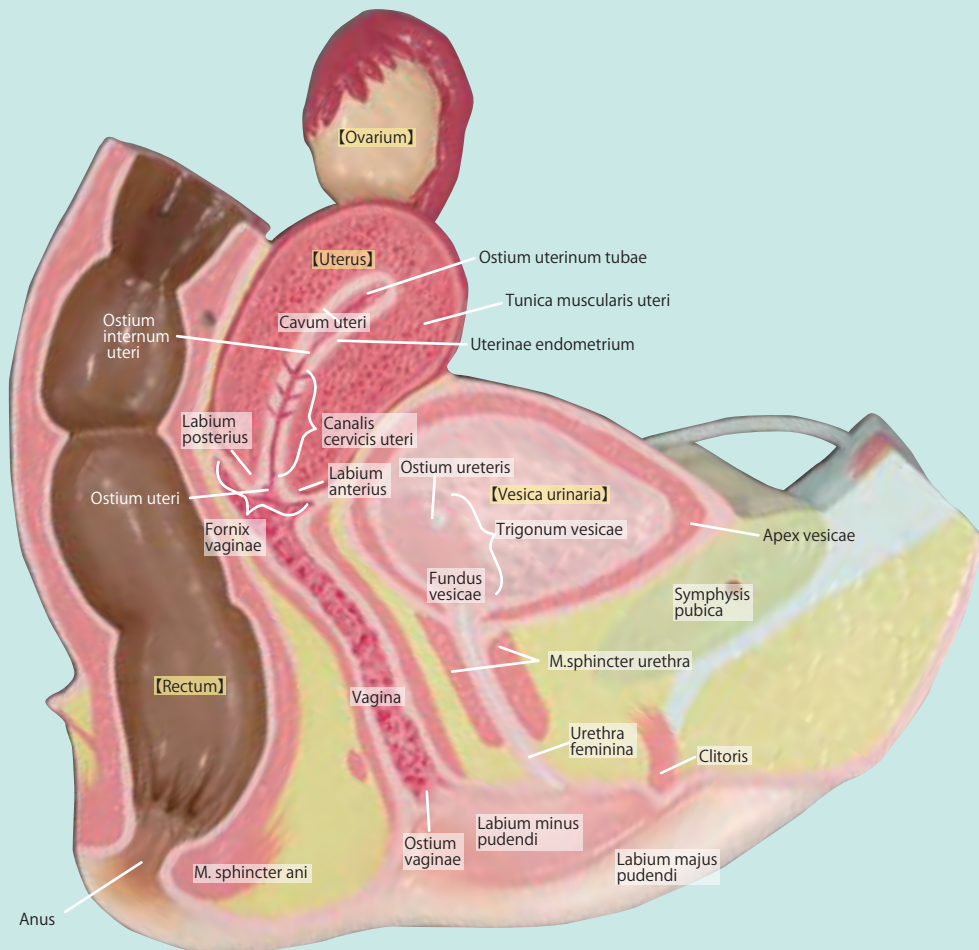
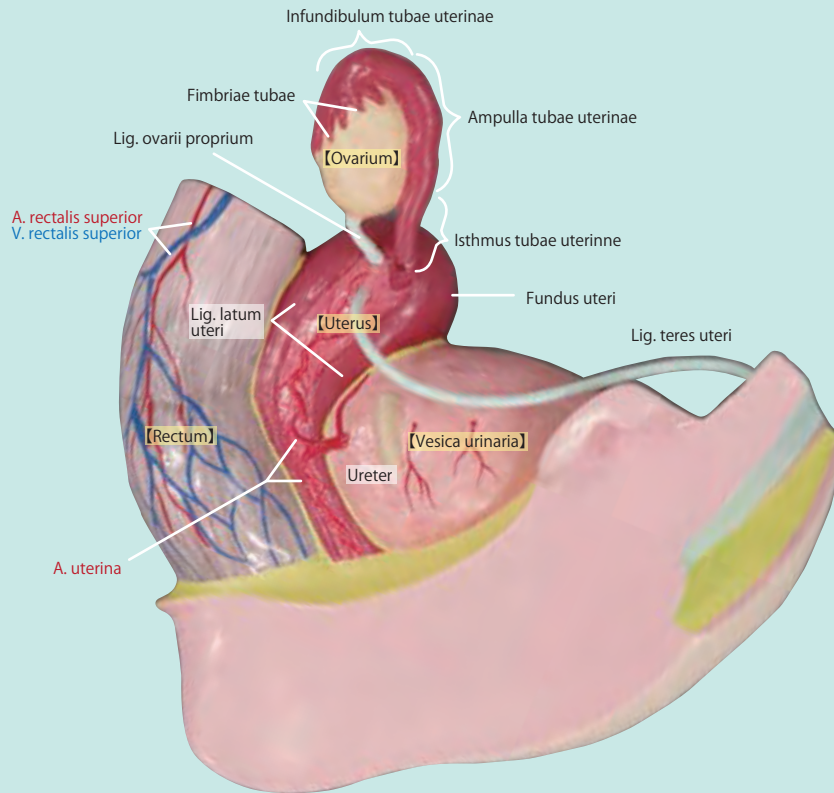
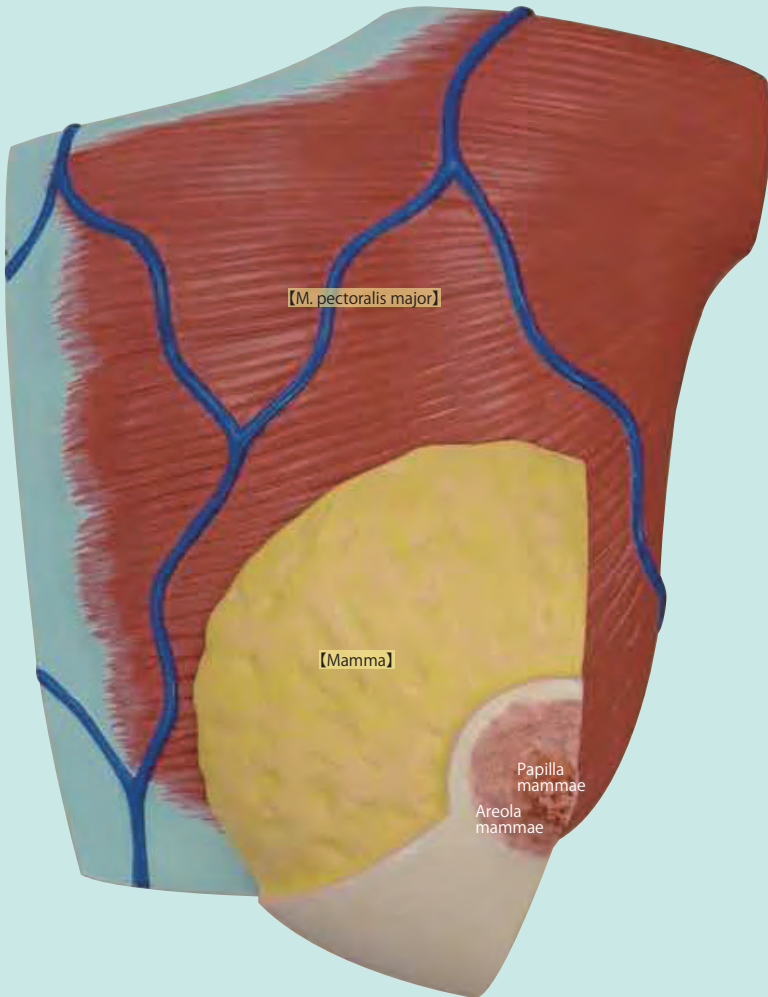
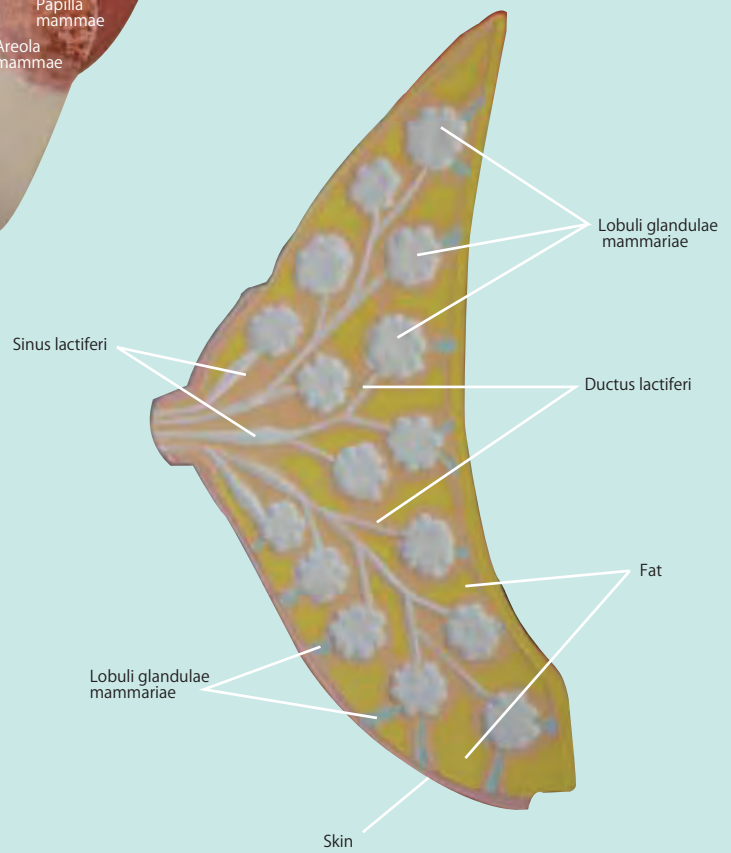
Caution

Do not mark on the model and other components with pen or leave any printed materials contacted on surface. Ink marks on the product cannot be removed.

For inquiries and service, please contact your distributor or **KYOTO KAGAKU CO., LTD.**

The contents of the instruction manual are subject to change without prior notice.

No part of this instruction manual may be reproduced or transmitted in any form without permission from the manufacturer. Please contact manufacturer for extra copies of this manual which may contain important updates and revisions.

Please contact manufacturer with any discrepancies in this manual or product feedback. Your cooperation is

WORLDWIDE INQUIRIES & ORDERING

 **KYOTO KAGAKU co.,LTD**

<http://www.kyotokagaku.com>
rw-kyoto@kyotokagaku.co.jp

● **KYOTO KAGAKU HEAD OFFICE & FACTORIES**

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